


version 7.1.0 rev 13 April 2016

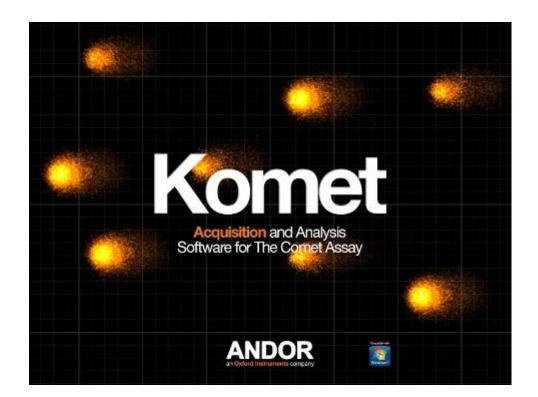

Software Guide



COPYRIGHT

Andor Technology 2005, 2016

Kinetic Imaging Ltd, 1992, 1993 ,1994, 1996, 1998,1999, 2000, 2001, 2002, 2003, 2004.

All rights reserved. No part of this publication may be reproduced, transmitted, or translated into any language in any form without the written permission of Andor Technology. Andor Technology is an Oxford Instruments company.

TRADEMARKS

Microsoft, Microsoft Windows, Excel and Word for Windows are trademarks of The Microsoft Corporation.

Andor Technology hereby reserves rights over the name Komet[®] for use with regard to the analysis of SCGE image data by digital image processing. The name Komet[®] has been reserved as a trademark in the United States of America.

This manual was correct at the time of printing, but may not include all developments of the product, due to our policy of continually upgrading and improving Komet and our other products.

Contact

For customer support please contact: www.andor.com/contact us/support request

Manual Revision History

08 Dec 2014	Amended information and updated presentation throughout for Komet 7.0.1 (all sections).
13 Apr 2016	Updated for Komet 7.1.0.



Contents

SECTION 1: Welcome to KOMET!
1.1 Computer Specifications
1.2 Installation of Software14
1.3 New Features in Komet15
1.3.1 New Imaging Interfaces and Digital Camera support15
1.3.2 Database Viewer
1.3.3 Printing Datasets in Database Viewer16
1.3.4 LED Light Source Recommendation17
1.4 Other Andor Products
SECTION 2: Komet and the Comet Assay18
2.1 An Imaging Protocol for 'KOMET' Analysis18
2.2 Background Effects
2.3 Comet Visibility, Electrophoresis Time and Tail Length19
2.4 Analogue CCD Imaging19
2.5 Digital CCD and sCMOS Cameras20
2.6 Hints on Microscopy21
2.7 Light Source, Lamp-house and Epi-fluorescence path22
2.7.1 Fluorescence Filter Set and Microscope Objective23
2.7.2 Beam-splitter and C-mount Adapter25
2.7.3 Summary
2.8 User Management and Logon Procedure26
2.8.1 Running Komet While Logged Onto a Domain27
2.8.2 Configuration Manager
2.9 Repeatability Measurements
2.10 The Benefits of Komet
2.10.1 Flexibility, Protocol Setup and Data Integrity29
2.10.2 Calibration - Simple and Infrequent29
2.10.3 Overlays – Region of Interests (ROIs) and Calipers
2.10.4 Dose Statistics in a Secure Access Database
2.10.5 Suspending and Resuming Experiments
2.10.6 Conclusion
SECTION 3: Getting Started



3.1 Starting Komet	.33
3.2 The Komet User Interface	.34
3.2.1 The Image Window	35
3.2.2 The Comet Toolbar	36
3.2.2.1 Live/Store	36
3.2.2.2 Interactive	36
3.2.2.3 Delete Last	36
3.2.2.4 Express	36
3.2.2.5 Experiment/Abort	37
3.2.2.6 Open	.37
3.2.3 The Key Results Panel	.37
3.2.4 The Komet Control Window	.38
3.3 Working with Live and Stored Images	.39
3.3.1 Live Images	.39
3.3.2 Stored Images	.40
3.3.3 Notes on Single versus Multi-Cell Scoring	.41
3.3.4 Free and Experiment Modes	.41
3.3.4.1 Free Mode	.41
3.3.4.2 Experiment Mode	43
3.3.5 Live Image Experiment Sequence	.44
3.3.6 Live Image Experiment Sequence With Database	.45
3.3.7 Stored Image Experiment Sequence	.46
3.3.8 Stored Image Experiment Sequence with Database	.47
3.3.9 Data Management	.48
3.4 Calibrating the System	.49
3.4.1 Calibration Hints	56
3.4.1.1 Haemocytometer	56
3.4.1.2 Graticule	56
3.4.1.3 Binning	57
3.5 How to Setup Comet Protocols	58
3.5.1 Creating a New Comet Protocol	59
3.5.2 Protocol Calibration Options	60
3.5.3 Camera Settings and Adjustment	.61
3.5.4 Palette or Look Up Tables (LUTs)	C 2



3.5.5 Comet Options Settings	65
3.5.5.1 Comet Options: Processing	65
3.5.5.2 Comet Options: Experimental	66
3.5.5.3 Comet Options: Operational	66
3.5.6 Editing a Comet Protocol	67
3.5.7 Comet Protocol File Management	68
3.5.8 Sharing Comet Protocol Files	68
3.6 Running an Experiment with a Camera	69
3.6.1 Starting an Experiment	70
3.6.2 Acquire an Image	71
3.6.3 Set the Cell / Background ROI	72
3.6.4 Scoring Data	73
3.6.5 Experiment Mode - User Prompts	74
3.6.6 Running an Experiment on Stored Images	75
3.6.7 Starting an Experiment	75
3.6.8 Set the Cell / Background ROI	77
3.6.9 Scoring Data	78
3.6.10 Experiment Mode - User Prompts	78
3.6.11 Running an Experiment Using a Database	79
3.7 Starting an Experiment	80
3.7.1 Acquire an Image	82
3.7.2 Set the Cell / Background ROI	82
3.7.3 Scoring Data	83
3.7.4 Experiment Mode - User Prompts	83
3.7.5 How to Suspend an Experiment	86
3.7.6 How to Resume an Experiment	87
SECTION 4: Komet Function Reference	89
4.1 The Image Window	89
4.2 The Image Window Menus	90
4.2.1 Image Window File Menu	90
4.2.1.1 File Menu: Open Image	90
4.2.1.2 File Menu: Save Image	92
4.2.1.3 File Menu: Save Display (BMP) / Save Display (JPEG)	93
4.2.1.4 File Menu: Print	94



4.2.1.5 File Menu: Exit	95
4.2.2 Image Window Edit Menu	95
4.2.3.1 Edit Menu: Copy to Clipboard	95
4.2.3.2 Edit Menu: Calibration	95
4.2.4 Image Window View Menu	96
4.2.4.1 View Menu: LUT	96
4.2.4.2 View Menu: Mapping	97
4.2.4.3 View Menu: AutoMap	99
4.2.4.4 View Menu: Toolbars	99
4.2.4.1 Toolbars Menu: Toolbar	100
4.2.4.2 Toolbars Menu: Status	
4.2.4.5 View Menu: Image Background	
4.2.5 Image Window Image List Menu	
4.2.6 The Comet Toolbar	
4.2.6.1 Live/Store	
4.2.6.2 Interactive	
4.2.6.3 Delete Last	
4.2.6.4 Express	
4.2.6.5 Experiment/Abort	
4.2.6.6 Open	
4.3 The Key Results Panel	
4.4 The Komet Control Window	
4.5 The Komet Control Window Menus	
4.5.1 Komet Control Window File Menu	
4.5.1.1 File Menu: Save Image	
4.5.1.2 File Menu: Set Save Path	
4.5.1.3 File Menu: Settings	
4.5.1.4 File Menu: Exit	
4.6 Komet Control Window QC Menu	
4.7 Komet Control Window Wizards Menu	
4.7.1 Wizards Menu: Calibration	
4.7.2 Wizards Menu: Comet Protocol	111
4.8 Komet Control Window Help Menu	112
4.8.1 Help Menu: Help	112



4.8.2 Help Menu: Product Activation
4.8.3 Help Menu: About
4.9 The Current Context Panel114
4.10 The Analyse Tab
4.10.1 Intensity Profiles
4.10.2 Individual Analysis Parameters
4.10.3 Comet Protocol Tab
4.11 Comet Options Settings
4.11.1 Comet Options: Processing
4.11.1 .1 Head Threshold %
4.11.1.2 Smoothing Value
4.11.1.3 Tail Break Length121
4.11.1.4 Tail Threshold/1000
4.11.1.5 Maximum Head Diameter
4.11.1.6 Background Height122
4.11.1.7 Saturation Limit
4.11.1.8 Saturation Maximum
4.12 Comet Options: Experimental124
4.12.1 Cells per Slide
4.12.2 Replicate Slides
4.12.3 Samples
4.12.4 Retain Images in Memory
4.12.5 Use Automatic Data Directories125
4.12.6 Use Database
4.13 Comet Options: Operational
4.13.1 Fluorescence
4.13.2 Auto Background
4.13.3 AutoExpress on Store
4.13.4 Stop on Saturation
4.13.5 Floating Live Button
4.13.6 Head Ellipse On
4.13.7 Maximum Head Diameter129
4.13.8 Head Symmetry129
4.13.9 Tail From Centre of Cell



4.13.10 Interactive Tail
4.13.11 Floating Open Button130
4.13.12 Flip On Open
4.13.13 Highlight Scored Cells
4.14 Komet Option Switches Tab131
SECTION 5 Concepts in Komet GLP A Reference Guide
5.1 Introduction
5.1.1 Managing Validated Komet GLP workstations133
5.2 Continuity of Core Algorithms with QC Checking
5.3 Other Services
5.4 Core Concepts and Terminology134
5.4.1 Two-level Logon and Multi-mode Access
5.4.2 Studies, Datasets, Scorers and Protocols135
SECTION 6 Administrator Functions
6.1 Creating Users in KGLP – Windows Access Issues
6.2 System Configuration by the Local Machine Administrator (LMA)137
6.3 Running KGLP in System Setup Mode - Local Machine Administrator (LMA)137
6.4 Security Menu
6.4.1 Personnel Access Control
0.4.1 Personner Access control
6.4.2 Timeout Delay provides a lockout when KGLP is left open
6.4.2 Timeout Delay provides a lockout when KGLP is left open141
6.4.2 Timeout Delay provides a lockout when KGLP is left open
6.4.2 Timeout Delay provides a lockout when KGLP is left open
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset143
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password143
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware143
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware1436.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts144
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware1436.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts1446.6 View Audit Trails Menu tracks Local Machine Administrator and System QC activity146
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware1436.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts1446.6 View Audit Trails Menu tracks Local Machine Administrator and System QC activity147
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware1436.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts1446.6 View Audit Trails Menu tracks Local Machine Administrator and System QC activity147SECTION 7 Study Manager Functions148
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware1436.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts1446.6 View Audit Trails Menu tracks Local Machine Administrator and System QC activity1466.7 Wizards Menu147SECTION 7 Study Manager Functions1487.1 Studies, Datasets, Scorers and Protocols148
6.4.2 Timeout Delay provides a lockout when KGLP is left open1416.5 Study Password Change Frequency1426.5.1 Study Password Reset1426.5.2 Dataset Password Reset1436.5.3 Change Setup Password1436.5.4 Setting up Hardware1436.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts1446.6 View Audit Trails Menu tracks Local Machine Administrator and System QC activity1466.7 Wizards Menu147SECTION 7 Study Manager Functions1487.1 Studies, Datasets, Scorers and Protocols1487.2 Create New Study – Name, Password Entry – Use Template Option149



7.6 Review and/or Edit Established Datasets
7.7 Adding New Scoring Protocols – Protocol Name, Calibration, Camera and Comet Options 154
7.8 Use Existing Study - Study Manager
7.9 The Study Menu - Existing Studies
7.10 Close Study
7.11 E-Signatures and Dataset Integrity161
7.12 Check Dataset
7.13 Scoring Datasets – Study Managers
7.14 On Completion of a Dataset163
7.15 Review Closed Studies Mode
7.16 View Audit Trails
7.17 Reopen the Study166
7.18 Write Dataset to Standalone CD
SECTION 8 KOMET GLP for Scorers
8.1 Preparing to Score Datasets in Studies
8.2 Scoring Datasets
8.3 Key Results and Current Context
8.4 When the Current Slide is Complete175
8.5 Image Gallery for Current Sample and Printing177
8.6 Deleting Last and Interactive Scoring
8.7 Suspending the Scoring Session
8.8 Resume Scoring of a Dataset
8.9 Writing Stand Alone CD - On Completion of a Dataset182
PROTOCOL FOR THE APPLICATION OF THE pH>13 ALKALINE COMET ASSAY TO THE DETECTION
OF DNA DAMAGE IN MAMMALIAN CELLS
SECTION 9 Comet Assay184
9.1 Preparation of Reagents
9.2 Preparation of Slides for the Comet Assay
9.3 Cell Isolation
9.4 Electrophoresis of Microgel Slides191
9.5 Evaluation of DNA Damage
9.6 Model Experiment: Human Leukocytes and Hydrogen Peroxide194
APPENDIX 1 INSTALLATION OF SOFTWARE
A1.1 Evaluation license



A1.2 Installation of Sentinel Drivers197
A1.3 Initializing your Dongle with an Upgrade Code198
A1.3 Installing your Komet Software199
APPENDIX 2 Hardware Devices
A2.1 Camera Control Interfaces
A2.2 Virtual Camera and Analogue CCD Cameras205
A2.4 QImaging Cameras
A2.4.1 The QICam
A2.4.2 The Device Tab211
A2.4.3 The Camera Tab212
APPENDIX 3 Configuration Manager
A3.1 Run Configuration manager216
A3.2 Configuration Manager Menus217
A3.2.1 Settings Menu
A3.2.1.1 Settings Menu: Global Configuration Settings217
A3.2.1.2 Multi-User Access
A3.2.1.3 Image Disk Configuration
A3.2.2 Settings Menu: Software Protection Device (Dongle)220
A3.2.3 File and Edit Menus
A3.2.4 Tools Menu
A3.2.4.1 Tools Menu: Register/Unregister Any ActiveX Component
A3.2.4.2 Tools Menu: Register/Unregister Plugin Components
A3.2.4.3 Tools Menu: File List Generator
A3.2.5 About Menu
A3.2.6 Exit Menu
A3.3 Create a New Configuration
A3.4 Open Configuration
A3.5.1 Edit an Existing Configuration231
A3.5.2 Deleting an Existing Configuration
A3.6 Configuration Manager Summary234
APPENDIX 4 Komet Measurement Definitions
A4.1 Integrated Intensity Profiles and their Analysis235
A4.2 Komet Analysis - A Mathematical Description238
A4.2.1 Head and Tail Discrimination239



A4.2.2 Ensuring Intensity Profile Equivalence
A4.3 Komet Measurements and Parameter Definition242
A4.4 Comet Parameters
A4.5 Head Parameters
A4.6 Tail Parameters
A4.7 Global Comet Parameters
A4.8 Conclusions
APPENDIX 5 Komet Repeatibility Measurements
A5.1 Measurements Made248
A5.2 Summary of Results249
APPENDIX 6 Troubleshooting
A6.1 GENERAL PROBLEMS
A6.1.1 No Image appears when I click on Live254
A6.1.2 No Image appears when I select On-CCD
A6.1.3 Driver Problems
A6.1.4 User Name System Set-up Lost255
A6.1.5 Saturation Detection with the Meteor or Meteor 2255
A6.1.6 Regions Resizing and Disappearing



SECTION 1: WELCOME TO KOMET!

Welcome to the world of Andor Technology's scientific analytical imaging, and in particular our unique Comet Assay Software. Komet has been developed to image and analyse the samples produced using the Single Cell Gel Electrophoresis (SCGE) Assay, first developed by Singh *et al.* (1988). There is a laboratory protocol for the SCGE Assay included in this manual (Section 6), with permission from Ray Tice of Research Triangle Park, North Carolina. Andor has a keen interest in the development of the assay and offers services which include distribution of publication lists and the Comet Newsletter.

Note: Kinetic Imaging was acquired by Andor Technology in 2004. Andor Technology became an Oxford Instruments company in 2014.

We are delighted that you have purchased Komet, and hope it will prove both powerful and easy to use. This manual is designed to provide you with descriptions and discussions of the following:

- 1) Komet analysis functions and measurement definitions.
- 2) Komet user interface.
- 3) The considerations you should take into account when matching laboratory protocol to image analysis.

In addition you will find a user guide specifically for the **Andor Database Viewer** product which is now supplied with Komet and allows for flexible and powerful review, presentation, summarization and distribution of data from Komet studies.

Should you have a need to perform GLP studies, there is a GLP version of this popular software, which provides audit trails, and extensive functionality aimed at meeting FDA 21 CFR part 11 compliance. Section 5 of this user guide describes GLP features and configuration.



1.1 Computer Specifications

Komet requires the following minimum computer configuration:

- Windows 7 Professional
- Minimum Ram 2GB Mb, recommended 8GB
- 1TB Hard Disk (Solid State Disk preferred)
- Graphics adapter with a minimum resolution of 1280 x 1024 and 16K colours.
- A suitable display. We recommend at least a High Definition 24" monitor with resolution set at 1920 x 1080.
- Mouse or other suitable pointing device.

A number of sensitive Charge Coupled Device (CCD) and Zyla sCMOS cameras are supported, which will require installation of specific drivers and in some cases installation of hardware.

To run the full version of Komet, you will need to plug a USB software protection key (dongle) into your PC.

We are always pleased to hear from users with information on their applications, suggestions and ideas. Please feel free to contact us:

www.andor.com/contact_us/support_request



1.2 Installation of Software

Insert the Andor Installation CD-ROM and run the program **Start.exe** from the root folder. This will start the installation program. Select **Perform a Full Installation of Komet** and follow the steps below:

- 1. Andor Software Protection Key drivers for Windows 7. (Appendix A1.2)
- 2. Dongle should be supplied already initialized for Komet Software. (Appendix A1.2)
- 3. Camera and acquisition interface software: Andor Camera drivers (refer to the supplied installation documentation) include three models which support almost any type of camera. You will need to make sure you follow instructions to install this software, for example:
 - Andor cameras are typically operate via USB2 or USB3 interfaces
 - Video CCD cameras (previously via Matrox frame grabbers) require USB-Vid I/F adapter
 - IEEE1394 or "Firewire" drivers are required for Q Imaging CCD cameras.
 - All other cameras require Andor Virtual Camera (AVC) which captures images from the native application Live camera Window.
- 4. Komet or Komet GLP software
- 5. Database Components.

A separate Database Viewer Installation CD is supplied and instructions on the installation are provided in the Database Viewer User Guide.

For a detailed description of the installation procedure, please refer to **Appendix 1**. If this does not help, our technical support team will be glad to assist. To report a problem please contact us at <u>www.andor.com/contact_us/support_request</u>, or contact your supplier who will then liaise with us directly.

(Appendix A1.2 and Section 5)

(Appendix A1.2)



1.3 New Features in Komet

Komet 7 is certified Windows 7 compatible and provides improvements to dataset management as well as a number of bug fixes and increased robustness measures in the overall software platform. To ensure backward compatibility we have retained the analysis algorithms and options as presented in version 6. Komet 7GLP has been extended to ensure compliance with data handling and reporting requirements of OECD TG 489, mainly around the treatment of GLP Datasets and in the Database Viewer software.

A new range of cameras and camera interfaces is supported offering both new opportunities to speed scoring and improved ease of use of microscope by offering LED (light emitting diode) illumination options. As always we provide you with flexibility to keep working your existing hardware or to upgrade your system with the advance of Windows and imaging technology.

1.3.1 New Imaging Interfaces and Digital Camera support

To keep pace with the rapid developments in technology we have improved the Komet software to support the wide variety of analogue and digital cameras available today. Andor is known for excellence in scientific cameras and our most recent technology is showcased in Zyla 5.5, a camera whose field of view is almost as large as the microscope eyepieces, making on-screen scoring of comets even faster and more comfortable. With its highly convenient USB connection, Zyla offers exceptional quality and sensitivity not only for comet assay, but for a multitude of other applications.

At the same time we recognize that budgets are under pressure in many laboratories and that backward compatibility with existing hardware is critical. Therefore with Komet 7 and GLP we introduce the Andor Virtual Camera (AVC) which supports older video CCD cameras via cost-effective Video-USB adapters (VID2-USB2 from Star and DFG/USB2Pro). If you own a Matrox Metor or Metor II card, this card is now obsolete and must be replaced with the Star or DFG products.

Moreover, AVC supports Komet analysis from ANY camera running under Windows 7 and able to deliver an image into a "live window"; this is achieved by capturing an image directly from the screen and analysing it. We recommend that you setup a second display monitor for this live image window so that Komet does not have to share screen space during scoring.

1.3.2 Database Viewer

Database Viewer (DBV) was first introduced with Komet 5 and rapidly became a most popular feature due to its ease of use and power. Uniquely DBV saves an image of every comet scored which



is linked to the associated analytical data. Data obtained from a Komet experiment is now saved by default into the Database format which we refer to as a Dataset. In GLP versions, Komet even captures comets that were scored but rejected along with the reason for this action.

In version 7, DBV has been made even more robust and performance improved. A number of formatting bugs have been addressed e.g. Date and time have been changed to a universal international format (e.g. 20160206: YYYMMDD meaning 6th Feb 2016). Audit Trail date formatting has been made consistent with this and its redundancy reduced.

Efficiency has been markedly improved in reading large data sets. A speed up of approximately ten times has been achieved with typical in-vivo comet studies involving 36 exposure units and approximately 5400 data points.

Additional functionality has been provided to ensure full compliance with OECD Test Guideline 489: In-Vivo Mammalian Alkaline Comet Assay. This includes the option to "log transform" raw data before summaries are produced and optional reporting of the median data for each individual slides.

As before DBV provides fast and easy to use tools for critical functions such as QA data review, specimen decoding (essential for GLP studies) and creation of data summary tables. These functions speed data review and preparation for statistical hypothesis testing and are key to the DBV's popularity. Another benefit is that <u>DBV is a free distribution product</u> so that you can share Datasets with colleagues and sponsors while maintaining security and data integrity via password and electronic signature protection.

For researchers who are not working under GLP or submitting their data for Regulatory Approval, DBV provides a powerful tool for quality assurance of data in the research context and offers unique tools to explore and contrast data in a teaching and research environment. DBV allows students to explore Datasets, understand the concepts of decoding and statistical presentations in bar charts, 3D histograms and dose response curves. Where analysis not included with DBV is required, then DBV offers an export to Microsoft Access file format. Free distribution means that the benefits can be shared throughout the lab or department.

1.3.3 Printing Datasets in Database Viewer

In some laboratories, recoding of electronic records requires printable formats to be stored. DBV now supports efficient printing and storing of any Microsoft-compatible format including PDF and XPS. This mode is available for Image Galleries, Data Tables and Audit Trails.



1.3.4 LED Light Source Recommendation

Over many years of experience we have found that one of the main technical issues for nonspecialist microscope users is the arc lamp. Because the arc lamp gets very hot, emits intense light with UV content and usually contains mercury vapour it is hazardous and must be handled with great care. Moreover standard lamps must be replaced every few (2-3) hundred hours, disposal must be managed and their intensity and stability drops rapidly as they approach the end of life. Alignment of the lamp is also critical for good quality (uniform) illumination, otherwise image quality and data analysis can suffer. Fortunately, Light Emitting Diodes (LEDs) are now universally available and have been adapted to microscopes.

Andor recommends the use of LED light sources due to their longevity, safety, cost of ownership, uniformity of illumination and the fact that no specialist knowledge is required to obtain optimal performance. Typical lifetime of an LED is 20,000-25,000 hours, meaning that with an average working week of 40 hours, you should only need to replace them every 10 years if you use your microscope eight hours a day and 5 days a week! Moreover, LEDs are highly efficient, have virtually zero warm up and cool-down times, produce little heat and are "optically quiet". This means they do not produce high frequency fluctuations inherent to arc lamps, leading to lower variance in your image analysis. We now offer a range of LED light sources suited to your microscope, budget and experimental needs. Please let us know if we can advise you on such a device.

1.4 Other Andor Products

We hope you are aware of the range of Andor products for quantitative fluorescence imaging and microstructure quantification. Having entered the field with Komet, these other products might also help in your application of scientific imaging. Further information can be seen at www.andor.com. Full details will be sent on request. Andor Imaging was acquired by Andor Technology, an Oxford Instruments company in 2004 and is now Andor Bio-imaging Division. Komet remains under the Andor Imaging name for historical purposes.

 Andor iQ - Our flagship Live Cell Imaging application. Andor iQ provides powerful multidimensional acquisition and analysis for applications such as intra-cellular ion imaging, timelapse fluorescence for GFP and developmental studies. Andor iQ offers extensive support for CCD and sCMOS cameras, filter wheels, stages and motorized microscopes. Andor iQ provides powerful tools for specialized analysis of dynamic image data. It offers users flexibility in image acquisition, processing, analysis and storage.

Andor Bio-imaging Division has extensive development facilities and expertise in both fluorescence imaging hardware and software. Custom solutions can be developed.



SECTION 2: KOMET AND THE COMET ASSAY

2.1 An Imaging Protocol for 'KOMET' Analysis

When using the Komet Analysis system it is important to remember that this is an instrumental tool and requires an appropriate and reliable protocol for its use. This is true of any analytical software or instrument. Komet analyses the distribution of photometric data gathered through the image sensor of a camera. If reliable quantitative results are to be obtained, then you should use the suggested protocol in the following sections, or develop a similar one to suit your own requirements. If care is taken, the system will reward the researcher with an abundance of useful information. For a detailed explanation of the Komet analysis algorithms see **Appendix 3**.

Note: Cells MUST be imaged with the tail extending to the right of the screen.

2.2 Background Effects

In working with a number of research groups we have found that the uniformity of the gel preparation is an important factor in analysis, because an uneven gel may lead to varying or structured background intensity in the Region of Interest (ROI). Since the background is integrated and used to correct all intensities in the adjacent comet cell, it is important that there is a uniform and slowly varying background function. If your specimens show such varying or structured background intensity, the gel should be thickened a little to overcome some of the image structure produced by the ground glass slide. Better still, set up the dried slide protocol, which uses plain slides and exhibits low background intensity as described in **Section 9**.

The background effect will be even more apparent if Ethidium bromide (or Propidium iodide) solution shows signs of crystallization. Crystals will be apparent as very bright spots in the field of view. The starting solution should be filtered regularly to avoid this, or new aliquots made up frequently.



2.3 Comet Visibility, Electrophoresis Time and Tail Length

Clearly the electrophoresis time is a critical factor in affecting the resultant comet image. To enable the best performance of the Komet system it is important to aim for comet cells that can be seen fully in a single image frame. This is normal when viewing with a 16X or 20X objective, but in some cases where extensive DNA damage has occurred and electrophoresis time is long, the tail may be extremely long, causing analysis problems.

Also bear in mind that a comet tail is only of value if a reliable analysis of its characteristics can be achieved. Thus it will be clear that the longer a tail, assuming a fixed quantity of DNA, the less intensity it will show in a given elemental area and the more difficult it will be to image. A very bright head and a very dim tail require a wide dynamic range of the detection system. Normally a video system will deliver 8-bit dynamic range, while a cooled digital camera will provide 12-16-bits. Technically, Komet can support any of these cameras, but cost considerations may preclude further investment. It will be clear that a little effort expended in matching the comet cell preparation to limit tail length and brightness range will be well spent.

2.4 Analogue CCD Imaging

It should be stressed at this point that the Comet Cell preparation is not a major problem, but one of matching the preparation protocol to the analyser. It is more likely to be troublesome with older microscopes and in cases where 100% of the image cannot be directed to the CCD Camera. If you have sensitivity problem please contact your local supplier for details about the options we recommend

The **Brightness** and **Contrast or Exposure** controls in the Camera Settings dialogue should also be setup and fixed to ensure the most suitable imaging performance. The best approach to use for these settings is to present typically bright and typically dim comet cells to the system and adjust the imaging controls, as far as possible, to obtain a good quality image,. This normally involves setting the Brightness to offer a dim but visible background and adjusting the Contrast to ensure little saturation of the system. Refer also to Mapping Controls in the Komet Function Reference, Section 4.



2.5 Digital CCD and sCMOS Cameras

For users with 12 or 16-bit digital cameras: We normally recommend utilizing the binning feature of these camera only if sensitivity is an issue. Select a binning value of 2 and adjust the exposure time to provide a bright image of approximately 2000 grey levels (for 12 bit cameras) and 15000 (for 16 bit), taking care to avoid saturation.

If Gain and Offset are available, we generally recommend setting both to a mid-range value. Many digital camera control interfaces have an Auto-Exposure feature, which can be used to obtain a suitable exposure time.

sCMOS Cameras (such as the Andor Zyla series): largest field of view almost matches the microscope. This enables more multi-comet scoring in each acquired image, speeding up the scoring process and delivering data faster than ever.

.



2.6 Hints on Microscopy

Comet slides are normally viewed on a fluorescence microscope, because most convenient labels for DNA are fluorescent e.g. Propidium iodide, Ethidium bromide, DAPI, SYBR green. In some studies, DNA stains have been used which are visible in the microscope in brightfield mode e.g. silver stain, Feulgen. It is perhaps obvious, but often overlooked, that correct setup of the microscope is of central importance in obtaining good quality results from the comet assay. The following comments are intended to help you get the best from your instrument.

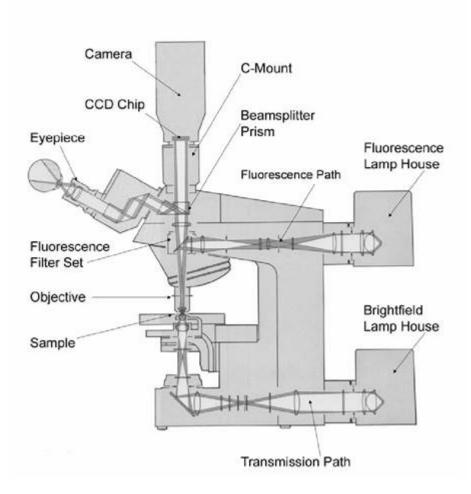


Figure 1. Microscope Schematic (after Zeiss) Showing Fluorescence and Brightfield Paths



2.7 Light Source, Lamp-house and Epi-fluorescence path

A schematic of a microscope with both epi-fluorescence and transmission illumination is shown in Figure 2.1. We will focus on the epi-fluorescence light path. The light source will be for example a Mercury arc or Xenon arc lamp. These lamps rely on maintaining an electrical discharge (arc) between metal electrodes in a mercury or Xenon vapour-enriched atmosphere, which is encapsulated in a Silica envelope. The lamps get very hot in operation and must be handled with great care. Do not re-ignite a lamp within 15-20 minutes of extinguishing and do not handle for the same period. Ozone-free lamps are generally preferred and less hazardous, because they do not emit very short wave UV. Even these lamps, however, must never be viewed directly or there is a serious risk of damage to your eyes.

Note: Andor recommends the use of LED light sources due to their longevity, safety, cost of ownership, uniformity of illumination and the fact that no specialist knowledge is required to obtain optimal performance. Typical lifetime of an LED is 20,000-25,000 hours, meaning that with an average working week of 40 hours, you should only need to replace them every 10 years if you use your microscope eight hours a day and 5 days a week! Moreover, LEDs are highly efficient, have virtually zero warm up and cool-down times, produce little heat and are "optically quiet". This means they do not produce high frequency fluctuations inherent to arc lamps, leading to lower variance in your image analysis.

However, if you do own a conventional arc lamp, here are some tips on how to set align them for optimal performance.

Arc lamps should be replaced after approximately 300 hours of use. Lamps will occasionally last for longer or shorter periods and users should be alert to the symptoms of ageing. These include reduced intensity, as metal from electrode erosion is gradually deposited on the silica envelope, and instability in the arc, which is seen as intensity fluctuations and can lead to apparent random variations in image intensity.

When replacing an arc lamp, be sure to use gloves to avoid leaving deposits of oil from your hands on the Silica envelope. Such deposits can lead to hotspots and lamp failure by cracking of the Silica or, in severe circumstances, explosion! Always wear UV goggles to protect your eyes and always follow the instructions in your microscope guide to ensure correct replacement procedure. If in doubt, seek help because this is a potentially hazardous operation.

Once the lamp has been replaced, it is important to follow the correct procedure for alignment. A badly aligned lamp will result in reduced and uneven illumination intensity, resulting in a poor fluorescence image as seen by your eye and the CCD camera. The arc lamp-house has a reflector (mirror) in it. When the illuminator is correctly aligned the image of the arc itself and the reflected



image will be adjacent to one another in the illuminated field (Figure 2) and the image pair will be centred in the field. A centred pair will result in uniform and bright images, provided the collector lens focus is correctly set. Once again, you should follow your microscope setup procedure.

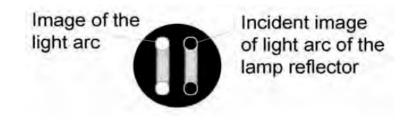


Figure 2. Focusing a fluorescence lamp

After the arc lamp collector lens has been correctly focused, it is important to make sure that the rest of the fluorescence illuminator train is centred and correctly adjusted. The train contains an aperture or "field stop". Make sure this is open sufficiently to just illuminate the whole required field-of-view and that it is correctly centred.

2.7.1 Fluorescence Filter Set and Microscope Objective

The light from the source is projected onto the specimen by way of the fluorescence (interference) filter set, specifically the excitation filter (Figure 3). The fluorescence filter set performs a crucial job in ensuring an image with high intensity and low background. This is important for effective imaging and analysis of the comet cell data. The filter set must be matched to the fluorescent label used. This means making sure of the following points:

1) **Excitation filter** – the peak transmission wavelength should closely match the peak excitation wavelength of the fluorescence label.

2) **Dichroic filter** – the transmission cut-on edge wavelength should be close to the minimum overlap between the excitation and emission spectra of the fluorescence label.

3) **Emission filter** – the transmission wavelength should be close to the peak emission of the fluorescence label.

If using a microscope inherited from previous research, or even a new instrument, make sure that the correct filter set is available. If this is not the case, investment in a tailored set is worthwhile. Several companies make these, including Chroma. Andor Bio-imaging Division is an agent for Chroma and can provide all such filters.



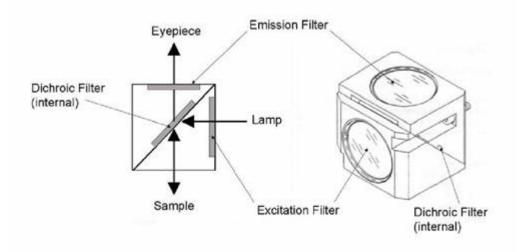


Figure 3. A typical fluorescence filter cube schematic

Please note that interference filters are made by evaporation of very thin layers of dielectric materials onto glass or Silica substrates and consequently should not be handled without gloves. Deposits of oil or other surface contamination will adversely affect the performance of the filters.

Some older microscopes were designed primarily for observation by eye and may use a long-pass emission filter, which transmits infrared (IR) light. Some CCD cameras exhibit substantial sensitivity in the IR region which may result in a very low contrast image. The solution is to obtain a band pass emission filter, or, where suitable, fit an IR cut filter to the CCD camera. The latter will reduce sensitivity, so the former is preferred when performance is limited by light throughput.

Furthermore, since both the excitation and emission light pass through the microscope objective, it is important to make sure that the lens is capable of a high level of transmission at both wavelengths. Fluar and Plan-Neofluar lenses are best suited to UV excited labels such as DAPI, but others are satisfactory for longer wavelength operation, e.g. Propidium iodide (green excitation and red emission). Obvious, though often overlooked, is the fact that lenses must be kept scrupulously clean. Use only proper lens tissues and alcohol for cleaning. Dry lenses must be kept free of immersion oil and oil lenses must be cleaned with alcohol after use. Dirty lenses can result in image distortion, low light levels and generally poor imaging performance.



2.7.2 Beam-splitter and C-mount Adapter

Fluorescence light from the specimen is actually emitted in all directions, but in the epi-fluorescence microscope, the light returning to the filter set along the excitation path is collected for image formation. It makes its way through the dichroic filter, by virtue of the fact that the filter transmits the emission wavelength. The emitted light is further selected by the emission filter, to produce a low background image of the specimen. The beam-splitter/prism directs the image of the specimen to the eyepiece or to the CCD camera, and in some cases, to both. The best condition for imaging results is when 100% of the light from the specimen is directed to the CCD camera. In older instruments this is especially important, because light levels are inherently lower. In some instruments, stray light from the eyepieces and elsewhere in the instrument can find its way through the prism into the camera, creating a poor contrast image. For this reason, fluorescence microscopy is often carried out in a darkened room and the eyepieces may be covered. This makes viewing the image on the computer screen even more desirable.

The fluorescence image, which is finally projected onto the CCD camera, must arrive in the correct optical plane for a sharp image. This plane should be "parfocal" with the eyepiece image to avoid refocusing between the eyepiece and camera. To achieve this and mechanical compatibility with most CCD cameras, a C-mount adapter is used. The C-mount is the almost universal CCTV connection and has been adopted by the majority of digital CCD camera manufacturers. The standard C-mount has a one inch (25mm) aperture and a 32 TPI thread.

In many cases the camera will have a CCD chip which is much less than one inch (commonly twothirds, half or third inch CCD chip) and so the field of view it sees will be much less than that projected into the eyepieces and the camera image plane. Hence, some C-mounts have demagnifying optics built in to increase the camera field of view. Common C-mount magnifications are 0.63x, to match two-thirds inch CCDs, and 0.5x, to match half inch CCDs. A 1x C-mount will not adversely affect imaging however, and provided the user understands the issues it is not essential to match the adapter to the CCD chip. Note that scientific CMOS such as our Zyla 5.5 camera also uses a 1X C-mount adapter. The sensor diagonal is 22 mm and images almost the entire field of view of the microscope.

2.7.3 Summary

The points above should be borne in mind when setting up the imaging system for the comet assay and indeed any microscopic imaging of fluorescent samples. If problems are encountered with imaging specimens, the material can be used as a checklist for troubleshooting and problem solving.



2.8 User Management and Logon Procedure

Andor software products now have a 2-level logon procedure. The first level is the Windows logon and the second is the Komet logon when the software starts. The Komet Research logon involves selection of a Configuration that is essentially a hardware configuration, along with associated Protocols and Channels. All Users (i.e. Windows users) share these hardware Configurations on a system. Komet GLP (see Sections 5, 6 and 7) has a more sophisticated user management system because FDA 21CFR Part 11 requires this. Please refer to that section for further information.

Configurations are created by the LOCAL system Administrator using the Configuration Manager utility (Section 6). Within the Configuration, each Windows User has his/her own data paths, preferences, etc. After the Configurations are set up by the system Administrator, it is recommended the software be run by the Administrator to ensure that hardware settings are correctly configured (e.g. COM ports, board settings, calibrations). The Configuration Manager also allows the Administrator to control Global Settings, which relate to ImageDisk organization and whether Multilevel User Access is enforced by the Komet software. When enabled, Multi-level User Access allows different permissions to users, as determined by the system Administrator. If Multi-level User Access is OFF, then all users will have full access to the software.

Members of Windows Local Machine Administrators and Power Users groups will have full access to the software, but Restricted Users will only be able to use Channels and Protocols set up by other users.

It is important therefore that Komet software users log on to Windows as themselves. It is also important that Users maintain their access integrity if access control is important. For example, this will be true where the system is used for GLP studies.



2.8.1 Running Komet While Logged Onto a Domain

Some customers may wish to run Komet on a computer while logged onto a remote domain. In this case, Windows grants limited access to certain areas of the local machine. Therefore, it is very important that the LOCAL Administrator runs Komet first and sets up the hardware device settings.

By default, from Win2000/XP onwards, all Domain Users are treated as restricted users on the local machine. In order to allow Domain Users full access to the Software, e.g. to be able to create, edit and delete Protocols and Channels, if required, those Users should be made local Power Users (Windows On-line Help).

Note to Administrators: To find out more about Domain User access, go to the Windows Start Menu\Settings\Control Panel\Users and Passwords and use Windows on-line help. (For Users with Windows NT 4, go to Start Menu\Programs\Administrative Tools\User Manager).

The Access Levels provided in Komet depend on Local Status and are summarized in the following table:

	Full+Configuration	Full	Restricted
Local Admin	Y		
Local Power User	Ν	Y	
Local User	Ν	N	Y
Domain Admin	Ν	N	Y
Domain Power User	Ν	N	Y
Domain User	Ν	N	Y
Domain+Local Admin	Ν	Y	
Domain+Local Power	Ν	Y	
Domain+Local User	Ν	N	Y

Table1. Access Levels in Komet with Multi-Level Access ON



2.8.2 Configuration Manager

Refer to Appendix 2.

2.9 Repeatability Measurements

In order to evaluate and prove the repeatability of the Komet analysis system, we have undertaken extensive testing and evaluation. Each new release undergoes the same rigorous testing regime before we supply it to our clients. The results summarized in Appendix 4 are included to establish the reliability of the system and its measurements. Both static tests, for ensuring algorithm correctness, and live tests, to establish noise performance, have been carried out. The results show that even under noisy conditions, with a little attention to preparation and imaging conditions a repeatability of a few percent can be achieved for even the most sensitive parameters, which are generally those associated with the tail. We hope that you find these of interest, since they establish the Komet system as a reliable tool for analysis. They also act as examples of how you can undertake a similar study in your own lab, with your own samples, to identify the significance of various parameters to your scientific endeavours!



2.10 The Benefits of Komet

2.10.1 Flexibility, Protocol Setup and Data Integrity

A major benefit of the current version of Komet is Protocol management. Using the Setup Wizard (**Section 3.3**), the users are freed of the detail of settings, once a Protocol has been created. Every Komet Protocol contains Comet Options, Calibration and Camera Settings selected for a given imaging and data analysis protocol. These can be set up once for a given experimental procedure and re-used by simply selecting from the Comet Protocol Manager. This reduces settings errors and makes changing settings mid-experiment more difficult, therefore, adding enhanced control of data obtained during scoring.

Protocols also facilitate recovery of an experiment following interruption, thus ensuring that the system is driven by the user's requirements. To interrupt an experiment simply press the **Abort** button on the Comet Toolbar, then when asked if you wish to suspend the experiment, select **Yes** and close the system down (**Section 3.3.4.2**). On start-up, Komet will recover the current user settings. When you press the **Experiment** button you will be asked whether you wish to resume an experiment (**Section 3.7.6**). The data will be stored in the same automatic data structures. Ensure **Use Automatic Data Directories** is selected in the Experimental Comet Options (**Section 4.12**).

2.10.2 Calibration - Simple and Infrequent

Once the system has been calibrated through the Calibration Wizard with a given microscope and objective, the calibration will be saved to file. Every time the system is started, Komet will recall the calibrations without further user-intervention. If the objective is changed, simply create another calibration from the Calibration Wizard as described in **Section 3.4**. These calibrations are then available when creating or editing Komet Protocols.



2.10.3 Overlays – Region of Interests (ROIs) and Calipers

A live video overlay is used to show the cell Region of Interest (ROI) while Live. Thus, the ROI is always visible during specimen sampling and the user need only locate a new discrete cell in the ROI before clicking on **Store**, or clicking the right mouse button on the image window to complete each analysis. However, if other cells in the field are to be scored, the ROI can be picked up and positioned with the mouse. Analysis proceeds by a right mouse button click.

When Analysis delivers its estimates, e.g. after an **AutoExpress On Store, Right mouse click** or **Interactive analysis (Section 4.2.6)**, Komet presents head and tail extents, indicated by red lines (vertical calipers) in the image screen overlay. Figure 4 is a comet cell image showing the Cell ROI, automatic background ROI and analysis calipers, delineating the start of Head, end of Head and end of Tail.

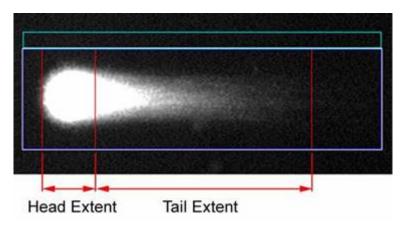


Figure 4. FA Comet Cell ROI after analysis



2.10.4 Dose Statistics in a Secure Access Database

In the latest version of Komet the data can be stored in a secure database, by selecting **Use Database** in the Comet Options Experimental section of your Protocol (**Section 4.12**).

Komet is distributed with a Database Viewer application, where the Comet data can be viewed and analysed. Histograms, Bar Charts and Dose Response curves can be produced and printed out from the Database Viewer. An Image Gallery displays all the comet cells scored in the selected dose, together with the calculated data for each cell.

2.10.5 Suspending and Resuming Experiments

It is now possible to interrupt an experiment before its completion. The current state of the experiment is remembered and at a later date you can continue from the same point. To **Suspend** an experiment press the **Abort** button on the Comet Toolbar and you will be asked whether you wish to **End** or **Suspend** the experiment.

If you choose to suspend the experiment, you will be asked to enter a name for the suspended experiment and notes describing the experiment. If you have many suspended experiments, entering notes can be useful for quick reference (Section 3.7.5).

To **Resume** an experiment that has previously been stored, press the **Experiment** button on the Comet Toolbar. If there are no suspended experiments, a new experiment will start as normal. If there are suspended experiments, you will be asked whether you wish to start a new experiment or to select a suspended experiment to resume (**Section 3.7.6**).



2.10.6 Conclusion

We hope you will benefit from this new release of Komet and we look forward to continuing to support the scientific community in the future. Please let us know of any further suggestions and enhancements that you may wish to see.

Back issues of the Comet Newsletter edited by Dr Luc Verschaeve, VITO, Belgium, which has been produced for all users of the SCGE assay, are available on the Andor Imaging website (See bottom of page at the following link) <u>http://www.Andorimaging.com/Komet.html</u>. The publication of the newsletter is currently suspended due to the large number of other publications relating to the method.



SECTION 3: GETTING STARTED

3.1 Starting Komet

When Komet is started, a User Registration dialogue is presented (Figure 5). The User Name, which you enter, will be saved in the Registry. When you exit from Komet the current system setup is saved in the Registry for that User Name.

Registration	X
Enter User Name:	
Test	•
ΟΚ	Cancel

Figure 5. User Registration Dialogue Box

A System Setup includes the parameters for Window Size and Position, File Save paths and Comet Protocols. Data in the Analysis Table will also be automatically saved to the file **LastSession.Xld** on exiting Komet. This file will be saved in two directories: 1) The directory from which you are running Komet and 2) The directory:

C:\Andor\Komet CAMERA**Data\UserID**, where **UserID** is the User Name entered on starting Komet. If this directory did not previously exist, it will automatically be created.

Note: This file directory can be changed by selecting **Set Save Path** from the Komet Control Window File menu.

The next time Komet is started, if you enter a User Name that already exists, the previous system setup for that user will be loaded from the Registry, and data loaded from the file **LastSession.Xld**. Komet will therefore be setup the same as when that user last logged off. If a new User Name is entered, the last used system setup will be loaded and a new entry created for the new user.



3.2 The Komet User Interface

KOMET offers a powerful user interface for comet cell analysis within the Microsoft Windows 98, NT, 2000, XP and onwards environments. This means that not only is the product attractive and easy to use, but you can also use a vast range of other compatible software products within the Windows environment.

The Interface consists of two floating dialogues, which can be positioned and resized as required (Figure 6).

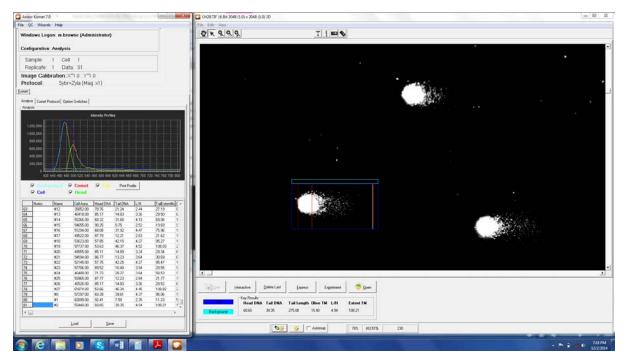


Figure 6. The Komet User Interface Image Window and Komet Control Window

.



3.2.1 The Image Window

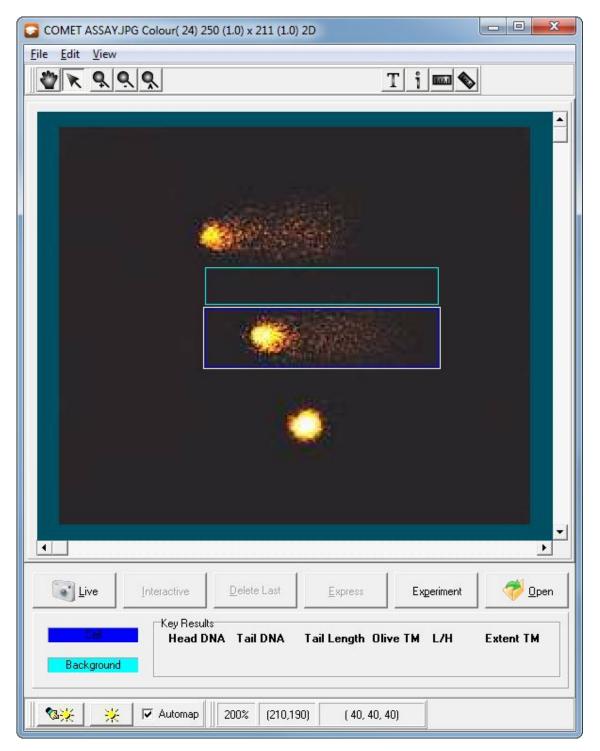


Figure 7. The Komet Image Window

The Image Window is where you view Comet cells (Figure 7). Open the images into the Image Window or go Live, position the Cell Region of Interest (ROI) around a Comet cell and then perform a comet analysis.

-0



3.2.2 The Comet Toolbar

The Comet Toolbar (Figure 8) supplies a rapid and convenient means of controlling the system both during an Experiment (**Section 3.6 – 3.7**) and in free mode.



Figure 8. The Comet Toolbar

3.2.2.1 Live/Store



Clicking on **Live** switches the system to live image grabbing mode and changes the button label to **Store**. When you have a specimen suitably positioned within the cell ROI, click on the **Store** button or more conveniently click the Right mouse button (**Section 4.2.6**).

3.2.2.2 Interactive



Cells may exhibit fragmented comet tails or unusual morphology after certain treatment. This can lead to inaccurate estimation of the Comet parameters. The **Interactive** function enables the user to define the Tail ROI under such circumstances (**Section 4.2.6**).

3.2.2.3 Delete Last



Successful analysis of a comet will insert a row of data into the analysis table in the Komet Window (**Section 4.2.6**). Pressing the Delete Last button will remove the last row of data from this table.

3.2.2.4 Express



Press the Express button or right click on the image window to analyse a Comet Cell. This will compute all the parameters and post them to the table on the Analyse tab (Section 4.2.6). Note that clicking the Right Mouse Button while the cursor is on the Image Window will achieve the same action.



3.2.2.5 Experiment/Abort



The **Experiment** button will switch Komet into Experiment mode (**Sections 3.6** – **3.7**), in which it keeps track of how many cells have been scored and prompts you when to change slides, save data and begin a new sample.

<u>S</u>uspend

When you click on the **Experiment** button, the button label changes to **Suspend**. Click on the **Suspend** button if you wish to cancel the Experiment before its completion, or Suspend it to Resume at a future date (**Section 3.7.5-3.7.6**).

3.2.2.6 Open



The Open button is used to load a stored image file. This has the same function as **Open Image** on the **File Menu. (Section 4.2.1.1)**.

3.2.3 The Key Results Panel

C AN	Head DNA	Tail DNA	Tail Length	Olive TM	L/H	Extent TM
Background	66.13	33.87	83.02	9.27	5.13	28.12

Figure 9. The Key Results Panel on the Image Window

Key results from the latest Experiment / Express Analysis are displayed in the Key Results dialogue for quick reference (Figure 9), as well as in the results table on the Analysis dialogue (Figure 10).

The Dark Blue ROI in the Image Window defines the Cell region.

The Light Blue ROI in the Image Window defines the Background region.



Komet[®]

3.2.4 The Komet Control Window

🖸 Andor Komet 7.0 -	Andor Virtu	ial Camera			X
<u>F</u> ile <u>Q</u> C <u>W</u> izards	<u>H</u> elp				
Windows Logo	n: j.webb				
Configuration:	vcam				
Sample:	1 Cell:	1			
Replicate:	1 Data	: 4			
Image Calibration:X*1.0 : Y*1.0					
Protocol:			l (Mag: xi	1)	
Komet					
Analyse Comet Pro	tocol Option	Switches			
Analysis	· ·				
		Intensity Pr	ofiles		
7,000					
6,000		\wedge			
5,000		/÷i-∖-`.			
4,000					
3,000	<mark>.</mark>		4		
2,000	<mark>/</mark>	<mark>A</mark>			
1,000					
0			1		
	430 440 450	460 470 480	490 500 510 9	520 530 540	550 560 570
R Backgr	ound 🔽	Comet		Print Pro	file
I Cell		Head			
Notes	Name	Cell Area	Head DNA		L/H
1	#0	5616.00	50.92	49.08	3.36
2	#1 #2	5130.00 5728.00	56.40 69.72	43.60 30.28	3.32 2.27
4	#3	5549.00	52.36	47.64	2.37
Image: A state of the state					•
	L	oad	Save		

Figure 10. The Komet Control Window

The Komet Control Window shows the Analyse, Comet Protocol and Option Switches tabs (Figure 10). Analysis data is displayed on the Analyse tab in both graphical and numerical format (Section 4.10). The comet options, system calibrations and camera settings are set up from the Comet Protocol tab prior to analyses (Section 3.5.2). The Option Switches tab allows certain parameters to be changed during an experiment without invalidating your data (Section 4.8.2).



3.3 Working with Live and Stored Images

In Komet you can score comet cells viewed live through a microscope, or from images saved to disk.

3.3.1 Live Images

When working with Live images, you should calibrate your microscope objective lenses as described in **Section 3.3**. In the first step of the Calibration Wizard the **Live Image** option is selected for the **Type of Calibration Required**.

When setting up a Comet Protocol (**Section 3.5**) the first step in the Wizard is to select a calibration. If you have calibrated all the objectives, select the calibration for the objective in use. The Comet Setup Wizard will then move on to the Camera settings. With the correct objective in place select **Live** to image a comet sample. The Camera settings can then be adjusted to obtain a good quality image (**Section 2**).

The final step of the Wizard is to specify the Comet Options for your experiment (Section 3.5.5).

Note: If **Floating Live Button** is selected in the Comet Options a **Live** button is displayed next to the mouse cursor after each score, which greatly reduces the amount of mouse movements between acquisitions.

You are now ready to run a Live experiment (Section 3.6).



3.3.2 Stored Images

When working with stored images you have the option of either using the calibration stored in the image headers, or applying a calibration previously defined in Komet. Some formats, such as Bitmap (BMP), do not store calibration information, so Komet must apply a calibration to them.

With **Apply selected calibration on Open Image** selected in the first step in the Comet Protocol Setup Wizard (Figure 12), the calibration selected will be applied to all opened images, overriding any calibration that might be present in the loaded image.

Note: This is only valid if the calibration applied has been created from an image of a calibration device produced under the same experimental conditions as the images being processed.

If **Apply selected calibration on Open Image** is NOT selected, the calibration will be taken from the loaded image, therefore great care must be taken to ensure all images have been acquired with a valid calibration (software dependant). The Wizard will then show the Camera Settings.

Note: This will not be present if the version of Komet does not support image acquisition. This part of the Wizard can be ignored when setting up a stored image configuration.

In the Comet Options Operational section, the **Floating Open Button** feature can be used (**Section 4.2.6**). This displays an **Open** button near the mouse cursor after scoring each comet. In Experiment mode you can specify a source image folder for each dose. The next image is automatically loaded on pressing **Open**.



3.3.3 Notes on Single versus Multi-Cell Scoring

The most repeatable performance will be achieved by using a fixed cell region of interest (ROI) box size and background regions, but this is not essential as shown in the repeatability studies in **Appendix 4**. Maintaining the position of the cell ROI will generally improve the overall precision of the measurements, especially with arc lamp sources where illumination intensity varies throughout the field, and this directly affects the resultant fluorescence intensity in the cell. LED and other light-guide coupled sources generally behave much better in this respect. When scoring multiple cells per field make sure that the fluorescence illumination has been set up correctly (**Section 2.6**) and show good uniformity across the field of view.

3.3.4 Free and Experiment Modes

The system has two modes of operation, Free and Experiment.

3.3.4.1 Free Mode

In Free mode, data is gathered and can be handled for storage, analysis or evaluation. Start by loading an image or going Live using the microscope camera, use the system default settings and then size the comet cell region of interest (ROI) large enough to enclose a full comet area; head plus tail (Figure 11). The ROI can be moved around the screen easily, using the mouse, to score several cells per field.

The system expects the comet tail to extend to the right of the image, so make sure you present the cells in this direction and make sure you select a suitable objective if you are live imaging.

We recommend you leave **Auto Background** on, if possible, but if you have switched it off in the Comet Options or Option Switches dialogue you will need to set up the background region. The system will warn you if you do not do this. If you have selected **Auto Background**, the system will place the background ROI immediately above or below the cell ROI. As with all background selections, you should avoid analysing cells where the background ROI contains other cells or substantial debris, as this will generate an underestimation in the measured cell intensity parameters.



Note: When going **Live** from the Comet Toolbar, the cell ROI is overlaid on the live image. This means that once set, you can simply navigate the microscope slide until the next cell for analysis is within the ROI and then click on **Store**.

Alternatively for scoring from file, open an image from your chosen directory by clicking on the **Open** button. Position and size the cell ROI. The Background ROI should follow this and be positioned immediately above or below the cell. To complete an analysis, click the **Express** button on the Comet Toolbar. Express integrates the cell and background in the regions specified, subtracts the background, smoothes the data and then proceeds with analysis. Alternatively, a Right Mouse click will perform the same Express action. The analysis results will be displayed in the Komet Window.



3.3.4.2 Experiment Mode

In Experiment mode (Section 3.6-3.7) the system will prompt the user:

- Each time a slide is complete: The correct number of cells has been measured, defined by the **Cells per Slide** parameter in the Comet Options.
- At the end of a sample: When the correct number of cells has been measured for all **Replicate Slides** in the sample.
- When data is saved.
- At the end of the Experiment: When all the **Samples** have been measured.

Cells Per Slide, Replicate Slides and Samples are specified in the Experimental section of the Comet Options (Section 4.12).

When you are ready to start work in earnest and embark upon an experiment, click the **Experiment** button on the Comet Toolbar (**Section 4.2.6**). The Results buffer will be cleared and you can begin your measurements, as described in **Section 3.3.4** (this assumes that **AutoExpress on Store** is selected in the Comet Options). The system will prompt you at the message points in the following procedure.

See **Section 4.2.6** on how to use the Comet Toolbar during an experiment.

Note: At any point in an Experiment the **Abort** button on the Comet Toolbar can be pressed. The user is then given the option of ending or suspending the experiment (**Section 3.7.4**).



3.3.5 Live Image Experiment Sequence

The following is a pseudo-code presentation of the sequence of Cell, Slide (Replicate) and Sample (Specimen or Donor) cycle defined within a Live Image KOMET experiment.

Start Experiment

Message "This action will delete your current data. Save it first?"

Message "Do you want to Resume an experiment?" (If any are present) If Use Automatic Data Directories is selected, enter Experiment ID

For M Samples

If Use Automatic Data Directories is selected, enter Sample ID For N Slides

For P Cells

- Select 'Live'
- For scoring single Cells:

Position the comet cell within the Cell ROI with the head towards the left and preferably, close to centre of the image screen

• For Scoring Multiple Cells:

Select the Comet Cell to score, then drag the ROI over the cell to be scored

• Select 'Store' or click the Right mouse button on the image window

Next Cell

Message "Select the next Replicate Slide, OK to continue, Cancel to repeat last measurement"

Next Slide

Message "Sample complete, data will now be saved, OK to continue,

Cancel to repeat last measurement"

If Use Automatic Data Directories is NOT selected, enter a directory and filename for the data to be saved

Next Sample

Message "Experiment complete" End Experiment.



3.3.6 Live Image Experiment Sequence With Database

Start Experiment

Message "This action will delete your current data. Save it first?"

Message "Do you want to resume an experiment?" (If any are present) Enter Experiment ID (entering other setup parameters is optional)

Message informing the user of the path to the database For M Samples

If Use Automatic Data Directories is selected, enter Sample ID

For N Slides

For P Cells

- Select 'Live'
- For scoring single Cells:

Position the comet cell within the Cell ROI with the head towards the left and preferably, close to centre of the image screen

• For Scoring Multiple Cells:

Select the Comet Cell to score, then drag the ROI over the cell to be scored

• Select 'Store' or click the Right mouse button on the image window

Next Cell

Message "Select the next Replicate Slide, OK to continue, Cancel to repeat last measurement"

Next Slide

Message "Sample complete, data will now be saved, OK to continue, Cancel to repeat last measurement"

If Use Automatic Data Directories is NOT selected, enter a directory and filename for the data to be saved

Message "Do you want to view the image gallery?" Next Sample

Message "Experiment complete" End Experiment.

Komet[®]



3.3.7 Stored Image Experiment Sequence

The following is a pseudo-code presentation of the sequence of Cell, Slide (Replicate) and Sample (Specimen or Donor) cycle defined within a Stored Image KOMET experiment.

Start Experiment

Message "This action will delete your current data. Save it first?"

Message "Do you want to resume an experiment?" (If any are present) If Automatic Data Directories is selected, enter Experiment ID

For M Samples

If Use Automatic Data Directories is selected, enter Sample ID

If Floating Open Button is selected, the Browse for Images Folder dialogue is displayed to select a source folder for images to be used in the analysis.

For N Slides

For P Cells

- Select 'Open' to open an image for analysis
- Position the Cell ROI around the Comet Cell
- Select 'Express' or RIGHT mouse click

Next Cell

Message "Select the next Replicate Slide, OK to continue, Cancel to repeat last measurement"

Next Slide

Message "Sample complete, data will now be saved, OK to continue, Cancel to repeat last measurement"

If Use Automatic Data Directories is NOT selected, enter a directory and filename for the data to be saved

Next Sample

Message "Experiment complete" End Experiment.



3.3.8 Stored Image Experiment Sequence with Database

Start Experiment

Message "This action will delete your current data. Save it first?"

Message "Do you want to resume an experiment?" (If any are present) Enter Experiment ID (entering other setup parameters is optional)

Message informing the user of the path to the database For M Samples

If Use Automatic Data Directories is selected, enter Sample ID

If Floating Open Button is selected, the Browse for Images Folder dialogue is displayed to select a source folder for images to be used in the analysis.

For N Slides

For P Cells

- Select 'Open' to open an image for analysis
- Position the Cell ROI around the Comet Cell
- Select 'Express' or RIGHT mouse click

Next Cell

Message "Select the next Replicate Slide, OK to continue, Cancel to repeat last measurement"

Next Slide

Message "Sample complete, data will now be saved, OK to continue,

Cancel to repeat last measurement"

If Use Automatic Data Directories is NOT selected, enter a directory and filename for the data to be saved

Message "Do you want to view the image gallery?"

Next Sample

Message "Experiment complete" End Experiment.



3.3.9 Data Management

If **Use Automatic Data Directories** is selected in the Comet Options, on completing each sample the data is automatically saved. The directory and filename are saved as follows:

C:\Andor\Komet CAMERAData\UserID\Date\ExperimentID\SampleID.XLD

where **UserID** is the User Name specified on starting Komet, **Date** is in the format, year-month (e.g. 2014-07), **ExperimentID** is the ID given by the user to the experiment, when prompted after clicking on the Experiment button. The **SampleID** is the ID given to the Sample by the user, when prompted at the start of that slide and **CAMERA** is the type of camera your version of Komet supports; Meteor, PCO, etc.

The default directory can be changed by selecting the **Auto Save Path** via the **Browse** button on the Comet Options interface or the Set Save Path option on the Komet Control Window File Menu (**Section 4.5.1.2**). Experimental Data will then be saved to the following destination:

Auto Save Path \UserID\Date\ExperimentID\SampleID.XLD where **Auto Save Path** is the destination path selected. If **Use Database** is selected the database will be saved in the following folder:

C:\Andor\KometCAMERAData\UserID\Date\ExperimentID\ ExperimentID.MDB

If **Use Automatic Data Directories** is not selected, a standard **Save Data File** dialogue window will be opened, in which you can specify a directory and filename to which to save your data. The data is saved as a tab delimited *.XLD file, which can be read in Microsoft Excel. If **Use Database** is selected the database will be saved to the folder specified in the Comet Options: **Database Path**\ DatabaseName.MDB

By default, Automatic Data Directories is selected.



3.4 Calibrating the System

Before you use Komet to analyse images from file or live from a microscope, you should calibrate the system. This will ensure that data referring to DNA migration are computed in real distances, i.e. micrometers. If you do not calibrate the system, these distances will be computed in units of pixels, making comparison between labs more difficult.

If you are scoring from file make sure you use an image file of a calibration device, such as a graticule or haemocytometer, which has been captured with the camera settings you are using for image capture.

If you are scoring live from the microscope, you must use a calibration device (graticule or haemocytometer) with the system set up to image with the correct objective in place.

When running Komet for the first time you will notice that there is a Default calibration in the Calibration Manager. This cannot be deleted or edited, but will be replaced with the new calibration you setup for your system. From here you can create (**New** Button), **Edit**, **Delete** and **Close** calibrations (Figure 11).

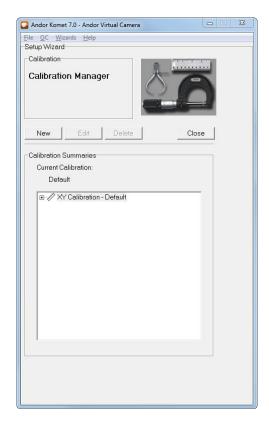


Figure 11. The Calibration Manager



The Calibration Wizard guides you through the process of creating calibrations. It is accessible from the Wizards Menu on the Komet Control Window, or through the Comet Protocol Manager when creating or editing a comet protocol.

The first step of the calibration wizard prompts the user to enter a name for the new calibration and to select the type of calibration required (Figure 12). Enter a name for the Calibration that clearly describes the objective that is being calibrated into the text box, as this calibration will be available to all users

🖸 Andor Komet 7.0 - Andor Virtual Camera	
<u>File QC Wizards</u> <u>H</u> elp Setup Wizard	
TXY Calibration	Personal start and the first data persons
Set up calibration settings:	
<< Back Next >>	Cancel
-Setup New Calibration	5
Enter a name for the new calibration:	
This name must be unique as it will be used filename when the calibration is saved.	as the
Type of calibration required:	
Live Image	•

Figure 12. Setup New Calibration





There are three methods of calibration available through the Setup Wizard (Figure 13).

Still Image		
Import From	Calibrated Image	_
Still Image		

Figure 13. Select Calibration Type

1. Still Image: Use an image which has been captured with the same camera and optical configuration that you will use for scoring. This will commonly be an image of a haemocytometer or graticule (Section 3.4.1).

2. Live Image: Use an image, Live from the microscope with the same camera and optical configuration that you will use for scoring. This will commonly be an image of a haemocytometer or graticule (Section 3.4.1).

3. Import from Calibrated Image: Use an image that you have previously calibrated and saved as a Andor TIFF file. The Andor TIFF format retains calibration information in the file (**Section 3.4.1**).

Press **Next** to continue.



The Calibration Setup Wizard then prompts you to setup the Calibration Region of the image. The image for calibration should contain a discrete feature of known size along the X and/or Y dimensions. A graticule, haemocytometer or cell counting chamber is usually a good choice.

In the Image Window, define a Region of Interest (ROI), so that it encloses a feature of known X and/or Y dimensions (Figure 14).

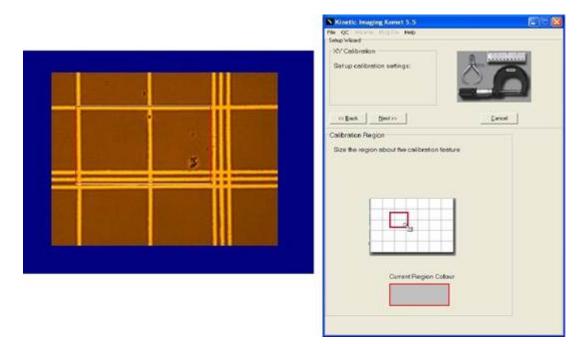


Figure 14. Define the Calibration Region

Note: Clicking on the Current Region Colour rectangle at the bottom of the

Wizard allows you to select a different colour for the region, if it clashes with features on the image (Figures 14).

Press Next to continue.



In the **X** and **Y Distance** text boxes, type the known X and Y dimensions of the feature defined in the previous step, then select the appropriate units (µm or mm) from the drop down list box. An added feature is for the calibration of square pixel CCDs, which can be facilitated from the **Type** drop down list box at the bottom of the tab. When a Square Pixel Type is selected, the user is only allowed to enter a value in one of the distance text boxes, depending upon whether the X or the Y calibration was selected. The value of the other distance is taken to be identical (Figure 15).

S Andor Komet 7.0 - Andor Virtual Camera	×
<u>File QC W</u> izards <u>H</u> elp	
-Setup Wizard	
-XY Calibration	
Set up calibration settings:	
<< Back Next >>	Cancel
-Actual Distance	
Enter the actual distance of the region in X and Y	
X Distance: Units: um Type: XY Calibration	



Press Next to continue.



The final step in the Setup Wizard is to provide a summary of the Calibration Settings (Figure 16).

🖸 Andor Komet 7.0 - Andor Virtual Camera	
<u>File QC Wizards Help</u>	
Setup Wizard	
	R LANGALANCE
Set up calibration settings:	
<< Back <u>N</u> ext >>	Cancel
-Calibration Summary of calibration se	ettings:
Name: NewCal	
Settings	
Type: Live Image	
Description	
Objective Lens	
Lens Magnification:	
Numerical Aperture: Lens Refrac 1.515	tive Index:
Immersion medium	
None RI: 1	
X Resolution	
	ration:
50 200 4	um
Y Resolution	
	ration:
50 200 4	um

Figure 16. Calibration Summary

Enter a **Description** for the calibration into the text box. The XY Calibration will be available to all users, so the description should include any information that will be valuable to the users. Enter the **Lens Magnification** and **Numerical Aperture** of the objective lens you are using into the text boxes.



The **Lens Refractive Index** is set at the default 1.515, but can be changed to correspond to the type of **Immersion Medium** you choose from the drop down list (Figure 16).

Press Next to Continue.

The XY Calibration is now complete and all calibrations are listed at the bottom of the dialogue. Clicking on the plus sign to the left of the calibration name will display summary information for that calibration (Figure 17).

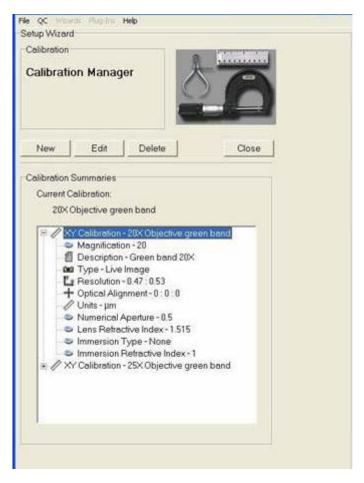


Figure 17. Calibration Summary overview

All of these calibrations are now available to every user when creating a new Comet Protocol. Press **Close** to exit the Calibration Manager.



3.4.1 Calibration Hints

3.4.1.1 Haemocytometer

In most laboratories, a haemocytometer is available for cell counting. A haemocytometer makes a very good calibration device. In the centre of the device are a series of small squares, which can be used to calibrate your microscope/imaging system, whether using live or file based scoring. You will need to image the device in transmission on the central small squares, select XY Calibration and make sure you align the sides with the X and Y axes of the image (Figures 14 and 15). If the device is marked 1/400 mm² (or 0.0025 mm²) then these small squares have a side of 50 μ m.

Note: See the *Directions for Use Sheet* that came with your haemocytometer to verify the ruling pattern, as not all haemocytometers are the same.

3.4.1.2 Graticule

If you have only a linear graticule, then image in transmission and align with the X axis (Figure 18) Select the **Square Pixel (Calibrate X)** option in the calibration type drop down list box (Figure 16). Then enter the real X dimension of the region you have selected on the graticule. The Y calibration will be filled in and assumes square pixels, which is correct for most CCD cameras. For measurements relating to DNA migration, the X-axis is the most important to calibrate precisely.

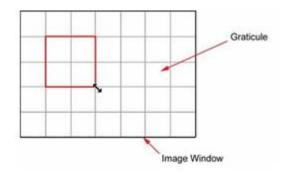


Figure 18. Defining a Calibration Region on the image



3.4.1.3 Binning

If using a digital camera with binning 2x2, you must calibrate using the image graticule at the same binning.

Note: You will need to use a shorter exposure time or turn down the illumination while in transmission.



Komet[®]

3.5 How to Setup Comet Protocols

A Comet Protocol is created from the Comet Protocol Manager. From the **Wizards** menu on the Komet Control Window select **Comet Protocol** or press the **Modify Comet Protocol** button from the Comet Protocol Tab (Figure 19).

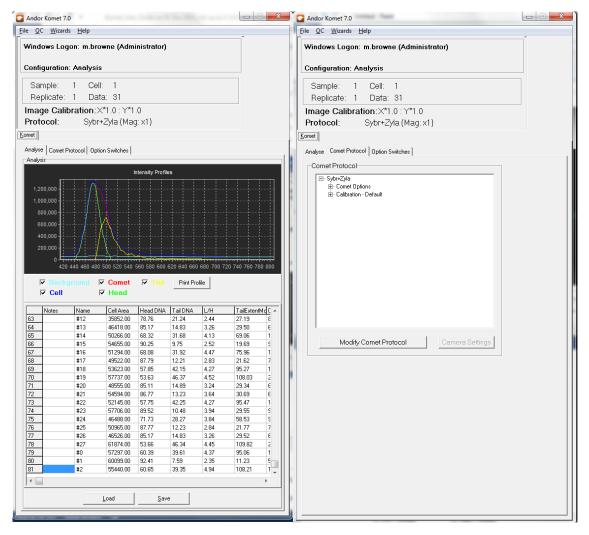


Figure 19. Access of Comet Protocol Manager through Wizards or the Comet Protocol Tab

When running Komet for the first time you will notice that there is a Default Comet Protocol in the Comet Protocol Manager. This cannot be deleted or edited. From here you can create (**New**), **Edit**, **Delete** and **Select** Protocols (Figure 20).





Andor Komet 7.0	J
<u>File QC Wi</u> zards <u>Help</u>	
Setup Wizard	
Comet Protocol	L
Comet Protocol Manager New Edit Delete Select	
Protocol Summaries	
Sybr+Zyla	
(Individual wizard pages can be edited by right-clicking on the component in the tree. Default Protocol cannot be edited)	
⊞- Komet Protocol - Sybr+Zyla	

Figure 20. The Comet Protocol Manager

3.5.1 Creating a New Comet Protocol

Press the **New** button in the Comet Protocol Manager to start. The first step in the Wizard is to give the new protocol a name (Figure 21). It is important to have some sort of naming convention for protocols, to make it easy for users to select a protocol from the manager, for the specific experiment they are about to undertake. This page of the Wizard also requires information about the system calibration. The options available are: **Use existing**, **Edit existing** and **Create new (Section 3.5.2)**.

S Andor Komet 7.0	
<u>File QC Wizards</u> <u>H</u> elp	
Setup Wizard	
Comet setup Enter Protocol name:	
Protocol001	
<< Back Next >> Cance	el
-Select Calibration	
Stored Images	
Apply selected calibration on Open Image	
	is opened, the selected calibration is ap
Select Calibration	1
Zyla_20X	
Calibration Wizard	
Click Calibration Wizard to Create and Edit calibration files.	
and Luit canoration mes.	

Figure 21. The Protocol Name and Calibration Selection



3.5.2 Protocol Calibration Options

Use existing	Select this option and one of the existing calibrations from the drop down list box.
Edit existing	Select this option and one of the existing calibrations you wish to edit from the drop down list box.
Create new	Select this option to create a calibration.
Apply selected calibrat	tion on Open Image
	If checked, the calibration selected in this Comet Protocol will be applied to images that are loaded from the image list or file. If not checked, loaded images will retain their existing calibrations.
Next	This will take the Wizard onto the next step, which will depend
	upon which calibration selection was made. Selecting Edit
	existing or Create new will step through the Calibration Wizard
	(Section 3.4). Selecting Use existing will proceed to the next
	step in the Protocol Setup Wizard.
Cancel	This will take the Setup Wizard back to the start of the Comet
	Protocol Manager.



3.5.3 Camera Settings and Adjustment

Depending upon which camera you are using this part of the Setup Wizard will be slightly different. The **Image Options** are common to all cameras (Figure 22).

Solup Wittend	
Comet setup Enter Protocol name: SYBR Gald	
< Bed >> Qencel Select Camera Settinge Image Options Image Options Image Name: Defoult Polete: Polete: Furnace Image Options	Select Camera Settings Image Options Image Name: Default Palette: Fumace •
Carriers Binning Gain and Other Control $G = 2x^2$ Gain Gain and Other Control $G = 2x^2$ Gain Set to Factory Default Auto Exposure Exposure Time Target 3 Subaration $G = x^2$ $C = 75$ $C = 55$ Set Set Set Set	


Figure 22. Camera Settings and Adjustment Image Options

All acquired images will be stored in memory with an **Image Name** using the prefix entered here (Figure 22). In addition, all acquired images will have the **Palette** or Look Up Table (LUT) selected here, applied to them (Figure 23). This can be altered after acquisition (**Section 4.2.3.1**). We recommend using the Furnace LUT, because it is good for image visibility by the human eye.



3.5.4 Palette or Look Up Tables (LUTs)

Image Name:	Default	
Palette:	Grey	•
Image Control	Negative Camera Red Green Blue Bandit	~
	Furnace Binary	~

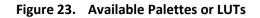

Once one of these Palettes has been chosen from the dropdown list, it is attached to the image currently being displayed and will always be used with that image, unless another palette is later assigned to it. If none of the palettes are selected, the default palette in the current Comet Protocol is used.

Note: The Mapping Range selected will affect the displayed rendering, but not the captured data.

The following standard Palettes are available.

Note that w shape.	ve recommend either Grey or Furnace for good visualization of comet head and tails
Grey	The standard grey level Palette, in which each grey level has an equal proportion of Red (R), Green (G) and Blue (B) components. Thus black (grey level 0) has 0,0,0 and white (grey level 255) has 255,255,255. This is a standard method of visualizing monochrome images.
Negative	The inverse of the grey Palette, where black appears white and vice-versa. Obtained by the "1's complement" of image grey levels.
Camera	An input Palette from a camera, where intensity is proportional to the video signal level, i.e. this is a normal grey scale. However, grey level zero (black) is set to RED=255, B=G=0, while grey level 255 (white) is set to GREEN=255, R=B=0. Use this Palette when setting up image grabbing conditions to detect immediately when saturation or lack of image brightness occurs.



Note: Meteor users must set the image mapping limits to 0 - 253 in order to see the benefit of this LUT.

- **Red** A monochrome palette using only the Red component (intensity 255 the maximum colour brightness). This provides an image shown in a manner which may be reminiscent of viewing a fluorescent image with the appropriate fluorochrome.
- **Green** A monochrome palette using only the Green component (intensity 255 the maximum colour brightness). This provides an image shown in a manner which may be reminiscent of viewing a fluorescent image with the appropriate fluorochrome.
- **Blue** A monochrome palette using only the Blue component (intensity 255 the maximum colour brightness). This provides an image shown in a manner which may be reminiscent of viewing a fluorescent image with the appropriate fluorochrome.
- **Bandit** A colour-banded Palette, ("the banded one") in which groups of 16 grey levels is visualized as the same colour. This is useful for pseudo-colouring and for visualizing and enhancing structural and intensity variation in images. Valuable for comparing changes to the image after a processing operation.
- **Furnace** This Palette is a yellow-brown table in which intensity is matched to brightness along the yellow-brown colour scale. Furnace provides enhancement to the visibility of low contrast image structure.
- **Binary** This palette is not available for Komet 5 or 6.

From the camera settings interface you can adjust the acquisition parameters to obtain the best image quality for the selected calibration. This is done by clicking the **LIVE** button to view a prepared slide in the image window, then adjusting camera settings (Figure 24).





Comet setup Editing Protocol		
SYBR Gold		- 41
LWE		LIVE
<< Back New >		Garcel
elect Camera Setting hage Options	15	
Image Name: 0)etault	
Palette:	umace	<u>•</u>
Intege Control Brightness cu Contrast cu Exposure (ms) [40	146 160 	

Figure 24. Test Camera Settings Using Live

Depending upon which camera is being used, the available controls will differ. Generally for analogue CCD cameras we would recommend an exposure time between 30 to 40 ms and for digital CCD cameras, exposure times in the range of 5 to 40 ms with binning set to 2 x 2. For more details about the Camera control interface, see **Appendix 2**.



3.5.5 Comet Options Settings

Komet®

The Comet Options Settings part of the Protocol Setup Wizard is split into three sections: Processing, Experimental and Operational (Figure 25).

Andor Komet 7.1 - Andor Virtu	al Camera		-	Х
	lelp			
Setup Wizard				
Comet setup Enter Protocol name:				
Protocol001	-			
To test settings click Right Mouse Button in Image Win	dow.			
Then see Key Results field				
<< Back Next >>		Cance		
			51	
Comet Options				
Processing Head Threshold% 10				
	Dealers and Listel	00		
Smoothing Value 1	Background Height	20		
Tail Break Length (um) 5	Saturation Limit	25		
Tail Threshold/1000 10	Saturation Maximum	254		
Experimental Cells per Slide 25	Replicate Slides	2		
	Retain Images in Me			
Samples 2		anory		
C:\Kinetic\KometData\G.J				
Use Database				
(Automatic Data Directory	used)			
Operational			_	
Fluorescence	🔲 Max Head Diamete	ſ		
Auto background	Head Symmetry			
AutoExpress on Store	Tail from Centre of I	Cell		
Stop on Saturation	Interactive Tail			
Floating Live Button	Floating Open Butto	on		
 Head Ellipse On Highlight Scored Cells 	🦳 Flip on Open			

Figure 25. Comet Options Settings

3.5.5.1 Comet Options: Processing

These settings determine how the head and tail regions are defined by the software. They also control how the software will react to saturation levels in the acquired images; therefore indicating when camera adjustments are needed (**Section 4.11.1**).



3.5.5.2 Comet Options: Experimental

These parameters determine the duration of the experiment, whether data is automatically saved to the selected default folder at the end of each Sample, whether acquired images remain in memory where they can be accessed from the Image List, and if the data is automatically stored in a secure database (**Section 4.11.1**).

3.5.5.3 Comet Options: Operational

These determine which processing options will be implemented, and how the experiment will proceed. Ergonomic features such as **Live/Open Floating Buttons** on the interface and **AutoExpress on Store** can increase your experimental throughput (Section 4.11.1).

After specifying the Comet Options, right click on the image to carry out an Express analysis, in order to test the settings. Select **Next** to complete the Comet Protocol Setup Wizard. You should now see the Protocol you set up in the Protocol Summaries (Figure 26).

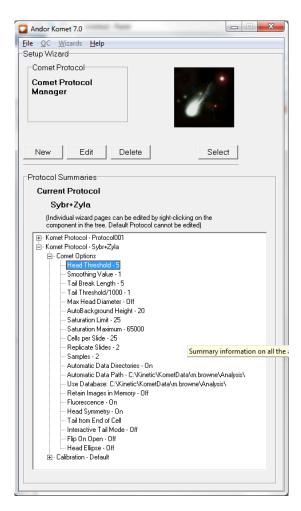


Figure 26. Protocol Summaries



3.5.6 Editing a Comet Protocol

To edit a comet protocol, start the Comet Protocol Manager from the Wizards menu or select the Modify Comet Protocols button under the Comet Protocols tab (Figure 26). Select the Protocol that you wish to edit in the list and press the **Edit** button. This will take you through the whole Setup Wizard, where changes can be made accordingly.

If you wish to edit the Comet Options, Calibration or Camera Settings, then just expand the tree to display the part of the configuration you wish to change. Clicking the right mouse button will now display an **Edit** button which will take you straight to the appropriate page in the Setup Wizard to edit the selected options (Figure 27).

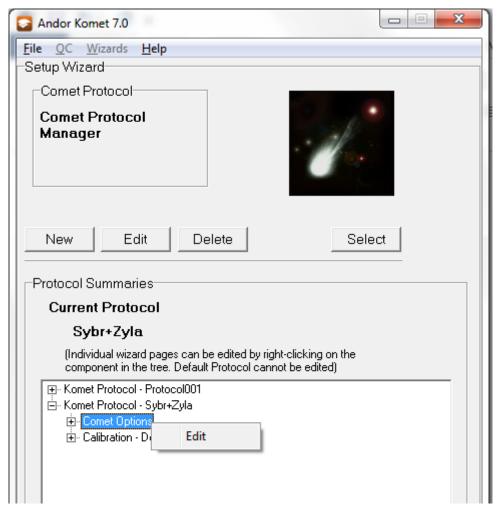


Figure 27. Editing a Comet Protocol



3.5.7 Comet Protocol File Management

Protocols are made available through the Comet Protocol Manager. This allows easy cataloguing and reproduction of experimental configurations without tedious manual recording of all settings. These Protocols are only available to the User that has created them, unlike Calibrations that are available to all Users. However, the procedure in **Section 3.5.4** explains how to share Protocols among Users.

3.5.8 Sharing Comet Protocol Files

To share a Protocol the user will have to copy the **.CAM** and **.CFG** files from one user directory to another using the windows explorer.

For Example:

- Copy C:\Andor\KometCAMERAData\UserIDME \TEST.CFG
- To C:\Andor\Komet CAMERAData\UserID YOU\ TEST.CFG



3.6 Running an Experiment with a Camera

The normal operations involved in the analysis of a Comet cell are described in this section. The description assumes that you have:

- 1) Calibrated the microscope objective lens (Section 3.4).
- 2) Set up the Imaging Options (Section 3.5).
- 3) Set up the Comet Options in the Comet Protocol Setup Wizard (Section 3.5).

Image saturation can be detected and used to halt the analysis, since it will result in estimation errors and affect the quality of data you gather with any imaging system. We recommend the selection of **Stop on Saturation** (Section 4.13.4) in the Comet Options dialogue in all circumstances, unless there is some strong reason to over-ride in a particular case.



3.6.1 Starting an Experiment

To begin an experiment, click on the **Experiment** button on the Comet Toolbar (Figure 28). You will first be asked whether you wish to save the current data in the Komet Control Window. If the data has been saved, either automatically or manually, then select No. The results list will then be cleared of all data.

Live	Interactive	<u>D</u> elete Last	<u>E</u> xpress	Experiment	<u>O</u> pen
------	-------------	---------------------	-----------------	------------	--------------

Figure 28. The Comet Toolbar

If **Automatic Data Directories** is selected in the Comet Options dialogue, you will be asked to enter an ID for the Experiment (Figure 29).

Database Information	
Please complet	e for Database labelling
Project ID:	Test1
Group ID:	1
Individual:	1
Compound or Exposure Regime	EMS
Cell Type:	Hela
*Experiment ID:	,
(^)	Required field)
ОК	Cancel

Figure 29. Experiment ID Text Entry Dialogue Box

When you have entered an Experiment ID and clicked **OK**, a dialogue will be opened asking for an ID for the next Slide for scoring in the Experiment (Figure 30).



Slide ID (max 4 chars)		
Enter an ID for the next Slide		
ЗВ		
ОК	Cancel	

Figure 30. Sample ID Text Entry Dialogue Box

When you have entered an ID for the first Slide of the Experiment in the text box, click on the **OK** button to continue.

In the **Current Context Panel** the Sample, Replicate and Cell values are set to 1, and the Data value to 0. These values are highlighted in RED and are continually updated throughout the experiment (Figure 31). If **Use Automatic Data Directories** is selected in the Comet Options dialogue, the Experiment ID and Sample ID are also displayed.

S Andor Komet 7.0		
<u>File QC W</u> izards <u>H</u>	lelp	
Windows Logon:	m.browne (Administrator)	
Configuration: Ar	alysis	
Sample: 1	Cell: 19 Exp. ID: EMS	
Replicate: 1	Data: 19 Sample ID:3B	
Image Calibration:X*1.0 : Y*1.0		
Protocol:	Sybr+Zyla (Mag: x1)	

Figure 31. The Current Context Panel

3.6.2 Acquire an Image

With your camera connected to the imaging board, click the **Live** button on the Comet Toolbar or the floating **Live** button, if the option has been selected in the current Protocol. A live image will be displayed in the image window.


The **Live** button always uses the conditions set in the Camera Settings current configuration, so once the parameters have been specified there is no need to alter them in the current work session.




3.6.3 Set the Cell / Background ROI


Steer a Comet cell into the Cell Region of Interest (ROI) or drag the ROI box over a cell. The user is able to size and move the Cell ROI manually using the mouse. The boundaries of the Cell ROI should align with the known rectangular boundaries of the cell (Figure 32).

If **Auto background** is selected in the Comet Options dialogue, the background ROI is automatically positioned on the image immediately above or below where the user positions the Cell ROI. If **Auto background** is not selected, the user will have to size and move the Background ROI manually (Section 4.13.10).

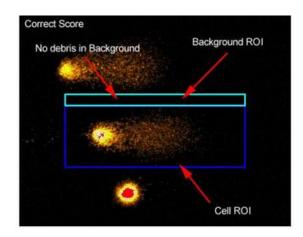

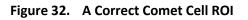

Note: Avoid analysing cells where the Background ROI contains other cells or substantial debris, since this will generate an underestimate in the measured cell intensity parameters (Figure 33).

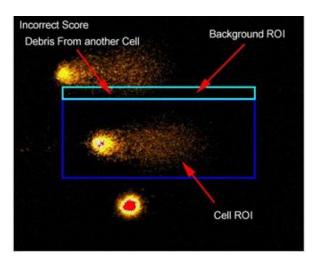





Komet®



3.6.4 Scoring Data

After clicking the **Live** button on the Comet Toolbar, the label changes to **Store**. When you have the cell located in the Cell ROI, click on the **Store** button or right mouse click on the image window to digitize the video image.

If **AutoExpress on Store** is selected in the Comet Options the analysis will be carried out automatically, and the result posted to the Analysis Parameters table. If the result is unacceptable, clicking the **Delete Last** button in the Comet Toolbar deletes the last result in the table. If **AutoExpress on Store** has not been selected, use the **Express** button on the Comet Toolbar or click the right mouse button over the image to perform the analysis.

If the morphology of the cell(s) has some unusual features, click on the **Interactive** button on the Comet Toolbar. You will be prompted to define the tail region using the mouse, after which an express analysis will occur.

Note: If **Interactive Tail** is selected in the Comet Options, you will be prompted with a message below the image window, to manually define the comet tail for each cell (Figure 34). After analysing you can proceed to the next cell.

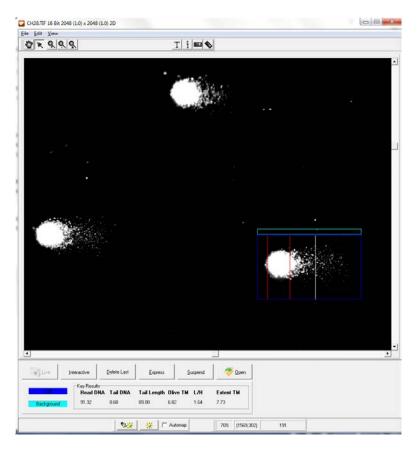


Figure 34. Interactive Definition of the Comet Tail



3.6.5 Experiment Mode - User Prompts

When the system detects that you have scored the number of cells, set in the **Cells per Slide** setting of the Comet Options dialogue, Komet will prompt you to change to the next slide, if the **Replicate Slides** setting is greater than one, and continue the data capture phase. This continues until the correct number of cell data is in the Results list. When all Replicates for the Sample are complete the system will Save the Sample data file. However, this depends on whether or not you have selected **Use Automatic Data Directories**.

If **Automatic Data Directories** is selected in the Comet Options, on completing each Sample the data is automatically saved as follows:

C:\Andor\Komet CAMERAData\UserID\Date\ExperimentID\SampleID.XLD

where **UserID** is the User Name specified on starting Komet, **Date** is in the format year-month (e.g. 2014-07), **ExperimentID** is the ID given by the user to the experiment, when prompted after clicking on the Experiment button. The **SampleID** is the ID given to the Sample by the user, when prompted at the start of that slide and **CAMERA** is the type of camera your version of Komet supports; Meteor, PCO, etc.

The default directory can be changed by selecting the **Auto Save Path** via the **Browse** button on the Comet Options interface or the Set Save Path option on the Komet Control Window File Menu (**Section 4.5.1.2**). Experimental Data will then be saved to the following destination:

Auto-Save Path UserID\Date\ExperimentID\SampleID.XLD where **Auto-Save Path** is the destination path selected.

If **Use Automatic Data Directories** is not selected, a standard **Save Data File** dialogue window will be opened, in which you can specify a directory and filename to which to save your data. The data is saved as a tab delimited *.XLD file, which can be read in Microsoft Excel.

Komet®



3.6.6 Running an Experiment on Stored Images


The normal operations involved in the analysis of a Comet cell are described in this section. The description assumes that you have:

- 1) Calibrated the microscope objective lens (Section 3.4).
- 2) Set up the Comet Options in the Comet Protocol Setup Wizard (Section 3.5).

3.6.7 Starting an Experiment

To begin an experiment, click on the **Experiment** button on the Comet Toolbar (Figure 35). You will be asked whether you wish to save the current data in the Komet Control Window. If data has been saved, either automatically or manually, then select No. The results list will be cleared of all data.

	Live	Interactive	Delete Last	<u>E</u> xpress	Experiment	<u>O</u> pen
--	------	-------------	-------------	-----------------	------------	--------------

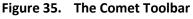

If **Automatic Data Directories** is selected in the Comet Options dialogue, you will be asked to enter an ID for the Experiment (Figure 36).

Database Information					
Please complete for Database labelling					
Project ID: Test1					
Group ID: 1					
Individual: 1					
Compound or EMS					
Cell Type: Hela					
*Experiment ID:					
(* Required field)					
OK Cancel					

Figure 36. Experiment ID Text Entry Dialogue Box

When you have entered an Experiment ID and clicked **OK**, a dialogue will be opened asking for an ID for the next Sample in the Experiment (Figure 37).



Slide ID (max 4 chars)				
Enter an ID for the next Slide				
3B				
ОК	Cancel			

Figure 37. Sample ID Text Entry Dialogue Box

When you have entered a Sample ID click on the **OK** button to continue. If you have the **Floating Open Button** Option selected in your current Comet Protocol (**Section 4.13.11**), the **Browse for Folder** dialogue will prompt you to select the folder containing your images for the first Sample (Figure 38). On pressing the **Open** button, the next image in the folder will automatically open for analysis.



Figure 38. Browse for Images Folder Dialogue Box

In the **Current Context Panel** the Sample, Replicate and Cell values are set to 1, and the Data value to 0. These values are highlighted in RED and are continually updated throughout the experiment (Figure 39). If **Use Automatic Data Directories** is selected in the Comet Options dialogue, the Experiment ID and Sample ID are also displayed.


S Andor Komet 7.0	the free				×	
<u>File QC W</u> izards <u>H</u> elp						
Windows Logon: m.browne (Administrator)						
······································						
Configuration: Analysis						
Sample: 1 Cell: 23 Exp. ID: EMS						
Replicate: 1						
Image Calibration:X*1.0:Y*1.0						
Protocol:	Sybr+Zyla (M	lag:x1)				

Figure 39. The Current Context Panel



3.6.8 Set the Cell / Background ROI

Position the cell ROI around the comet cell making sure the boundaries of the Cell ROI aligns with the known rectangular boundaries of the cell (Figure 40). The user is able to size and move the Cell ROI manually using the mouse.

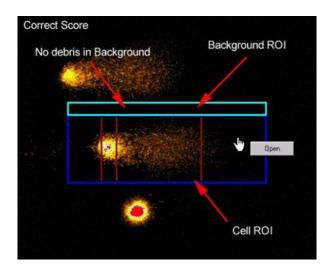



If **Auto background** is selected in the Comet Options dialogue, the background ROI is automatically positioned on the image immediately above or below where the user positions the Cell ROI. If **Auto background** is not selected, the user will have to size and move the Background ROI manually.

Avoid analysing cells where the background region of interest contains other cells or substantial debris, since this will generate an underestimate in the measured cell intensity parameters.

.

Komet®



3.6.9 Scoring Data

Press the **Express** button on the Comet Toolbar or click the right mouse button over the image for the analysis to be carried out. The results will be posted to the parameters table in the Analysis Dialogue. If the result is unacceptable, click on the **Delete Last** button on the Comet Toolbar deletes the last result in the table.

If the morphology of the cell(s) has some unusual features that you wish to score, click on the **Interactive** button in the Comet Toolbar. You will be prompted to define the tail region using the mouse, after which an express analysis will occur (Figure 40).

Note: If **Interactive Tail** is selected in the Comet Options, you will be automatically prompted to manually define the comet tail for each cell.

After analysing the comet cell you can proceed with scoring, by moving the cell ROI, with the mouse, to the next comet in the image window.

3.6.10 Experiment Mode - User Prompts

When the system detects that you have scored the number of cells, set in the **Cells per Slide** setting of the Comet Options dialogue, Komet will prompt you to change to the next slide, if the **Replicate Slides** setting is greater than one, and continue the data capture phase. This continues until the correct number of cell data is in the Results list. When all Replicates for the Sample are complete the system Komet will Save the Sample data file. However, this depends on whether or not you have selected **Use Automatic Data Directories**.

If **Use Automatic Data Directories** is selected in the Comet Options, on completing each sample the data is automatically saved as follows:

C:\Andor\Komet CAMERAData\UserID\Date\ExperimentID\SampleID.XLD

where **UserID** is the User Name specified on starting , **Date** is in the format year-month (e.g. 2014-07), **ExperimentID** is the ID given by the user to the experiment, when prompted after clicking on the Experiment button. The SampleID is the ID given to the Sample by the user, when prompted at the start of that slide, and **CAMERA** is the type of camera your version of Komet supports; Meteor, PCO, etc.

The default directory can be changed by selecting the **Auto Save Path** via the **Browse** button on the Comet Options interface or the Set Save Path option on the Komet Control Window File Menu (Section 4.5.1.2).



Experimental Data will then be saved to the following destination:

**Auto-Save Path ** UserID\Date\ExperimentID\SampleID.XLDwhere

Auto-Save Path is the destination path selected.

If **Use Automatic Data Directories** is not selected, a standard **Save Data File** dialogue window will be opened, in which you can specify a directory and filename to which to save your data. The data is saved as a tab delimited *.XLD file, which can be read in Microsoft Excel.

3.6.11 Running an Experiment Using a Database

The operations involved in a database enabled live analysis of a Comet cell are described in this section. The description assumes that you have:

- 1) Calibrated the microscope objective lens (Section 3.4).
- 2) Set up the Imaging Options (Section 3.5).
- 3) Set up the Comet Options in the Comet Protocol Setup Wizard (Section 3.5).
- 4) Selected Use Database in the Comet Experimental Options (Section 3.5.1.3).

Image saturation can be detected and used to halt the analysis, since it will result in estimation errors and affect the quality of data you gather with any imaging system. We recommend the selection of **Stop on Saturation (Section 4.13.4**) in the Comet Options dialogue in all circumstances, unless there is some strong reason to over-ride in a particular case.



3.7 Starting an Experiment

To begin an experiment, click on the **Experiment** button on the Comet Toolbar (Figure 41). You will be asked whether you wish to save the current data in the Komet Control Window. If the data has been saved, either automatically or manually, then select No. The results list will then be cleared of all data.

Live Interactive Delete Last Expression	ss Experiment 🧖 Open
---	----------------------

Figure 41. The Comet Toolbar

If **Use Database** is selected in the Comet Options, you can assign ID values to several parameters for future reference (Figure 42). The Experiment ID is the only required field; the other fields are optional. The Experiment ID name is given to the database file, and if **Use Automatic Data Directories** is selected, the name is also used in the Data Path and .XLD data. The remaining fields are optional. This information is recorded on the Setup tab when viewing the database in the Database Viewer.

te rer e ande doo rabenning	Please complet
Komet 04	Project ID:
:	Group ID:
:	Individual:
	Compound or Exposure Regime
Liver	Cell Type:
Expt 01 LV BC 2-25-04	*Experiment ID:
Liver	

Figure 42. Experiment ID Text Entry Dialogue Box





If **Use Automatic Data Directories** is selected, after entering the Experiment ID you will be prompted to enter an ID for the next Sample in the Experiment (Figure 43).

iample ID	
Enter an ID for th	e next Sample
LV02BC	
OK	Cancel

Figure 43. Sample ID Text Entry Dialogue Box

Enter the Sample ID then press **OK** to continue. A message is then displayed with the information about the location of the database (Figure 44).

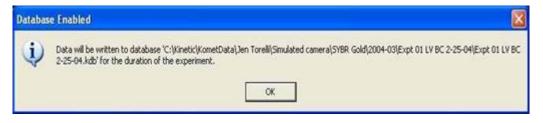


Figure 44. Database Path Message Box

In the **Current Context Panel** the Sample, Replicate and Cell values are set to 1, and the Data value to 0. These values are highlighted in RED and are continually updated throughout the experiment (Figure 45). If **Use Automatic Data Directories** is selected in the Comet Options dialogue, the Experiment ID and Sample ID are also displayed.

Andor Komet 7.0		ſ				
<u>File QC W</u> izards <u>H</u>	elp					
Windows Logon: m.browne (Administrator)						
······································						
Configuration: Analysis						
Sample: 1	Cell: 23 Exp. ID: EMS					
	Data: 23 Sample ID:3B					
Image Calibration:X*1.0 : Y*1.0						
Protocol:	Sybr+Zyla (Mag: x1)					

Figure 45. The Current Context Panel

.



3.7.1 Acquire an Image

With your camera connected to the imaging board, click the **Live** button on the Comet Toolbar or the floating **Live** button, if the option has been selected in the current Protocol. A live image will be displayed in the image window.

The **Live** button always uses the conditions set in the Camera Settings current configuration, so once the parameters have been specified there is no need to alter them in the current work session.

3.7.2 Set the Cell / Background ROI

Steer a Comet cell into the Cell Region of Interest (ROI) or drag the ROI box over a cell. The user is able to size and move the Cell ROI manually using the mouse. The boundaries of the Cell ROI should align with the known rectangular boundaries of the cell (Figure 46).

If **Auto background** is selected in the Comet Options dialogue, the background ROI is automatically positioned on the image immediately above or below where the user positions the Cell ROI. If **Auto background** is not selected, the user will have to size and move the Background ROI manually.

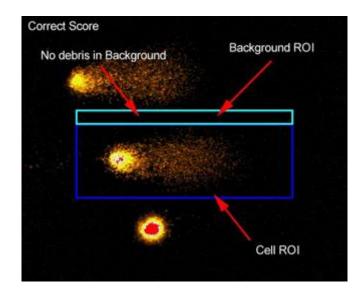


Figure 46. A Correct Comet Cell ROI

Avoid analysing cells where the background region of interest contains other cells or substantial debris, since this will generate an underestimate in the measured cell intensity parameters.

Komet®



3.7.3 Scoring Data

After clicking the **Live** button on the Comet Toolbar, the label changes to **Store**. When you have the cell located in the Cell ROI, click on the **Store** button or right mouse click on the image window to digitize the video image.

If **AutoExpress on Store** is selected in the Comet Options the analysis will be carried out automatically, and the result posted to the Analysis Parameters table. If the result is unacceptable, clicking the **Delete Last** button on the Comet Toolbar deletes the last result in the table. If **AutoExpress on Store** has not been selected, use the **Express** button on the Comet Toolbar or click the right mouse button over the image to perform the analysis.

If the morphology of the cell(s) has some unusual features, click on the **Interactive** button on the Comet Toolbar. You will be prompted to define the tail region using the mouse, after which an express analysis will occur.

Note: If **Interactive Tail** is selected in the Comet Options, you will be prompted with a message below the image window to manually define the comet tail for each cell (Figure 46). After analysing you can proceed with scoring.

3.7.4 Experiment Mode - User Prompts

When the system detects that you have scored the number of cells, set in the **Cells per Slide** setting of the Comet Options dialogue, Komet will prompt you to change to the next slide, if the **Replicate Slides** setting is greater than one, and continue the data capture phase. This continues until the correct number of cell data is in the Results list. When all Replicates for the Sample are complete the system will Save the Sample data file. However, this depends on whether or not you have selected **Use Automatic Data Directories**. If **Use Automatic Data Directories** is selected in the Comet Options, on completing each Sample the data is automatically saved as follows:

C:\Andor\Komet CAMERAData\UserID\Date\ExperimentID\SampleID.XLD

where **UserID** is the User Name specified on starting Komet, **Date** is in the format year-month (e.g. 2014-07), **ExperimentID** is the ID given by the user to the experiment, when prompted after clicking on the Experiment button. The **SampleID** is the ID given to the sample by the user, when prompted at the start of that slide, and **CAMERA** is the type of camera your version of Komet supports; Meteor, PCO, etc.

Komet®



The default directory can be changed by selecting the **Auto Save Path** via the **Browse** button on the Comet Options interface or the Set Save Path option on the Komet Control Window File Menu (**Section 4.5.1.2**). Experimental Data will then be saved to the following destination:

Auto-Save Path\ UserID\Date\ExperimentID\SampleID.XLDwhere Auto-Save Path is the destination path selected.

If **Use Automatic Data Directories** is not selected, a standard **Save Data File** dialogue window will be opened, in which you can specify a directory and filename to which to save your data. The data is saved as a tab delimited *.XLD file, which can be read in Microsoft Excel.

The data is also saved to a password protected Microsoft Access database. The **Andor Database Viewer** is the only method of reviewing the data contained in a database generated during a Komet analysis.

At the end of each treatment unit or individual (usually two replicate slides to improve statistics) the user is given the option to delete the last scored cell and score another with the following dialogue. This allows a poor measurement to be replaced even if it is the last score in the current unit.

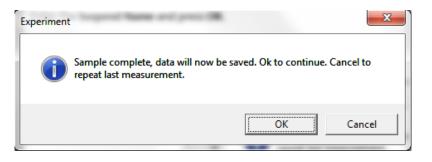


Figure 47. Continue or repeat experiment message box

Once the user is happy with the acquired data, the system provides an opportunity to review all the comets scored in the current sample.



Figure 48. View Image Gallery Message Box





Selecting **Yes** opens the Image Gallery (Figure 49). Exiting the Image Gallery continues the Experiment.

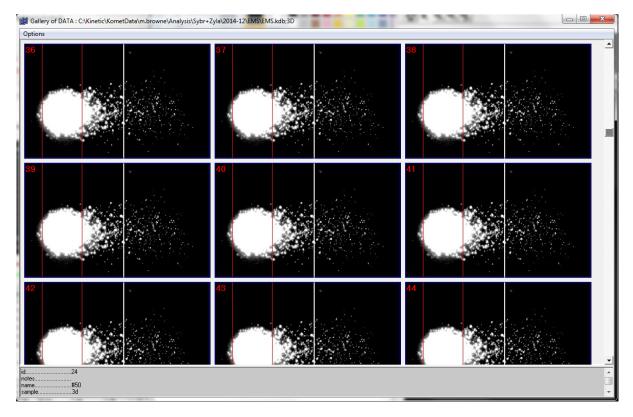


Figure 49. Image Gallery View in Database Mode. Note data per cell is visible at the bottom of the dialog in the Split panel.

When the Experiment Scoring is complete, Komet asks allows the user to write the Dataset directly to a CD/DVD for sharing or archive. If Yes make sure to insert a Clean writeable CD or DVD into the CD or DVD drive on the PC. Then Komet will write the dataset and the Database Viewer software onto the CD/DVD.

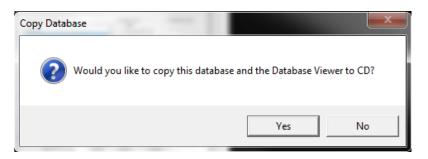


Figure 50. The Copy Database message window

.



Komet[®]

3.7.5 How to Suspend an Experiment

On pressing the **Abort** button on the Comet Toolbar during an experiment, the user will be shown the suspend dialogue (Figure 51).

E	Experiment	nt 📃	
	i	This action will end your current experiment! Do you wish to suspend the experiment for completion at a later date?	
		Yes No Cancel	

Figure 51. The Suspend / Abort Dialogue Box

Yes Suspend the experiment for completion at a later date. You will be asked to enter a name and description for the suspended experiment (Figures 52).

No Abort the current experiment.

Cancel Resume the current experiment.

If you have **Use Automatic Data Directories** selected in the Comet Options the Experiment ID will be suggested as a default name; you may enter a different name if you wish. Similarly, if **Use Automatic Data Directories** is not selected and **Use Database** is selected, the database name will be suggested (Figure 52, left).

Enter the Suspend Name and press OK.

Suspend Name	Notes
Enter name for the suspend state	Enter text to describe experiment
Lunch break	EMS test experiment
OK Cancel	OK Cancel

Figure 52. The Suspend Name and Notes Dialogue

The **Notes** dialogue box describing an experiment can be useful for quick reference (Figures 52, right). Enter the information in the Notes dialogue and press **OK**. The experiment is now suspended and you can close the software or even shut down your computer. The experimental settings will be remembered and you can continue at any time from the same point.



Komet[®]

3.7.6 How to Resume an Experiment

To **Resume** an experiment that has previously been suspended, press the **Experiment** button on the Comet Toolbar. If there are no suspended experiments, a new experiment will start as normal. If there are suspended experiments, you will be asked whether you wish to Resume a suspended experiment (Figure 53).

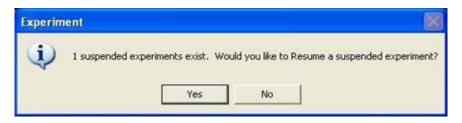


Figure 53. The Resume a Suspended Experiment Dialogue Box

- **No** A new experiment will begin.
- **Yes** You will be prompted to select the suspended experiment that you wish to resume from a dialogue list showing all suspended experiments (Figure 54).

n State	
on state to resume:	
25-04 2004 10:30:16 AM	
	on state to resume:

Figure 54. The Suspended Experiment List Box

.



Click on the experiment you wish to resume and press the **Select** button. You will be prompted to ensure you have selected the appropriate objective lens if acquiring live images (Figure 55).

Comet (ptions			
<u>.</u>	The magnification for this Proto	ocol is x20. Please e	nsure you have the	correct objective selected before going Live.
		ОК	Cancel	

Figure 55. The Magnification Prompt Dialogue Box

After pressing **OK** the experiment will resume at the point where it was suspended (Figure 56). The Comet Protocol used for that particular experiment will be loaded and scoring can resume.

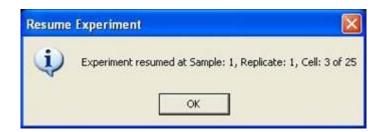


Figure 56. The Resume Dialogue Box

.





SECTION 4: KOMET FUNCTION REFERENCE

4.1 The Image Window

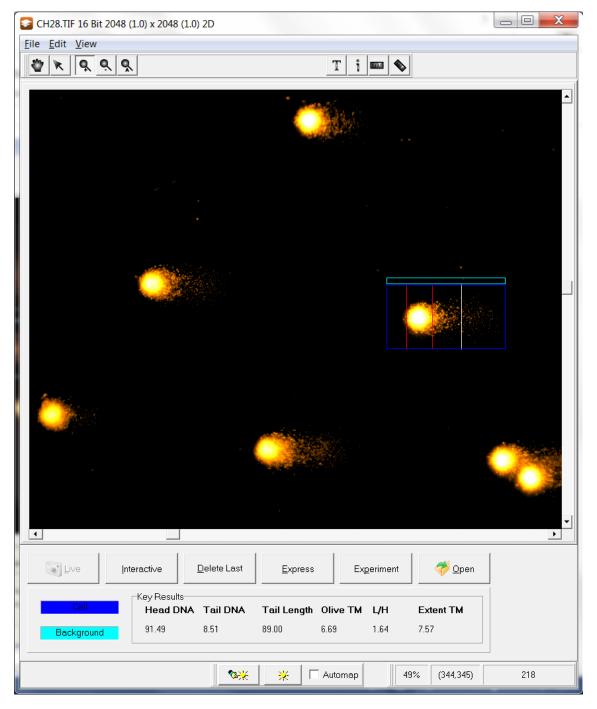


Figure 57. The Image Window

The Image Window is where you view Comet cells (Figure 57). Open the images into the Image Window or go Live, position the Cell ROI around a Comet cell and then perform a comet analysis.



4.2 The Image Window Menus

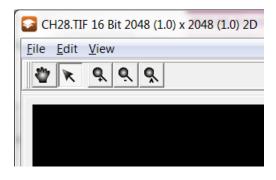


Figure 58. The Image Window Menu Bar

4.2.1 Image Window File Menu

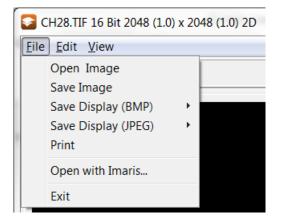


Figure 59. The Image Window File Menu Expanded

4.2.1.1 File Menu: Open Image

When **Open Image** is selected (Figure 59), a digital image is read into memory from disk from a standard Load File dialogue. It will be displayed in the Image Window, ready for analysis. This is equivalent to clicking on the **Open** button on the Comet Toolbar.

Komet has been equipped to Read and Write a wide range of proprietary and industry standard image file formats, as well as non-standard rasterised images. The image file to be read may have any of the following extensions. They are identified by the manufacturer who created the format and, in parenthesis, the product in which the format first appeared (to the best of our knowledge at time of going to press).

Komet®



File Types Supported by Komet are:

.AVI	Microsoft Windows Audio / Video Interleave (first frame)
PIC	Biorad MRC Confocal format
.BMP	Microsoft (Windows bitmap)
.OGB	Olympus GB200 confocal
.IMG	Andor (Fenestra), Kontron (IBAS),
.ICS	Image Cytometry Standard (open source)
.JPG	JPEG Compressed industry standard
.LSM	Zeiss early LSM format (circa 2001)
.MVX	Indec Systems (Microvoxel)
.ANI	Indec Systems (Microvoxel) (loads first frame)
.MIF	Micro-image format – originator unknown.
.PMI	Photometrics (Cooled CCD camera driver software)
.SIF	NASA Scientific Image Format (Andor supported)
.STK	Molecular Devices – Metamorph – limited
.TGA	Targa (Truevision)
.TIF	Tagged Image File Format
UserDef	Non-standard rasterised images
.IMG	Noesis (Visilog)

To select the file type to read, scroll through the options in the **Files of type** drop down list box. Images of the selected type that are in the current directory will be listed in the File list box (Figure 6.).





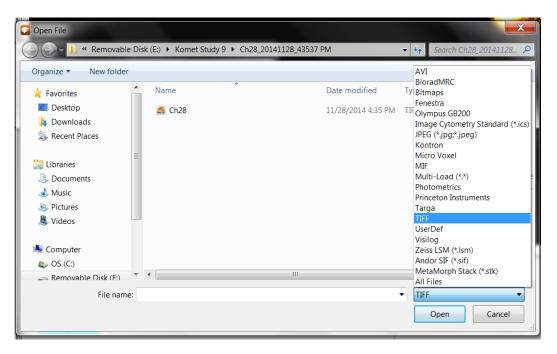


Figure 60. Standard Windows File Open Dialogue Box

4.2.1.2 File Menu: Save Image

When **Save Image** is selected (Figure 61), the current digital image being displayed in the image window can be saved to disk. A standard Windows Save dialogue is used to specify the file name and the desired destination folder. Images can be saved in the selected format from the following list: TIF, IMG, PIC, BMP, PMI, TGA, MVX, ANI, and AVI. Selection of the File Save format is performed in a way similar to Opening Image files, as described in **Section 4.2.1.1**. Saving in TIFF, JPG and BMP formats is recommended for export of images to any Desktop Publishing (DTP) or image archive system.

.



4.2.1.3 File Menu: Save Display (BMP) / Save Display (JPEG)

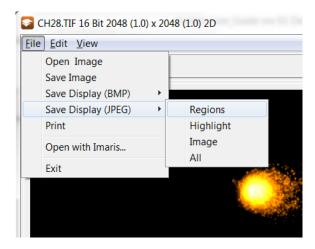


Figure 61. Save Display (BMP)/ (JPEG) Menu Box

The **Save Display (BMP)/ (JPEG)** menu item will save the display options in a Bitmap or Jpeg image format (Figure 61).

Regions	Saves the currently displayed regions (Cell, Background or Calibration) on
	the image; the rest of the image will be blank.
Highlight	Not applicable
Image	Saves the currently displayed image without regions.
All	Saves the currently displayed image and regions.

If the user chooses Save Display (JPEG) Image, he/she will be prompted with a Save Image View Dialogue box in which the file can be saved to a specific folder. Then the user is prompted with a JPEG compression dialogue box (Figure 62).

Select OK to save the image.

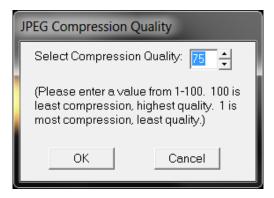


Figure 62. The JPEG Compression Quality Dialogue Box



4.2.1.4 File Menu: Print

Selecting Print from the File Menu (Figure 58), will display the Andor Print Preview dialogue box

(Figure 63).

TeeChart Print Preview		-			
Printer: \\NC-DC-01.ando	rtech.net\Brotł 💌	<u>S</u> etup	Print	Clos	se
Oriențation: Portrait Landscape Margins (%) 15 ÷ 20 ÷ 20 ÷ 15 ÷					
Reset <u>M</u> argins					
Proportional					

Figure 63. The Print Dialogue Box

Printer	A drop-down list of the printers available on your system.
Setup	Displays the Printer Setup dialogue of the selected printer.
Print	Press this button to print the contents of the main window.
Close	Closes the Print Preview Dialogue.
Orientation	The page can be printed in either Portrait or Landscape
	format.
Margins (%)	Page Margins can be adjusted as a % of the page size.
Reset Margins	If margins have been altered they are reset to defaults.
View Margins	Toggles the visibility of the margins (Not Printable).

Komet®



Detail Sets the dpi of the printout.

Proportional Retains the aspect ratio of the image when checked.

4.2.1.5 File Menu: Exit

Selecting **Exit** from the File Menu (Figure 58) will ensure all files used during the work-session are properly closed. The data currently in the Results table in the Analysis dialogue is saved to the file LastSession.XLD. The data is saved to the folder from which you are running Komet and also to the Komet User folder:

C:\Andor\Komet CAMERAData\UserID.

The current system setup is also saved to the Registry. It is reloaded the next time that user starts Komet, so that the system setup is the same as for the previous session.

4.2.2 Image Window Edit Menu

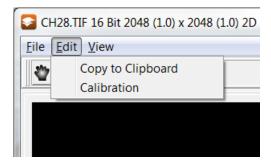


Figure 64. Copy to Clipboard

4.2.3.1 Edit Menu: Copy to Clipboard

Choosing **Copy to Clipboard** (Figure 64) copies the current contents of the image window to the clipboard, where it can be pasted into other windows applications, for example word processing or DTP documents.

4.2.3.2 Edit Menu: Calibration

Choosing **Calibrations** provides a list of Calibration settings to apply to the current image.



4.2.4 Image Window View Menu

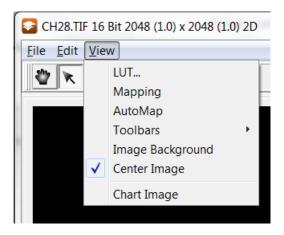


Figure 65. The Image Window View Menu Expanded

4.2.4.1 View Menu: LUT...

Selecting **LUT...** gives the user the option of changing the Look Up Table (LUT) or palette of the image being displayed in the Image Window (Figure 66).

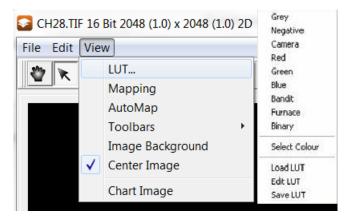


Figure 66. The View Menu LUT Expanded

Note: This does not change the setting of the LUT or palette in the Comet Protocol.

.



4.2.4.2 View Menu: Mapping

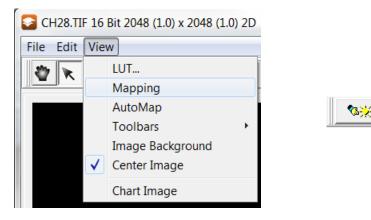


Figure 67. Figures 4.11 The View Menu Mapping option and Mapping shortcut key

Select **Mapping** from the View Menu (Figure 67, left) or select the **Mapping** shortcut key (Figure 67, right) to open the Image Contrast dialogue, in which you can manually set the image mapping (Figure 68). Using the **Max** and **Min** scrollbars, select the intensity values within the image to be better contrasted and press the OK button to accept the current mapping values for all images subsequently opened or grabbed. The example in Figure 68 reflects settings for an 8-bit image.

🗠 Image Contrast	
Options View	
Gamma Gamma Exclude	■ 12360.00 ÷ Max ■ 462.0000 ÷ Min 1.000000 ÷ Step
Gamma Coefficien ,	0.50
Smooth Use Region of Interest	Ok Cancel

Figure 68. The Image Contrast Dialogue Box

This modifies the display of image pixel intensities in the Image Window and is used to improve the contrast of the image to the operator, or to view specific image properties. Only the Image Window display is changed; analysis is not impacted as the underlying image pixel intensities remain the unchanged.



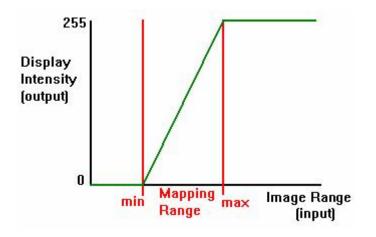


Figure 69. The Image Mapping Function

For display, the intensity at each image pixel is converted to a display intensity according to the mapping. Pixel intensities less than the Mapping Range minimum are displayed as value zero, while pixel intensities greater than the Mapping Range maximum are set to 255. Figure 70 is an example of linear contrast mapping for 8-bit data.

In practice it is recommended that you set the Mapping Range from 0 - 255 (0 - 254 for Meteor Controlled Cameras) while grabbing an image and then experiment with different brightness and contrast mappings in the Camera Settings dialogue (**Appendix 2**) to find the one that best contrasts image features of interest.

For 12 and 16 bit data Imager Mapping is essential and because the display range of monitors is usually in the range of 8 bits per channel, mapping compresses the image grey scales into this range. This means in order to see the wide dynamic range of dim tail and bright head you may need to saturate the display of the comet head by adjusting the Max and Min values. Once way to help visualize wide dynamic range data is to use the Gamma function to provide a non-linear scale. So for example if gamma<1 then contrast of low intensities pixels are boosted while high intensity pixels are slowly increased as shown in Figure 68 where Gamma = 0.5.



4.2.4.3 View Menu: AutoMap

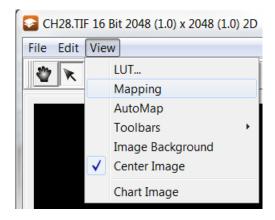


Figure 70. The View Menu AutoMap option and AutoMap shortcut key

Select **AutoMap** from the View Menu (Figure 70) or select the **AutoMap** shortcut key . This will automatically map the current image by analysing a randomized selection of 16K pixels from the current image and setting the display to the max and min of the pixel values.

4.2.4.4 View Menu: Toolbars

The **Toolbars** menu has a Toolbar and Status option (Figure 71). If these are selected (check mark present), then the Image Window Toolbar (Figure 72) and the Image Window Status bar will be present.

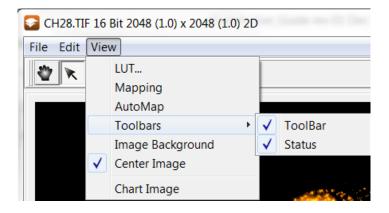


Figure 71. The View Menu Toolbars Expanded

4.2.4.4.1 Toolbars Menu: Toolbar

The buttons available on the **Toolbar** in Komet are the **Pointer** and **Zoom** tools (Figure 72)



Figure 72. The Image Window Toolbar

The **Pointer** tool is used when defining or moving a region (Comet Cell, Comet Background or Calibration Region) on the image window. Note the Hand tool has the same function as the Pointer.

To move a region, position the cursor over the Region of Interest (ROI) box. When the cursor changes to the shape of a hand, hold down the left mouse button and move the mouse. Releasing the mouse button will drop the ROI at that point.

To resize a region position the cursor over an edge or corner so that the scrolling arrows appear. Hold down the left mouse button and move the mouse, dragging the edge or corner in the direction of the scrolling arrow. When the ROI is in the desired position release the mouse button.

Note: When using the right mouse to initiate a Komet analysis the Pointer tool MUST be selected.

Q Q Q N The Zoom tools enables you to zoom, in out and fit the image displayed in the Image Window. This allows you to view the fine detail of features of interest more clearly. Selecting the **Zoom** icon will enable zooming in/out using the right and left mouse buttons. Right Mouse to zoom **OUT** and Left Mouse to zoom **IN** about the point clicked. The current magnification of the image is displayed in the text box to the right of the Mapping/AutoMap shortcut keys on the Status bar (Figure 73). For Example: 50%, 100%, 200% or 120 %.



4.2.4.4.2 Toolbars Menu: Status



Figure 73. The Image Window Status Bar

When **Status** is selected from the Toolbars menu options, a Status Bar will be displayed in the Image Window (Figure 73). The Status Bar displays the manual **Mapping** and **AutoMap** buttons, the image information, such as zoom factor, image position (X,Y coordinates) and pixel intensity value in grey or RGB format. When you position the cursor over the image displayed in the Image Window, the image position coordinates and pixel intensity value changes.

4.2.4.5 View Menu: Image Background

Selecting **Image Background** from the View Menu enables the user to select a background colour for the Image Window (Figure 74). This background colour surrounds the image being displayed in the Image Window.

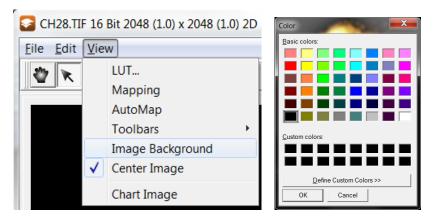


Figure 74. The View Menu Image Background

4.2.5 Image Window Image List Menu

Komet®

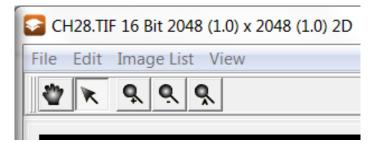


Figure 75. The Image Window Image List Menu

This Image List Menu (Figure 75) is only visible when the current Comet Protocol has Retain Images In Memory (Section 4.12.4) selected in Comet Options. Clicking the Image List menu item opens the Image Manager dialogue (Figure 76).

😡 Image Manager : Virt	ual			
File Edit Comments	Options			
Image Details				
Name : CH28.TIF Path : E:\KOMET STUDY 9 Type : 16 bit grey/pseudo Intensity : 1.00000 lux x : 2048 * 0.322500 : um y : 2048 * 0.322500 : um	I\CH28_20141	128_43537 PM\		
[Created] Date=28/11/2014 Time=4:33:13 PM [Created End] [Grab Parameters] Camera=Zyla Window Left=0 Window Top=0				
Image List				
Image Name	# Date	- Time	Туре	
CH28.TIF1	1 28/11	/2014 - 4:33:13 PM	16 bit grey/ps	seudo
CH28.TIF	2 28/11	/2014 - 4:33:13 PM	16 bit grey/ps	seudo
Close Refresh	Show	Rename Image	Save Image	

Figure 76. The Image Manager Dialogue Box

Acquired and opened images are handled by Andor's Image Manager and are displayed in the Image List. Images under Andor Memory Management are either stored in system RAM or onto a disk partition formatted with NTFS.

Komet[®]



Note: From Windows 98 onwards, users can only use system RAM to store images; therefore, exiting Komet will lose all retained images from memory. Ensure you save any images you wish to keep prior to exiting Komet.

The Image Manager displays Image Details, including the image name, type, intensity and dimensions. It is therefore an excellent method of reviewing recent acquisitions.

When you acquire images by pressing **Live** then **Store** on the Comet Toolbar (**Section 4.2.6**), they can be accessed from the Image List. By default they will be named with a prefix, specified in the **Image Name** parameter in the Comet Setup Wizard (**Section 3.5.1.2**).

Close

Exits the Image Manager.

Show

Displays the selected image in the Image Window.

Rename Image

You can rename the highlighted image (Figure 77).

Rename Image	Constants Constants
Enter new name for image:	1128_43537 PMCH28TIF
ОК	Cancel

Figure 77. The Rename Image Dialogue

Edit the name in the text box as required, and then press **OK**.

Note: This action does NOT save the image.

Save ImageOpens a standard Windows Save File dialogue from which you cansave the currently selected image to disk.

Delete Selected

ImagesAll selected images will be deleted from the Image Manager. Note: If the
image is also saved to disk, the saved image will NOT be deleted.



4.2.6 The Comet Toolbar



Figure 78. The Comet Toolbar

The Comet Toolbar (Figure 78) supplies a rapid and convenient means of controlling the system both during an Experiment (Sections 3.6 - 3.7) and in free mode.

4.2.6.1 Live/Store

Clicking on **Live** switches the system to live image grabbing mode and changes the button label to **Store**. Provided you have the microscope ocular beam splitter correctly set, as you move the specimen about under the microscope you will see a live image of the specimen. When a cell is suitably positioned and focused within the Cell Region of Interest (ROI) click on the **Store** button, or more conveniently click the Right mouse button and the image will be digitized and acquired. If **AutoExpress on Store** is selected in the Comet Options, the Express function will be invoked; equivalent to clicking the **Express** button on the Comet Toolbar (**Section 4.3.4**). The **Store** button will be re-labelled **Live**, indicating the next action it will instigate.

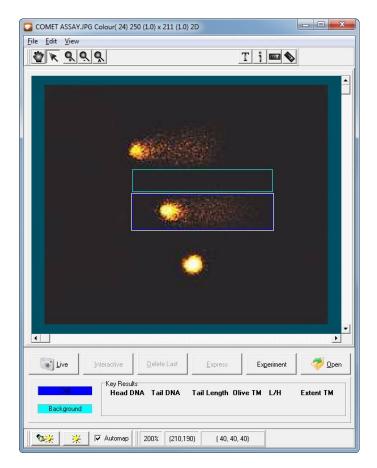


Figure 79. Comet Assay Live view



4.2.6.2 Interactive

Interactive

Cells may exhibit fragmented comet tails or unusual morphology after certain treatment. If the comet tail has a break greater than the value defined for the **Tail Break Length (Section 4.11.1.3)** on the Comet Options, this can lead to estimated tail lengths significantly shorter than the correct value. The **Interactive** function enables the user to define the Tail ROI under such circumstances. The user is prompted to define a ROI using the mouse from the start to the end of the tail (Figure 80). On releasing the mouse button, the interactive Comet Cell is analysed. Selecting the **Interactive Tail** function (**Section 4.13.10**) in the Comet Options will automatically invoke the interactive function for all analyses.

Note: The Pointer tool must be selected to initiate a Komet analysis using the Interactive function.



Figure 80. Defining an Interactive Tail

4.2.6.3 Delete Last

Delete Last

Successful analysis of a comet will insert a row of data into the analysis table in the Komet Window (**Section 4.10**). Pressing the **Delete Last** button will remove the last row of data from this table.



4.2.6.4 Express



Press the **Express** button or right click on the image window to analyse a Comet Cell. This will compute all the parameters and post them to the table on the Analyse tab. With **AutoExpress on Store** selected in the Comet Options, the **Express** button is rarely necessary when analysing live images, since when you click on **Store** (complement of Live) the Express function will be invoked.

4.2.6.5 Experiment/Abort



The **Experiment** button will switch Komet into Experiment mode in which it keeps a track of how many cells have been scored and prompts you when to change slides, save data and begin a new sample.



When you click on the **Experiment** button, the button label changes to **Suspend**. Click on the **Suspend** button if you wish to cancel the Experiment before its completion or Suspend the Experiment to be resumed at a later date

4.2.6.6 Open



The **Open** button can be used to load a stored image file. This has the same function as **Open Image** on the **File Menu.** (Section 4.2.1.1).

4.3 The Key Results Panel

Cel	Key Results Head DNA	Tail DNA	Tail Length	Olive TM	L/H	Extent TM
Background	66.13	33.87	83.02	9.27	5.13	28.12

Figure 81. The Key Results Panel

Results from the latest Experiment / Express Analysis are displayed in the Key Results dialogue for quick reference (Figure 81), as well as in the results table on the Analysis dialogue (**Section 4.11**). The Dark Blue ROI in the Image Window defines the Cell region and the Light Blue ROI in the Image Window defines the Cell region and the Light Blue ROI in the Image Window defines the Section 4.11).

When testing the Comet Options settings (**Section 3.5.5**) the text on the Key Results Panel will be displayed in RED, highlighting to the user the changes in the results.



4.4 The Komet Control Window

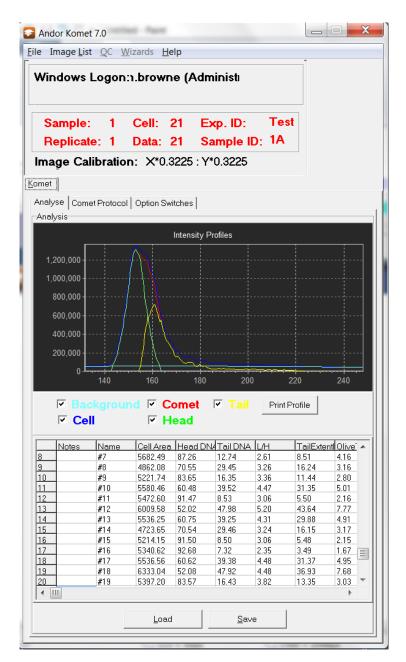


Figure 82. The Komet Control Window

From the Komet Control Window the Analyse tab, the Comet Protocol tab and Option Switches tab are available (Figure 82). See **Sections 4.11 to 4.14** for more information regarding each of the tabs.



4.5 The Komet Control Window Menus

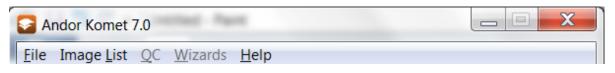


Figure 83. The Komet Control Window Menu

Note: The Komet Control Window Menu Bar will appear in two forms, depending on whether or not Retain Images in Memory is selected. See **Section 4.2.5** for more information regarding Retain Images in Memory.

4.5.1 Komet Control Window File Menu

4.5.1.1 File Menu: Save Image

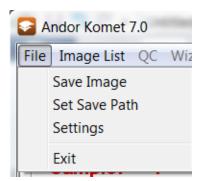


Figure 84. The Komet Control Window File Menu Expanded

4.5.1.2 File Menu: Set Save Path

Selecting **Set Save Path** (Figure 84) brings up a Select Data Path dialogue box, which allows the user to change the path for saving data (Figure 85). Clicking on the Browse button opens a dialog in which a user defined Folder can be chosen as the default path for data.



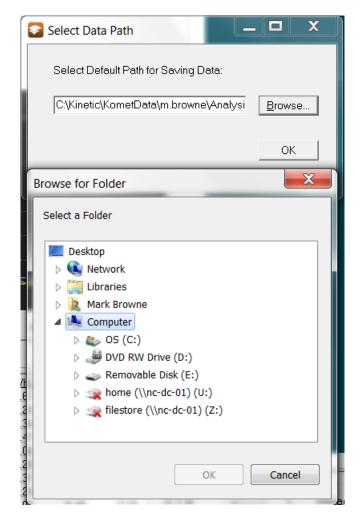


Figure 85. The File Menu Set Save Path Dialogue Box with the Browse Folder dialog open.



4.5.1.3 File Menu: Settings

Set Single File Size Limit	×
Single File Size Limit (1-2000 MB):
640	(MB)
ОК	Cancel

Figure 86. The File Menu Settings Dialogue Box

To set the size limit for a single file, if for example, saving to a CD, select **Settings** from the File Menu. The Set Single File Size Limit dialogue box (Figure 86) will appear and allow the user to select a size limit for the file(s) being saved.

4.5.1.4 File Menu: Exit

Selecting **Exit** from the File Menu ensures all files used during the work-session are properly closed. The data currently in the Results table in the Analysis dialogue is saved to the file LastSession.XLD. The data is saved to the folder from which you are running Komet and also to the Komet User folder:

C:\Andor\Komet CAMERAData\UserID.

The current system setup is also saved to the Registry. It is reloaded the next time that user starts Komet, so that the system setup is the same as for the previous session.

.




4.6 Komet Control Window QC Menu

File	QC W	zards P	lug-Ins	Help
-C	Run			
	User:	Jen To	orelli	(Admin
	Configu	ration	: Sim	ulated
	mage:	ком	ETST	ART.B

Figure 87. The OC Menu Expanded

The QC menu **Run** option (Figure 87) is only available in Komet GLP. This feature is particularly useful for those involved in GLP studies and can be purchased for an additional fee. Contact Andor for further information.

4.7 Komet Control Window Wizards Menu

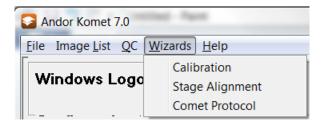



4.7.1 Wizards Menu: Calibration

Select **Calibration** from the Wizards menu to access the Calibration Setup Wizard (Figure 88). The Calibration Setup Wizard will guide the user through the process of calibrating the system (**Section 3.4**).

Note that Stage Alignment is not functional in Komet.

4.7.2 Wizards Menu: Comet Protocol

Select **Comet Protocol** from the Wizards menu to access the Comet Protocol Manager (Figure 89). A Protocol Setup Wizard will guide you through setting up a Comet Protocol (**Section 3.5**). The user will select a calibration, define a camera setup and select a set of Comet Options. These parameters are then used for subsequent analyses.



4.8 Komet Control Window Help Menu

Sandor Komet 7.0	
<u>File</u> Image <u>L</u> ist <u>Q</u> C <u>W</u> izards	<u>H</u> elp
	Help
Windows Logon:bro	Product Activation
	About

Figure 89. The Komet Control Window Help Menu Expanded

4.8.1 Help Menu: Help

Selecting **Help** from the Help Menu (Figure 89) opens the User Guide. The User Guide online Help is in the standard PDF format, familiar from many Windows application packages. This requires Adobe Reader or a third party PDF viewer. Use the bookmarks to menu to navigate through the User Guide.

4.8.2 Help Menu: Product Activation

The Product Activation function allows authorized personnel to activate the software for evaluation. A temporary license file can be created here (or provided by Andor) and placed on the Computer Desktop to enable full operation for a finite period. Selecting this menu will show the current status of the license file if any is shown. Note the license files can be shared with other Andor products e.g. Andor iQ – as will be recognized in the screen shot below.

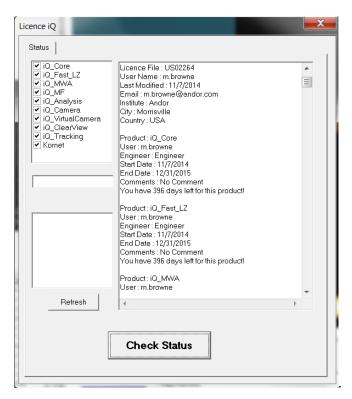


Figure 90. The iQ product Licence menu



4.8.3 Help Menu: About

Pressing **About** from the Help Menu (Figure 90) displays information about the version number and release date of Komet (Figure 91).

Ab	out		X			
	ANDOR an Oxford Instruments company	Andor Bioimaging Division Komet				
		Copyright : Andor Technology 1990-2014				
	Build Time : 14:44:55 No	v 7 2014 Build Number : 7.0.1.40				
	Corporate Headquarters 7 Millennium Way Springvale Business Park Belfast BT12 7AL United Kingdom					
	Tel: +44 (0)28 9023 712	6	APAC			
	row micros	copy_support@andor.com				
		www.andor.com				
		ОК				

Figure 91. The Komet About Information Box





4.9 The Current Context Panel

Andor Komet 7.0)		·				X	
File QC Wizard	s H	elp						
Windows Log Configuration			ie (Ad	lministra	ator)			
Sample: Replicate:						 		
Image Calit Protocol:	orat	ion:×*1	.0 : Y			 		

Figure 92. The Current Context Panel

The User ID, specified on starting Komet (Section 3.1), is displayed in the Current Context Panel, together with the Configuration (Figure 92). Experimental information such as the current Sample number, Cells and Replicates for an experiment is displayed here, together with the number of sets of Data that have been collected. These parameters are specified in the Comet Options Experimental section (Section 4.12).

Click the **Experiment** button in the Comet Toolbar (**Section 4.2.6**) to activate Experiment Mode. The Experiment ID and current Sample ID, input by the user, are also displayed and the experimental information is highlighted in red (Figure 93).

[Andor Komet 7.0	other from			x
	<u>File QC W</u> izards <u>I</u>	<u>H</u> elp			-
	Windows Logon	m.browne (Ad	dministrator)		
	Configuration: A	nalysis			
		-	Fxn ID:	EMS	
	Configuration: A Sample: 1 Replicate: 1	Cell: 23			

Figure 93. The Current Context Panel During an Experiment



4.10 The Analyse Tab

The **Analyse Tab** in the Komet Control Window (Figure 94) allows the user to view the Intensity Profiles and the Individual Analysis Parameters for each cell scored.

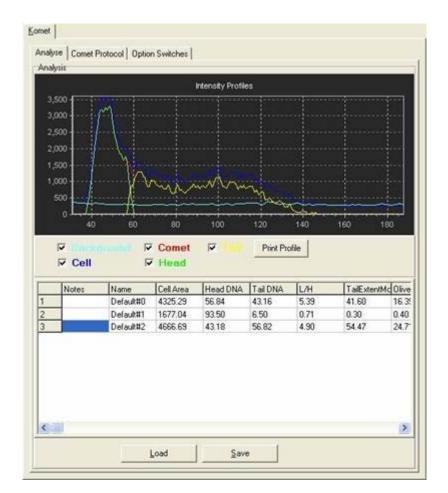


Figure 94. The Komet Analyse Tab



4.10.1 Intensity Profiles

Komet[®]

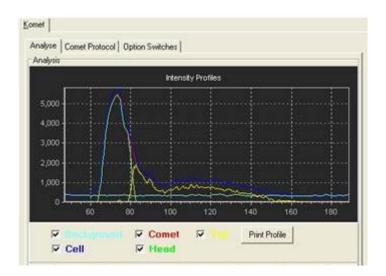


Figure 95. The Komet Analyse Tab Intensity Profiles

The **Intensity Profiles** are displayed as the analysis proceeds (Figure 95). The colour coded check boxes below the black graphic rectangle determine which of the profiles are displayed.

Colour coding is as follows:

Background - Light Blue Cell - Dark Blue Comet - Red Head – Green Tail - Yellow Print Profile – Prints the Comet profile (Section 4.2.1.4)

For a detailed description of these profiles and how they are used for DNA damage analysis see **Section 9**.

.



Komet[®]

4.10.2 Individual Analysis Parameters

The **Individual Analysis Parameters** for each cell is displayed in the results table. Double clicking in the corresponding cell under the **Notes** column enables the user to add comments about that cell (Figures 96).

Cancel

Figure 96. The Individual Analysis Parameters and Notes Dialogue Box

Users are able to Load their data at the beginning of a session and/or Save their data at the end of a session to reload it at a future date, by pressing the **Load** and/or **Save** button.

On exiting from Komet, the Data currently in the analysis table is saved to the file LastSession.XLD, in the directory from which you are running Komet. The data is also saved in the current user directory:

C:\Andor\Komet CAMERAData\UserID

where UserID is the user name entered on starting Komet. The next time that user runs Komet the data will be read from this file.

Data is saved in tab delimited spreadsheets, including Microsoft Comet Analysis Excel Macros interpretation of data (**Section 5**) format, so it can be read in Excel.

.



4.10.3 Comet Protocol Tab

Comet Protocols determine how the software defines the head and tail regions in the comet cell. The current Comet Protocol (Section 3.5) can be viewed from the Comet Protocol Tab (Figure 97). You can create (New), Edit, or Delete a Protocol from the Comet Protocol Manager, which can be accessed from the Modify Comet Protocols button on the Comet Protocol Tab or through the Wizards Menu on the Komet Control Window (Section 4.7).

) E	BR Gold Comet Opt Calibration Camera Sr	- 20K Objecti	ve green b	and	

Figure 97. The Comet Protocol Tab

On starting Komet, the Comet Protocol settings are loaded from the Registry with the values in use the last time the current user exited Komet. If the current user is a new user, the values from the previous Komet session are loaded.

For a full tutorial on creating a Comet Protocol from the Comet Protocol Setup Wizard see **Section 3.5.1**.

The **Camera Settings** button allows the conditions for grabbing and digitization to be established (Figure 97). On starting Komet, the Camera Settings parameters are loaded from the configuration saved to the Registry the last time the current user exited Komet. If the current user is a new user, default values are loaded. You can change the values in the Camera Settings dialogue at any time, but they will not be saved into the current Comet Protocol that is loaded (Figures 98). For more information regarding the camera settings see **Appendix 2**.

Komet®



loguisition/Auxiliary Devices	M		
Canesa Image Control Video Source Accumulation Contract 160 Brightness 128 Exposure (m) Frames [33.3 1 1] Set Window 540 : 430 Standard << Meteor-II Mi6 Version Dox	Acquisition/Auxiliary Devices Camera Image Control Video Source Accumutation Video Source Formal NTSC Video Source Formal Monochrome Device Camera Mode Field Grab Field Grab Channel (C) C 1 C 2 C 3 Sampling X 1 Y 1	Acquisition/Auxiliary Devices Carrere Image Control Video Source Accumulation	X
Standard Video with Meteor II/I	Standard Video with Meteor II/A	Standard Video with Meteor II/1	

Figure 98. The Camera Settings Dialogue: Image Control, Video Source and Accumulation

4.11 Comet Options Settings

Komet®

The Comet Options Settings are divided into three sections: Processing, Experimental and Operational (Figure 99).

	met Options			
Г	Processing			
	Head Threshold%	10		
	Smoothing Value	1	Background Height	20
	Tail Break Length (um)	5	Saturation Limit	25
	Tail Threshold/1000	10	Saturation Maximum	254
Γ	Experimental			
	Cells per Slide	25	Replicate Slides	2
	Samples	2	🔲 Retain Images in f	Memory
	C:\Kinetic\Kom			
	Use Database (Automatic Data		used)	
	Use Database			
	Use Database			
	Use Database Use Database (Automatic Database)	a Directory (used)	
	Use Database (Automatic Database) Operational Fluorescence	a Directory (used) Max Head Diame	
	Use Database (Automatic Database) Operational Fluorescence Auto background	a Directory (1 Store	used) Max Head Diame Head Symmetry	
	Use Database (Automatic Database) (Automatic Database) (Automatic Database) (Automatic Database) (Autobase)	a Directory (used) ☐ Max Head Diame ✓ Head Symmetry ☐ Tail from Centre of	of Cell
	Use Database	a Directory (used) ☐ Max Head Diame ✔ Head Symmetry ☐ Tail from Centre of ☐ Interactive Tail	of Cell

Figure 99. The Comet Options Dialogue Box

4.11.1 Comet Options: Processing



Figure 100. The Comet Options Processing Dialogue Box

.....



4.11.1 .1 Head Threshold %

The default value is 10%.

The **Head Threshold %** value is used to specify the intensity threshold of the head region for density measurements. Pixels with an intensity value above the background by this percentage of the maximum head intensity are included in the Head measurements. By selecting a separate Head Threshold and Tail Threshold (**Section 4.11.1.4**) you can provide tight calipers on the cell head (higher head threshold), while detecting very faint fragments in the comet tail (lower tail threshold). Values can lie in the range 0-90%, the default value being 10%.

4.11.1.2 Smoothing Value

The default value is 1.

The **Smoothing Value** is used to determine the smoothing applied to integrated intensity profiles prior to Head/Tail analysis. It is useful for noisy and/or low intensity data. The default value is 1, but for noisy data a higher value is recommended. Allowable values are in the range 0-20.

4.11.1.3 Tail Break Length

The default value is 5.

The estimation of comet tail length is probably the most difficult of all measurements undertaken in the analysis. It is such a variable characteristic of an intensity profile, as it depends on everything from the toxin under study to the type of cell, and of course, the execution of the experimental protocol.

Under some circumstances, treated cells may exhibit fragmented comet tails or there may be spot noise from crystals in the fluorophore solution. In this case, the intensity profiles used for analysis may show "breaks" (regions where intensity drops below background), even though more DNA fragments lie further from the cell. Breaks can give rise to under-estimation of tail length, while spot noise can have the opposite effect. This control allows you to decide above what limit a break should be considered sufficiently large to terminate a tail measurement. In this case, a 5 micron break will terminate the tail.

A **Tail Break Length** value of zero will use the first profile value below background to terminate the tail, while a value of 10 will allow breaks of up to 10 microns (or other user-defined calibration units; this is a distance calibrated by your system settings) and then resume Tail measurement. Values for the Tail Break Length can lie in the range 0-50, the default value being 5.



4.11.1.4 Tail Threshold/1000

The default value is 10 or 1%.

The **Tail Threshold/1000** value is used to specify the intensity threshold of the cell for area measurements and the tail region for density measurements. Pixels with an Intensity value above the background by this percentage of the maximum head intensity are included in the measurements. This parameter has a range of 0-900 which corresponds to 0-90%, thereby giving fine control over the tail threshold value. Values can lie in the range 0-90%, the default value being 10 or 1%.

4.11.1.5 Maximum Head Diameter

The **Maximum Head Diameter** can be set here, provided that the Comet Options Operational settings check box is selected (**Section 4.11**). This feature is used for damaged Comet cells, where it is hard to distinguish between the head and tail. You can specify a maximum size for the Comet head which will be used to restrict the search for a normal head morphology. This option is only available when the **Max Head Diameter** is selected in Comet Options Operational settings. By default this value is NOT selected.

4.11.1.6 Background Height

The default value is 20.

The default is OFF.

The **Background Height** option is only available when **Auto background** (Section 4.13.2) is selected in the Operational section of the Comet Options dialogue. It sets the height of the region used in automatic background creation. This parameter is provided to allow you to adapt to the cell density and background level of a given slide. It will not affect results significantly, except where noise is high or background shows large intensity fluctuations. In such cases, the background should be adapted to be as tall as possible, consistent with maintaining vertical similarity throughout the background, and consistent with the background in the cell region.

Values can be in the range 10-50, with the default being 20.



4.11.1.7 Saturation Limit

The default value is ON.

The **Saturation Limit** option is only available when **Stop on Saturation (Section 4.13.4)** is selected in the Operational section of the Comet Options. It determines the number of saturated pixels, specified by the **Saturation Maximum** parameter (**Section 4.11.1.8**), that will be accepted in the analysis of a comet profile. Saturation of the CCD camera can occur at either the black level (zero grey level, 0) or the white level (full scale grey level, 254 for 8-bit data and 4095 for 12-bit data and 65535 for 16-bit data) and results in an over- or under-estimation of damage. If the number of saturated pixels exceeds this setting, the analysis will terminate.

If **Stop on Saturation** is not selected, the user can override Saturation detection and score the data anyway, but must be aware of the resultant errors which may arise in the analysis. We recommend the **Saturation limit** be set to a low value, such as 25.

4.11.1.8 Saturation Maximum

The default value is ON.

The **Saturation Maximum** option is only available when **Stop on Saturation (Section 4.13.4)** is selected in the Operational section of the Comet Options.

For an 8-bit image: min grey level = 0, max grey level = 255 (i.e. 2^8 -1)

For a 12-bit image: min grey level= 0, max grey level = 4095 (i.e. 2^{12} -1), etc.

A pixel is saturated when: $g \le minimum$ grey level or $g \ge maximum$ grey level.

Komet is capable of loading non-standard rasterised Image types. The **Saturation Maximum** parameter is available so that the user can input the correct maximum grey level value for such an image.



4.12 Comet Options: Experimental

Experimental Cells per Slide	75	Replicate Slides	2	
Samples	Ø	🔲 Retain Images in	Memory	
Use Automatic I		ies h.browne\Analysis\		
Use Database V (Automatic	Data Directo	ry used)		

Figure 101. The Comet Options Experimental Dialogue Box

4.12.1 Cells per Slide

Komet[®]

The default value is 25.

The **Cells Per Slide** value is used to prompt the user when to change slides and so forms part of the experiment protocol settings. When the system is in Experiment Mode, selected by clicking the **Experiment** button on the Comet Toolbar, the user will be guided through the Slides and Replicates. In most studies, replicate data will be gathered from multiple slides. After scoring the specified number of cells you will be prompted to change slides. Recommended values are in the range 1-100, with the default value being 25. Note that OECD guidelines recommend scoring 75 cells per Slide.

4.12.2 Replicate Slides

The default value is 2.

The **Replicate Slides** value specifies the number of Slides to include per Sample. It is used when Experiment mode is active, selected by clicking the **Experiment** button on the Comet Toolbar, to guide the user through the data capture phase of an experiment. Recommended values are in the range 1-20, with the default value being 2.

4.12.3 Samples

The default value is 2.

The number of **Samples** in a given experiment is designated here. At the end of the data collection phase of each sample, the data for that sample is saved to the desired file. If **Use Automatic Data Directories** is selected, the data is saved automatically (**Section 4.12.5**). Otherwise the user is prompted to input a filename and directory to which to save the data. This data will be useful for Dose Response curves and other comparative analysis. The default value is 2. The typical in-vivo safety testing sample number would be 30, corresponding to 6 treatment or exposure groups and 5 units per group. If multiple organs or cell types are to be tested, then this number of samples would be measured per organ. In bio-monitoring studies or cancer research there may be many other experimental designs and sample cohorts. Komet provides the flexibility to deal with those cases.

Komet[®]



After the final replicate slide in a sample is complete, if **Use Automatic Data Directories** (Section **4.12.5**) has been selected, the data for the sample is saved automatically. Otherwise a standard **Save Data File** dialogue window is shown at this point (Figure 102).

Save Data Fi	ile		2 🛛
Save in:	Simulated camera		* III •
LastSession			
File name:	*XLD		Save
Save as type:	Data Files ("XLD)	•	Cancel

Figure 102. The Save Data File Dialogue Box

4.12.4 Retain Images in Memory

The default is OFF.

The default is

The **Retain Images in Memory** option can be toggled on and off by a mouse click. When on, the check box is labelled with a tick.

When this option is selected, acquired images are stored in memory. You can view their properties and open them in the Image Window from the **Image List** menu (**Section 4.2.5**), which is only visible if this option is selected.

By default, Retain Images in Memory is NOT selected.

4.12.5 Use Automatic Data Directories

If **Use Automatic Data Directories** is selected, on completing each sample during an experiment the data is automatically saved to a default Data Path directory and filename:

C:\Andor\Komet CAMERAData\UserID\Date\ExperimentID\SampleID.XLD where

UserID is the User Name specified on starting Komet, **Date** is in the format year-month (e.g. 2014-07), **ExperimentID** is the ID given by the user to the experiment, after starting the Experiment, and **SampleID** is the ID given to the Sample by the user.

Komet®



The default directory can be changed by selecting the **Auto Save Path** via the **Browse** button on the Comet Options interface or the Set Save Path option on the Komet Control Window File Menu (**Section 4.5.1**). Experimental Data will then be saved to the following destination:

Auto Save Path UserID\Date\ExperimentID\SampleID.XLD, where **Auto Save Path** is the destination path selected.

If **Use Automatic Data Directories** is NOT selected, a standard **Save Data File** dialogue window will be opened, in which you can specify a directory and filename to which to save your data. The data is saved as a tab delimited *.XLD file, which can be read in Microsoft Excel.

By default, Use Automatic Data Directories is selected.

4.12.6 Use Database

The default value is OFF.

When **Use Database** is selected, a database is generated when conducting a Comet Experiment. The database will contain ALL the Comet data acquired during the experiment and will also store thumbnails of all the cells scored.

The location of the database is set in the Comet Options Experimental section.

If **Use Automatic Data Directories** is selected, then the database will use the same location, otherwise you can browse to select the folder in which to save the data.

By default **Use Database** is NOT selected.

4.13 Comet Options: Operational

Figure 103. The Comet Options Operational Settings Dialogue Box

4.13.1 Fluorescence

Andor's Comet analysis software is designed to operate with brightfield as well as fluorescence images. The **Fluorescence** option is selected by default, since most researchers operate the SCGE assay with a fluorescence stain such as Propidium iodide. At least one study is under way to investigate use of a dual colour label, Acridine Orange, to differentiate between single- and doublestranded DNA. Imaging will be available to quantify the relative fractions of each. Silver staining has also been used for comet visualization and tail length estimation, though to our knowledge, not for densitometric quantification.

4.13.2 Auto Background

When the Cell Region of Interest (ROI) is specified, if the **Auto background** option is selected, the system will automatically identify a background region immediately above or below the cell region. Thus, the user need only specify the cell region and ensure that cells presented for analysis have a clear background region around them.

4.13.3 AutoExpress on Store

The **AutoExpress on Store** option ensures that the system automatically performs a "Comet Express" analysis on the cell within the Region of Interest (ROI) when the **Store** button is pressed. This is equivalent to clicking the **Express** button on the Comet Toolbar (**Section 4.2.6**). Thus in the normal course of an experiment when the Cell region has been specified, the user presses the **Live** button on the Comet Toolbar, locates the next cell in the ROI, clicks on **Store** and the analysis is completed automatically. The user then clicks on **Live**, locates the next cell and clicks on **Store** again. This reduces operator fatigue and speeds measurements.

4.13.4 Stop on Saturation

As mentioned in the description of the **Saturation Limit** (**Section 4.11.1.7**), when image saturation is detected in the cell or background region, then over or under-estimation of DNA damage can result. The Saturation Limit is therefore used to define the number of saturated pixels which can be

127



The default is ON.

The default is OFF.

The default is ON.

The default is ON.

Komet®



accepted before terminating an analysis. Termination will only occur if the **Stop on Saturation** check box is checked and the **Saturation Limit** is exceeded. We recommend that the **Saturation Limit** be set to a low value. When **Stop on Saturation** is selected, the **Saturation Limit** and **Saturation Maximum (Section 4.11.1.8)** selection parameters will be available in the Processing section of the Comet Options dialogue.

4.13.5 Floating Live Button

The default is ON.

When selected, a floating **Live** button will appear on the Image Window close to the mouse cursor after each analysis is complete, to minimize the number of mouse movements in the experimental procedure. This reduces operator fatigue and speeds the scoring process.

4.13.6 Head Ellipse On

The default is OFF.

The **Head Ellipse On** option creates an oval around the head region of the cell being scored, after the image has been stored and analysed. This does not affect the way the cell is analysed, but rather provides a visual for the user, to show what the system is designating as the head region of the cell.



4.13.7 Maximum Head Diameter

The default is OFF.

Selecting **Max Head Diameter** will display a dialogue box in the Processing section of the Comet Options (Section **4.11.1.5**) where you can enter the required size. This feature is used for damaged Comet cells, where it is hard to distinguish between the head and tail. You can specify a maximum size for the Comet head and this will be used to restrict the search for a normal head morphology. Care should be taken with this parameter if used since it can impact measurements especially in heterogeneous specimens.

4.13.8 Head Symmetry

The default is ON.

In previous versions of Komet, **Head Symmetry** has been assumed to facilitate the edge of the comet head in the presence of a continuous distribution which overlaps the comet tail (quite common). Although this works well in most cases, it can cause problems in cases where controls show DNA distribution skew towards the anode (right hand side of the image), or with cells of unusual morphology. A number of specific examples where this assumption stops viable scoring have been brought to our attention. For this reason the Head Symmetry option has been created. In cases of unusual morphology the user may wish to disable this option.

When **Head Symmetry** is selected, Komet will continue with the analysis as in previous versions. When **Head Symmetry** is NOT selected, Komet takes the head extent as defined by the fitting algorithm or the interactive settings, but uses the intensity of the profile it finds, without creating a mirror image of the cathode, to centre the profile. This has been tested on a wide range of cell morphologies and found to create acceptable analysis in both automatic and interactive analysis (**Section 4.13.3**).

4.13.9 Tail From Centre of Cell

The default is OFF.

When the comet tail is measured, it can be computed either from the edge of the extrapolated cell profile, or from the cell's centre. The **Tail from Centre of Cell** option allows the results to be expressed in either form. Recent work has shown that measurements from the centre are good for low dose sensitivity, while those from the edge provide excellent discrimination against background effects, where significant damage is present.

Revised 13 Apr 2016

4.13.10 Interactive Tail

When Interactive Tail is selected, the user is prompted to draw the extent of the comet tail before each comet analysis will proceed. This has the same functionality as the **Interactive** button in the Comet Toolbar (**Section 4.3.2**).

4.13.11 Floating Open Button

When selected, a floating **Open** button will appear on the image window close to the mouse pointer after each analysis is complete. This will greatly reduce the amount of mouse movement required between each analysis and produce a more ergonomic experimental process.

With the **Floating Open Button** option selected, once in Experiment Mode you will be prompted to select the folder at the start of each sample which contains all your images for that sample. On pressing the **Open** button, the next image in the folder will automatically open.

4.13.12 Flip On Open

The **Flip On Open** option is used to adjust the positioning of stored images. The Komet system is set up to analyse comets in which the head is on the left and the tail is on the right. Therefore, if images have been stored in which the comet tail has migrated towards the left, then select Flip on Open. The images will be flipped so that the tail is now on the right and in the correct position for Komet analysis.

4.13.13 Highlight Scored Cells

The **Highlight Scored Cells** option is used to keep track of the cells scored on a single frame. When using the live option the trail becomes active once the store or interactive button is clicked. Going live again will clear the list. When using Stored Image mode the list will be cleared each time an image is loaded. This feature can be useful where the user has acquired images from a large format camera and there are multiple cells to be scored within a single frame.





The default is OFF.

The default is OFF.

The default is OFF.

The default is ON.



4.14 Komet Option Switches Tab

The Komet **Option Switches** tab provides Comet Options that can be changed during an experiment (Figure 104).

Protocol Options	
Images	
Retain Images in M	emory
Operational	
🔽 Auto background	🔲 Interactive Tail
AutoExpress on Store	e 🔲 Floating Open Button
Floating Live Button	🗖 Flip on Open
🔽 Head Ellipse On	Highlight Scored Cells
Calibration	
Apply selected calib	ration on Open Image
Calibration: iXon_10	0Xoi
-	

Figure 104. The Komet Option Switches Tab

When you set up a Comet Protocol, changing most of the parameters during an experiment will invalidate your data. For this reason, you are not allowed to edit certain parameters while in Experiment mode. However, there are several parameters which control the Experimental flow without invalidating the calculated data and therefore can be toggled on and off during an Experiment. The default values for these parameters are taken from the current Comet Protocol. Adjusting the parameters on the Option Switches tab will change the Experimental flow accordingly, although the configuration file will remain unchanged. The default parameters from the Protocol can be reloaded by pressing the **Reset to Protocol settings** button (Figure 104).

.

Komet®



SECTION 5 Concepts in Komet GLP A Reference Guide

5.1 Introduction

Komet GLP (KGLP) has been developed to fulfil the increasingly important requirements for electronic data capture and management as outlined in the FDA 21 CFR part 11 guidelines. In order to deliver to the guidelines we have interpreted the documentation and worked with laboratories involved with GLP comet studies and evaluated several beta versions in real scoring environments.

This product release is a result of painstaking efforts to make a solution, which is not only rigorous in implementation and robust in performance, but also convenient and effective in use.

In the design of KGLP, we have considered aspects of its use in a managed IT environment as follows:

- System Administration and Integration into existing IT systems.
- Personnel Access management and security.
- Hardware configuration.
- Systems and Software QC.
- Study Manager tools for Study Creation and Management.
- User Interface and Data Management.
- Distribution and Presentation.

In order to obtain the best from the software and systems supplied with KGLP we recommend that you use the information below as a reference to the product as well as reading the Komet User Guide for details on specific Settings.

Training in the use of KGLP can be provided by Andor personnel, but the following will provide a reference, which we recommend you read and use to prompt your memory if an infrequent user. Many of the concepts introduced here will be familiar to those involved with GLP studies, but some of the terminology is particular in its use to KGLP and we recommend your detailed review of the following sections of this User Guide.



5.1.1 Managing Validated Komet GLP workstations

For use in GLP (Good Laboratory Practice) contexts, software must be validated in each individual laboratory and it is done in combination with standard operating procedures in the laboratory. According to FDA 21CFR part 11 regulations, Komet7-GLP can be considered an 'off-the-shelf' software for comet assay analysis. Proceeding with local validation will be significantly simplified if this designation is used.

Further, because all changes in the computer system (including the Windows operating system) may impact performance and require repeat validation, we recommend that the Komet7-GLP workstation is operated on a computer in which automatic updates of Windows are disabled. We also recommend that workstations are not connected to the internet due to the risk of infection by malware.

However, systems can be run in a managed IT environment if this is the preference of the local management. The following explains use of Komet GLP in both contexts.

Data can be readily shared from the machine using CDR or DVD disks, USB flash memory "sticks" or hard drives. Komet and Database Viewer have built-in CD/DVD recording capabilities to make this process straightforward. If USB memory sticks or hard drives are used then it is very important to ensure the machine has adequate virus protection software



5.2 Continuity of Core Algorithms with QC Checking

Komet GLP (KGLP) uses the same field-proven analysis algorithms which have formed the basis of Komet software, first established in 1991 and improved incrementally with various versions since then. Care has been taken at each step to ensure backward compatibility and consistency to allow for comparison of historical and current datasets. These algorithms have been tested with synthetic images of known intensity distributions to prove their correctness in the absence of noise. In real situations where electronic noise is present they have been shown to deliver mean values which reflect the correct score according to the Central Limit Theorem, widely used in statistical analysis.

The algorithms have been encapsulated into a library module, which is used by KGLP in both Scoring and QC modes. This structure ensures that exactly the same piece of code is QC tested and used for routine scoring, simplifying the validation of software function in our factory and in the customer laboratory.

5.3 Other Services

The experience we have gained over more than 20 years in the use and development of Komet as well as our other products, means that we can offer other services to help with "21 CFR part 11-compliance", data collection, analysis, management and IT systems. We have also developed an extensive network of collaborators, which allows us to recommend or offer other further services to clients involved in comet assay and Genetic Toxicology studies. Please enquire for details.

5.4 Core Concepts and Terminology

5.4.1 Two-level Logon and Multi-mode Access

In line with GLP and 21 CFR part 11 recommendations, we have created dual level password protection for the software. The first level relies on Windows user logon, which allows the host IT department to manage system wide access and resources. The second level is managed by KGLP itself and provides 3 modes of access depending on the user rights as determined by the company or institution and may or may not involve the local IT department. Minimal experience of Windows Administration is assumed or required.

These three user modes are:

• Local Machine Administrator (LMA):

The LMA has ultimate control of software, hardware, QC and other User access to the system.

• Study Manager (SM):

The SM can specify, create and manage Studies, which contain Datasets and Structures, Scoring Protocols and Scorer lists.



• Scorers:

Scorers can access Studies in which they are authorized to score specimens or samples.

For more details on Users, Windows access levels and User setup please refer to Section 2.1.

5.4.2 Studies, Datasets, Scorers and Protocols

A Study Manager (SM) has the authority to Create Studies. A Study in KGLP is a structure, which contains all the information required for gathering and reporting data from a Comet Assay study.

elect 9	Study
	o Administrator. You are a Study Director and an inistrator on this system.
Plea	se select the Study for this session:
٩U	se Existing (please select from list):
	CO-55654-040903
	reate New (please enter a name): Enter new study name here
C P	eview Closed Studies
C S	ystem Setup
	OK Cancel

Figure 105. Existing Studies are selected from a list at start up.

A Study contains:

- An Identifier or Name, which once created, appears in a list in the KGLP Startup dialog as shown in Figure 105.
- An access password.
- A list of Scorers, who are registered users of the Windows PC being used to score the Study. They can be Local or Domain users and are selected by the SM.
- The information for KGLP to create one or more Datasets, consists of a name (ID), password, sample type, number of samples, number of replicate slides per sample and number of cells to be scored per slide.
- When complete the Datasets contain images and data from every cell scored in the study, cells scored and "deleted" (excluded from analysis but still recorded) as well as scoring and system audit trails.

Datasets can be viewed and distributed with Andor Database Viewer (DBV) to allow Review, QC and or Training.



SECTION 6 Administrator Functions

6.1 Creating Users in KGLP – Windows Access Issues

As stated previously in Section 5, in line with 21 CFR part 11 recommendations, we have created dual level password protection for KGLP. The first level relies on Windows user logon, which allows the host IT department to manage system wide access and resources. This means that the LMA has to have some information about user accounts under the Microsoft Windows networking systems. Thus ultimate control of users and their access rights can be controlled by either the local system administrator or the facility IT department.

Three user modes exist:

• Local Machine Administrator (LMA):

The LMA has ultimate control of software, hardware, QC and other User access to the system. The local machine means the PC(s) on which KGLP runs. See Important Notes below.

• Study Manager (SM):

The SM can specify, create and manage Studies, which contain Datasets and Structures, Scoring Protocols and Scorer lists. There can be many Study Managers. See Important Notes below.

• Scorers:

Scorers can access Studies in which they are authorized to score specimens or samples. Results from scoring are stored in the Datasets created by the Study Manager.

There can be multiple LMA users. LMA users should be selected from individuals who understand how to setup the KGLP system and its associated hardware.

LMA, as implied in the name, must have Administrator rights on the local computer, but does *not* need Domain Administrator rights.

Important Notes relating to Windows and Domain Users:

- LMAs are the only users who can Create, Change or Remove Study Manager users in KGLP. SM Users are a class of Power User on the local computer created by KGLP, but may have different rights on the various managed Domains in the company networks. The names and rights of Power Users may differ in different versions of Windows, e.g. NT4, 2000, XP, 7 and 8.
- A group named Administrator has to be created. This group has to be named Administrator no matter what the local language is because the software will only recognize that particular spelling as having administrator functions.
- Administrator can create local Study Managers and Scorers when logged on to the local machine. These uses do not need to have Domain rights.



- In order to access Domain User Lists (e.g. to Create Study Manager users), each LMA needs to have a domain logon with the same user name and password as used on the local machine logon. Study Manager users must be already setup as local or domain users. Similarly, SM users can "Create Scorer" users. Therefore to access Domain User Lists each SM needs to have a domain logon with the same user name and password as used on the local machine, though Domain access level may be different.
- Make sure that all Users who Create Datasets i.e. Scorers and Study Managers, have rights to Access Folders where data will be stored.

6.2 System Configuration by the Local Machine Administrator (LMA)

Note: First run the Andor Configuration Manager...

Before KGLP can be used with a camera or other hardware the LMA first must run the Andor Configuration Manager, which provides a hardware profile (configuration name) for later use in KGLP and other Andor software products. See Appendix 2.

For GLP Operation, Multi-level Access must be enabled in the Configuration Manager Global Settings Page. We also recommend using the Virtual Image Disk Setting. In Komet Option Switches deselect Images in Memory.

6.3 Running KGLP in System Setup Mode - Local Machine Administrator (LMA)

The first time you run KGLP you must run in System Setup Mode, where the following settings will typically be configured or created:

- Hardware (e.g. camera) settings will be confirmed and configured other users have access only to limited camera settings.
- Study Managers can be Created, Removed and Changed.
- Scorers can be Removed from All Studies.
- QC Frequency is set.
- Password change frequency can be set for Studies and Datasets.
- System Timeout is set.
- GLP mode can be switched on or off. In non GLP mode KGLP reverts to standard Komet operation (no audit trails and different data entry requirement).

Komet®



Select Study	
Hello mbrowne. You are a Study Direct Administrator on this system. Please select the Study for this session	
C Use Existing (please select from lis	Password Entry
Enter new study name here	Please enter the password for Setup.
C Create New (please enter a name)	Landard Control of Comp.
Enter new study name here	posses
C Review Closed Studies	
System Setup	
	OK Cancel
Ok Cand	

Figure 106. . The Local Machine Administrator starting dialog for KGLP. The very first time KGLP is accessed in Setup Mode the LMA must define the Setup Password.

When the LMA logs onto Windows and starts KGLP, the starting dialog shows four options as shown in Figure 107. The LMA is the only user who can access the software in System Setup Mode. This option will not be offered to other users.

When the System Setup mode is selected (Figure 106), KGLP requires a password for access to LMA functions. The LMA sets this password the very first time KGLP runs the software in a Password dialog requiring password entry and confirmation. Thereafter the Setup Mode password can be managed/changed from the Security menu – see section 2.4.6.

Be sure to make a note of the Administrator Password for future reference!

KGLP then starts up and shows the Configuration selection dialog, which is shown in standard Komet operation. Here the Administrator selects a hardware configuration, which will be automatically selected for other users thereafter.

KGLP can be operated with a range of analogue and digital cameras and, as outlined in Chapter 1, the camera is specified in the Configuration which is setup by the LMA.



Please select the Configuration for this session: Virtual Camera Q Imaging Orca	
	•
Orca	^
Andor Zyla Video CCD	1 10 10
	~

Figure 107. Selection of the Hardware Configuration by the Local Machine Administrator. Once the camera is selected KGLP will start up with two Windows as described in detail in the Komet User Guide. These are referred to as the Control and Image Windows respectively, as shown in Figure 108.

Exercis Insurance Exercise 6.0.02P The Security OC Weekadd Itals without Pug-liss Hep-	Later former factors
Construction of connections waters inducts into the integration of connection of the integration of the inte	
System Setup Mode	Sec. 2 Million
Image Calibration: X*0.825 : Y*0.825 Protocol	
Ernet	
Analyse Devices	
Andyss	
Electricity Purples 5,000 4,000 1,000 0,000 1,	
Note: Name Cell Area Head DNA Tail Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Head DNA Tail Tr. Tr. 1 Image: Cell Area Tail Tr. Tr. Tr. Tr.	
Lord Jave	3.2 1005 2015

Figure 108. Control (left) and Image Windows (right) in Komet GLP.

KGLP shows a blue message in the Control Window indicating the current operating mode, in this case "System Setup Mode". In the Current Context section at the top of the Control Window, the

Komet®



User Profile, Configuration name and Calibration settings of the Image displayed in the Image Window is also shown.

The LMA can now specify the System Setup as desired. The following dialog descriptions explain and illustrate the features and settings which can be controlled.

6.4 Security Menu

6.4.1 Personnel Access Control

Personnel Changes	Create Study Director	
Time Out Delay Study Password Change Frequency Study Password Reset Dataset Password Reset	Remove Scorer from All Studies Remove Study Director Change Study Director	
Change Setup Password • GLP Mode		

Figure 109. The Security Menus

Personnel changes can be made from the Security menu. SDs can be Created, Changed or Removed and Scorers can be Removed from All Studies. This is an important and powerful security tool when an employee leaves the company. The Study Manager group is a subset of Power Users on the system and can be specified on the local machine or on selected local Network Domains.

Note: In the previous versions of Komet GLP Study Managers were referred to as Study Directors. Please substitute Study Managers if we have not updated text and you find further references to Study Director.



Add
Remove User Selection Select User to add to Study Directors group
Local Computer Domain Update
Administrator Guest HonyO JoeSD rey UserR

Figure 110. Clicking on Add or Remove in the first dialog provides the interface for Local or Domain users to be selected and their KGLP Access Rights controlled.

When Create Study Manager is selected the dialog in Figure 111 is shown. This allows Users, who must already be registered on the system or selected domain, to be granted SM rights. If a SM is Removed from all Studies then KGLP will request a substitute SM to be provided from existing SDs so that the Study does not become orphaned and therefore inaccessible.

6.4.2 Timeout Delay provides a lockout when KGLP is left open

Please select the ne out and re-logon:	w delay for time
Current TimeOut	1 mins
New TimeOut	1 mins
ок	Cancel

Figure 111. Selecting the Timeout Delay from the Security Menu allows the Timeout Delay to be modified. After the set number of minutes, KGLP will switch to a screen saver, which requires the user to log into Windows with a password again.




6.5 Study Password Change Frequency

Administrator Functio	ns:
Please select the new change frequency:	v study password
Current Frequency:	30 days
New Frequency:	30 days
OK	Cancel

Figure 112. Selecting the Study Password Change Frequency allows the current frequency to be changed to a new frequency.

6.5.1 Study Password Reset

Select Study		
Select the Study you wish to Reset the passwo	ard for	
CO-0127-09445	•	
Study Director: MARK-LAPTOP\Administrato Number of Datasets: 7 LIVER	r	
LIVER_RS1 LIVER_RS2	Reset P	assword
3	<u> </u>	Are you sure you wish to reset the Study password for Study 'CO-0127-09445'?
OK Cancel		Yes No

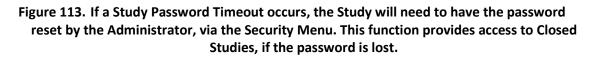

The next time the Study Manager opens the Study, KGLP will ask for a new password. If you are a Scorer, KGLP will ask you to contact your SM to update the password before scoring can proceed.



6.5.2 Dataset Password Reset

		Kinetic Imaging Komet 6.0 GLP
Select Study	Select Dataset	Bio Separity QC View-Budit Trails Woards Magitins Bolp Setup Waard Commet setup
Select the Study which contains the dataset.	Select the required Dataset	Dateset
Testing 123 Study Director D9×M9341\Jen Torelli Number of Datasets: 1 Basal001	Komet Dataset - Basal001	Based001 Cencel Sotup Dataset Current pasaword: Select pasaword:
	Selected Dataset Bosol001	Contirm password:
Ok Cancel	Select	NB: Database passwords have a maximum length of 14

Figure 114. A Dataset password can be Reset. Select Study, Dataset then enter the new password

6.5.3 Change Setup Password

Change Password	
Change password for	or Setup:
Old Password:	Autobala
New password:	Autodologi
Confirm New Passwo	rd: www.
ОК	Cancel

Figure 115. The System Setup

Access Password is changed by selecting from the Security Menu and then entering the required information in the dialog shown opposite.

6.5.4 Setting up Hardware

The Local Machine Administrator sets up the hardware devices for all users. Selecting the Devices Tab from the Control Window, shows the Device Setup dialog. In the example, shown in Figure 116,

.





the Andor dummy camera settings are shown. Each camera has its own interface and depending on the features, other controls will be revealed.

	Acquisition/Auxiliary Devices
Komet	Camera
Analyse Devices	Image Control
	Contract 160 Esposue (ms) Frames
	-Set Window 768 : 576
Device Setup	Advanced >>
	Kinetic Dummy Camera

Figure 116. The Administrator is the only user who can change the hardware settings in the camera or other devices controlled by a specific Komet GLP system. Other users will have access only to exposure time, brightness and contrast which are defined in the Protocol.

6.5.5 The QC Menu allows the Setup of QC Tests, Frequency and Reset after Lockouts

Setup QC Reset Lockout Run	
V Setup QC frequency	
Setup QC Procedure.	Setup QE frequency
Algorithm Validation Plug-In Stross Test Plug-In	Setup QC Procedure:
Please selected required QC plugins	☐ Algorithm Validation Plug-In ☐ Stress Test Plug-In
Report Directory : C\Documents and Settings\All Users\Application D	Please selected required OC plugins Frequency: 0 days Report Directory :
Backup Directory	C\Documents and Settings\All Users\Application Data\Kinetic Imaging\K
C\Documents and Settings\All Users\Application D	Backup Directory : C\Documents and Settings\All Users\Application Data\Kinetic Imaging\K

Figure 117. Selecting the QC Menu settings allows the Administrator to set the QC checks and the frequency of these QC checks.

Komet[®]



his is a system wide setting in which a QC check will be run every time KGLP starts and the frequency period has elapsed. The interface allows the Administrator to define where the QC report files will be logged. The QC reports will also be incorporated into the Study Datasets for the period during which the data was gathered. Thus, the Audit Trail will provide evidence that KGLP was functioning correctly during the duration of the data capture period.

The QC Test is suggested to be run by users with locally gathered and stored images. Results should be checked against standard values gathered during Validation tests. Komet facilitates manual test only, here the user should create a check list of their own choosing. The tests performed should validate that the system will create consistent images and measurements.

If a QC check fails three times in an execution then KGLP locks out all users until the Administrator Resets the Lockout. The system

QC should then be Run by the Administrator and any problems resolved. Further advice or service requirements can be addressed to Andor Imaging.



6.6 View Audit Trails Menu tracks Local Machine Administrator and System QC activity

1

QC Log Amended by W2000901FLVAdministrator Date 28/01/2004 ImmeDut Delag(min):1 Time:0ut Delag(min):1 Date 28/01/2004 Study Password Change Frequency(days) Lock Out:0 Date 28/01/2004 Amended by W2000901FLVAdministrator Date 28/01/2004 Date 28/01/2004 TimeOut Delag(min):1 TimeOut Delag(min):1 Study Password Change Frequency(days) In Use:0 Study Password Change Frequency(days) Date 28/01/2004 Frequency (days) In Use:0 Study Password Change Frequency(days) Ocketup report Directory: 16/05/2003 Amended by Amended by MARKLAPTOPVAdministrator Ocke: 23/07/2004 TimeOut Delag(min):1 Study Password Change Frequency(days) Ocke: 23/07/2004 TimeOut Delag(min):1 Number of DC:: 2 Appointm Validation PlugIn TimeOut Delag(min):1 In Use: 0 Study Password Change Frequency(days) Date: 23/07/2004 Time: 13/810/7 TimeOut Delag(min):1 In Use: 0 Study Password Change Frequency(days) Date: 23/07/2004 Time: 13/810/7 Date: 23/07/20	QC Setup	N Administrator Settings file		
FineDub Delay(min): 1 Study Pasaword Change Frequency(days): [current] Lock, Out, 0 Amended by MARK-LAPTOPVAdministrator Date: 29/07/2004 Time: 13/18/21 Date: 29/07/2004 Number of QCs: 2 Date: 29/07/2004 Number of QCs: 2 Date: 29/07/2004 In Use: 0 Time: 19/19/20 In Use: 0 GLP Mode: 1 Strate Test Plug-In In Use: Study Pasaword Change Frequency(days): Lock Out 0 Settings: Settings: Core Operation: Study Pasaword Change Frequency(days): 5 Operational Settings/All Users/Application Data/Kinetic Imaging/Komet/ Core Directory: C/Documents and Settings/All Users/Application Data/Kinetic Imaging/Komet/	QC Log	Amended by W2000901RLVAdministrator Date 28/01/2004 Time 11:11:57 GLP Mode 1 TimeDut Delag(min):1 Study Pasawood Change Frequency(days) Lock Out:0 Amended by W2000901RLVAdministrator Date 28/01/2004 Time 13:18:30 GLP Mode 1 TimeOut Delag(min):1 Study Parawood Change Frequency(days) Lock Out:0 Amended by MARK-LAPTOPVAdministrator Date 29/01/2004 Time 18:33:33 GLP Mode: 1 TimeOut Delag(min):1 Study Parawood Change Frequency(days) Lock Out:0 Amended by MARK-LAPTOPVAdministrator Date 29/01/2004 Time 18:53:43 GLP Mode: 1 TimeOut Delag(min):1 Study Parawood Change Frequency(days) Lock Out:0 Amended by MARK-LAPTOPVAdministrator Date 29/01/2004 Time 18:53:43 GLP Mode: 1 TimeOut Delag(min):1 Study Parawood Change Frequency(days) Lock Out:0 Interfield by MARK-LAPTOPVAdministrator Date 29/01/2004 Time 19:19:20 GLP Mode: 1 TimeOut Delag(min):1 Study Parawood Change Frequency(days) Lock Out:0	Note Date: 28/01/2004 Time: 11:10/43 Number of 0Cs: 2 Algorithm Validation Plughn in Use: In Use: 0 Frequency (days): 0 QC report Directory: MARK-LAPTOP-Vachimistrator QC backup report Directory: 16/05/2003 Amended by: MARK-LAPTOP-Vachimistrator QC backup report Directory: 16/05/2003 Amended by: MARK-LAPTOP-Vachimistrator Date: 29/01/2004 Time: 15/5/07 Number of QCs: 2 Algorithm Validation Plughn in Use: Stemss Test Plug/n in Use: D' backup report Directory: C-Documents and Settings/All Users/Application Data/Kine QC backup report Directory: C-Documents and Settings/All Users/Application Imaging/Komet/v (current) Amended by: MARK-LAPTOP-Vachimistrator Date: 29/01/2004 Time: 13/8/21 Number of QCs: 2 Algorithm Validation Plug/In In Use:	Data/Kinehc fic Imaging/Komel\

Figure 118. The View Audit Trails menu allows the LMA to i) Check the Administrator function activity ii) Check QC Setup activity and iii) Check QC tests and their reports.



6.7 Wizards Menu

System Calibration is performed by the Administrator for all Studies – New Objectives, Cameras or Microscopes can be configured for all users.



Figure 119. The Administrator calibrates the system and the SM selects the magnification at which the scoring is to be applied. The Calibration Manager provides a Wizard, which guides the Administrator or SM through the calibration process. Details are available in the Komet User Guide.

System Calibration must be completed prior to running the system for Study Creation and data collection.

Komet®

SECTION 7 Study Manager Functions

7.1 Studies, Datasets, Scorers and Protocols

A Study Manager (SM) has the authority to Create Studies. A Study in KGLP is a structure which contains all the information required for gathering and reporting data from a Comet Assay study. As listed below a Study must contain...

dministrator on thi		
hease select the 5	tudy for this session:	
Use Existing (ple	ease select from list):	
HX04-001		
HX04-001		
CH0x04-091		
Enter new stud	y name here	
Review Closed	Studies	
System Setup		

Figure 120. Existing Studies are selected from a list at start up.

- An Identifier or Name, which once created, appears in a list in the KGLP Startup dialog as shown in Figure 120. The Study list is filtered to make visible only those Studies for which the current user is authorized.
- An access password.
- A Scoring Protocol, which holds Comet Options, Calibration and Camera settings.
- A list of Scorers, who are registered users of the Windows PC being used to score the Study. They can be Local or Domain users and are selected by the SM.
- The information for KGLP to create a set of named and password protected Datasets, which includes sample type, number of samples, number of replicate slides per sample and number of cells to be scored per slide. When complete the Datasets contain images and data from every cell scored in the study.



7.2 Create New Study – Name, Password Entry – Use Template Option

Select Study		
Helio Administrator. You are a Study Direct Administrator on this system. Please select the Study for this session C. Use Existing (please select from list).	Select Study Do you want to use an existing file as a EX04-001	templote?
CO-0127-09445		Password Selection
Create New (please enter a name) Ho014-099 Review Closed Studies System Setup	Study Director WX20005DIRL(mbrow Number of Datasets: 2 001 002	Please create a password for Study HM04-039 Enter Password:
OK Cancel	Yes No	OK Cancel

Figure 121. Create New Study is Selected from the Start up Menu for Study Directors. The first dialog to follow Start up is the Template query as shown in Figure 121... if a previous Study has a similar structure then it can be used as to create a new Study. KGLP will ask the Study Manager to select an existing study to use as a template.

The same Scorers, Protocols and Datasets will be created. Any of these can be edited from the appropriate Manager as shown on the tabs in Figures 122 – 126. If the SM declines the Template then the Study details must be created from scratch.

The Password Setting dialog is then shown, which requires a duplicate password entry. Any alphanumeric format password is acceptable, provided more than four or more characters are used. After the password is accepted the Study is created and the SM is then able to create or edit Scorers,

Datasets and Scoring Protocols

Note: The SM can only use Studies created by him/herself as Templates.





Komet	Kinetic Imaging Komet 6.0 GLP
Scorers Datasets Scoring Protocols	Bit Right Weiwiguide Train Weiserds Plugstris Belo Setup Witard Scorers Study Scorer Manager Study Scorer Manager Study Scorer Manager
. Scorer - D9XM9341\Jen Torelli	Add Remove OK Cancel Scorers assigned to this Study: Current Scorer D9XM9341\Jen Torelli R. Scorer - D9xM9341\Jen Torelli

Figure 122. The SM, who creates a Study, is by default a Scorer. Selecting the Manage Scorers button shows the Study Scorer Manager, where Users can be Added or Removed from the Study.

At least one Scorer must be present. The SM is by default a Scorer in the given Study. To Add or Remove Scorers the User Selection dialog is shown and lists the Local Computer or specified Domain users (Figure 123) to be added or removed.

User Selection	
C Local Computer C Domain 2000-Serve	Current Scorer MARK-LAPTOP\HarryO
JoeSD ray UserR	E: Scorer - MARK-LAPTOP/Administrator E: Scorer - MARK-LAPTOP/Henvo - Full Name - Harry Olson
Full Name: Hany Olson Comment: Study Director Select Cancel	Last Logon to this Computer - Wed Jan 28 14:21:31 2004 Scorer - MARK-LAPTOP\JoeSD Scorer - MARK-LAPTOP\rey

Figure 123. Selecting a User, e.g. HarryO, from the list allows the User to be Added or Removed. Once selection is complete and the Study Scorer Manager dialog is closed, the Scorers Tab in the KGLP interface allows you to Review Scorer details, including recent activity.



7.3 Adding Datasets – Specify Name and Optional Folder

Elle Study QC View Audit Trails Wizards Plug-Ins Hel	p		
User: Jen Torelli (Study Director + Adm Study: Experiment 001 Image:	Kmetric (maging Komet 6.0 GLP Pile scath, gc, view facts the s. Witeries Setup Wand Comet setup Enter name for Dataset [Basal001]	Form Heb	
Please press Experiment to select a Dataset	Setup Dataset Select Detabase path.	Browse for Folder Select a Folder	থ× শ
Komet Scorers Datasets Scoring Protocols Anage Datasets	Date will be served to : \+K04-099(Basel001\	Guotes RECYCLED STSNCD Government Market RecycleD Government Market RecycleD Government Market RecycleR My Network Places My Network Places	
	Deta vill be served to UP04-029(Decal001),	Admin Files on Kinetinemal	

Figure 124. Selecting Manage Datasets shows the Comet Dataset Manager and the SM specifies a name for the Dataset, in this example Basal001.

The location to store the Dataset is specified using the Browse button and the Folder dialog shown above right.

Note: The optional Backup Copy of Dataset Option has now been removed for reasons of data integrity.

Important Note: Datasets Folders can be placed anywhere on the network, but SDs should be sure to locate them in Shared Folders, where ALL Scorers have access rights. Otherwise some Scorers will be unable to add to Datasets.

The other details of the Dataset, password and structure are added by selecting Next, as shown in Figures 125 and 126.

.



7.4 Adding Datasets - Password setup

Komet®

e Study <u>QC</u> View <u>A</u> udit Trails <u>W</u> oards etup Wizard	Plog-tris Help
Comet setup Dataset Basal001	1
< <pre><< Back Next >> </pre>	<u></u> ancel
Select password:	_
Confirm password:	

Figure 125. After selecting "Next" KGLP prompts the SM to assign a password to the dataset and confirm the password. Each Dataset can be given a different password, enabling the SM to assign and manage the Datasets among different Scorers. Make sure to take a note of the Passwords or use an agreed structure known to the whole team!

7.5 Adding Datasets- Dataset structure

Kinetic Imaging Komet 6J0 GLP	Kinetic Imaging Komet 6.0 GLP
File Study QC Merr Autit Trais Witerds Puppins Help	File Study QC Wen Acit Trais Weards Pupins Help
Setup Wizard	- Setup Wizard
Cornet setup Deteset	Comet Dataset
	Comet Dataset Manager
Besel001	
<< Back Next >>	Qancel New Edit Close
Describe Dataset	Dataset Summaries
	Current Dataset
Number of Samples. 30	Basal002
Slides per Sample	Komet Dataset - Basal001 S Komet Dataset - Basal002
	Database - C\\HX04-099\Basel002\Basel002\Basel002_R4W kdb Backup Database - C\Documents and Settings\AdministratorbD
Cells per Slide: 50	E Experimental Number of Samples - 30
Total Number of Slides = 60	• Number of Slides - 2
Sample Type: bladder	- 10 Cells per Slide - 50 - 10 Sample Type - bladder
Southe type. Introduct	

Figure 126. The Dataset structure is entered in the Wizard dialog by selecting 'Next".

Komet®



Number of Samples, Slides per Sample, and Cells per Slide are entered and Total Number of Slides for scoring is computed by KGLP. Enter Sample Type for the dataset and select "Next". The Comet Dataset Manager appears once again, where the structure of the Dataset(s) can be Reviewed using the '+' and '-', controls in the List and, if necessary, edited.

7.6 Review and/or Edit Established Datasets

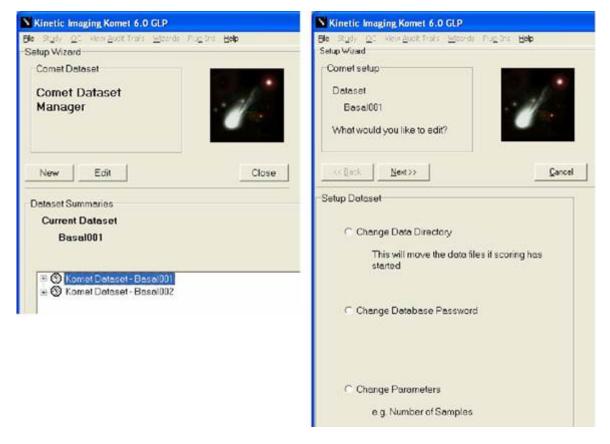


Figure 127. Once a Dataset has been added the Comet Dataset Manager allows Review and/or Editing of dataset details such as Change Data Directory, Change Database Password, and Change Parameters.

Edit the Dataset by selecting the Edit button or by a right mouse click on a dataset in the list. To Change the Data Directory (Folder), Database Password and/or Parameters follow the Wizard instructions and use the same dialog tools described in Figures 121 – 127.





Ele Study QC View Audit Trails Wittends F Setup Witzard	ing Ins Help
Comet setup	-
Doteset	· d
Basal001	Cancel
Setup Dataset Current Database path \\2000-SERVER\Server Filestore Select Database path:	NExperiment 001\Basal001
D.\	Browse
Data will be saved to : D:\\Experiment 001\Basal001\	

Figure 128. Datasets Folders can be placed anywhere on the network, but SDs should be sure to locate them in Shared Folders, where ALL Scorers have access rights. Otherwise some Scorers will be unable to add to Datasets.

When changing Dataset folder locations, make sure to update both Raw and Backup paths. If the Backup checkbox is not checked, then the Dataset will NOT be saved, even if previously checked and greyed out.

7.7 Adding New Scoring Protocols – Protocol Name, Calibration, Camera and Comet Options

omet		GLP	
Scorers Datasets Scoring	Protocols	Wizards	Plug-Ins He
			ation
Manage Protocols		Comet Protocol	
		Score	r Manager
F Komet Protocol - D	efault	Come	t Dataset

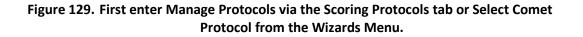








Figure 130. Then select New to create a new protocol or Import (Figure 131) to use a protocol from a previous study.

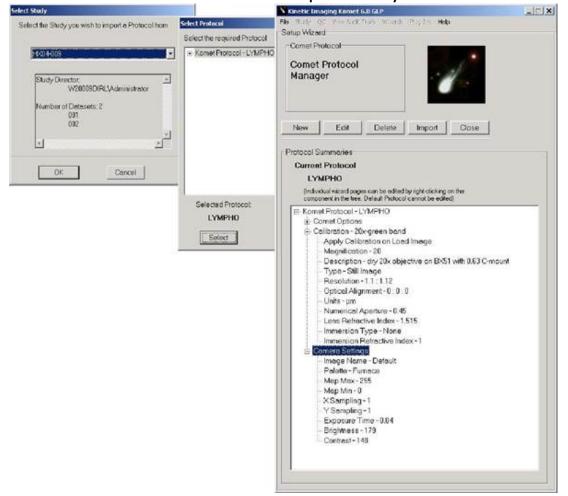


Figure 131. Selecting Import allows you to re-utilize a Protocol from a previous Study. Select the Study from the list shown and then Select the Protocol you wish to use from those listed within the Study. Once selected, the Protocol will appear in the Protocol Manager list. Right mouse clicking on the items provide Edit options.

Komet®



Select Calibration Stored Images Apply selected colibration on Open Image Select Calibration © Use existing: 20x Olympus © Et Olimaging 20X green band 20x Olympus © Create new General Select Calibration on Open Image Select Calibration on Open Image Camera Select Calibration (Camera Select Calibration (Camera	Solup Waad Comet setup Enter Protocol name: SVER Gold LIVE LIVE Select Comera Settings Image Options Image Name: Default Poloto: Funaco Select Comera Settings Image Options Image Name: Default Poloto: Funaco Set o Factory Defaults Ado Exposue Taget 2 Saturation C 100 C 50 C me: 40	Kinetic Imaging Komet 6.0 GLP		Kinetic Imaging Kome	1 6.0 GLP
Enter Protocol name: SYBR Gold Enter Protocol name: SYBR Gold LIVE LIVE Cancel Cancel Cancel Cancel Cancel Cancel Cancel Cancel Can	Enter Protocol name: SYBR Gold LIVE LIVE Cancel		Help		Trais Wizards Plug-Ins Help
VESC Net>> Select Calibration Cancel Stored Images < Apply selected calibration on Open Image Select Calibration © Use existing: 20x Olympus 20x Olympus Image Name: Centers Sinning Gain and Offset Control Gain 128 Offset Image Name: Cancel Select Calibration Concel Fundace Select Calibration Image Name: Original 20x Original 20x green band Image Name: Cancel Select Control Gain 128 Offset Image Name: Can	Cancel LIVE Cancel << Back Next >> Select Comera Settings Image Options Image Options Image Name Default Paloto: Fumace Paloto: Fumace • Binning Gain and Ottest Control Gain: 123 Ottest 0 Canera Binning Gain: 123 Ottest • Paloto: Fumace • • Canera Binning Gain: 123 Ottest • Canera Fundo • • • Otto foid foid • • • Otto foid foid foid • • Otto foid foid foid foid • Image Statistion foid foid foid foid foid				
Select Calibration Stored Images Apply selected calibration on Open Image Select Calibration Use existing: 20x Olympus Category C	Binning Select Comero Settings Image Name. Default Pelotic: Funace Senband Camesa Set brack Gain and Ottest Control Gain 128 Set brack Set to Factory Default Auto Exponse Exposse Time Target % Saturation Cimes Comesa Cimesa	SYBR Gold	· Mar		1
Stored Images Image Options Image Options Image Name: Select Calibration Image Name: Outse existing: 20x Olympus 20x Olympus Image Name: Camera Siming Gain and Ottest Control Gain 123 Ottest 0 20x Olympus Image Name: Create new Set to Factory Defaults Auto Exposuse Exposuse Time Target & Satuation Image 20 Image Options Image Name: Default Palots: Funace Image Name: Outset 0 Image Name: Outset 0 Image Name: Set to Factory Defaults Image Name: Image Name: Exposuse Time Target & Satuation Image Name: Image Options Image Name: Image Name: Exposuse Time Image Name: Image Name:	Image Options Image Name: Default Pelotto: Furnace Camera	Next Next >>	Cancel	<< Back Next>>	Cancel
Image Name: Default Paloto: Funace Select Calibration Image Name: Clock Olympus Image Name: 20x Olympus Image Name: 20x Olympus Image Name: Camera Select Control Camera Select Control 20x Olympus Image Name: Control Image Name: Default Paloto: Funace Image Name: Camera Select Control Gain 123 Othet: Image Name: Create new Set to Factory Defaults Auto Exposue Image X Satuation C 100 50 C 75 C 75	Ibrotion on Open Image Image Name. Default Paloto: Fumace Canera	Select Calibration			
Apply selected calibration on Open Image Select Calibration © Use existing: 20x Olympus 20x Olympus © EdQImaging 20X green band 20x Olympus © Create new	Peloto: Fumace Peloto: Fumace Camera Camer	Stored Images			
Imaging 20x Olympus Imaging 20x Green bond C EtQUmaging 20x green bond Imaging 20x Green bond C Create new Imaging 20x Green bond C Create new Imaging 20x Green bond Imaging 20x Olympus Imaging 20x Green bond	Sinning Gain and Offset Control Gain 128 Other 1 Set to Factory Defaults	C Apply selected calibration on Open	Image		a de la companya de l
20x Olympus Image: Control state of the contrel state of the contrel state of the control state of the control	Sinning Gain and Offset Control Gain 128 Other 128 Set to Factory Defaults	Select Calibration			
C EdQImaging 20X green band Binning Gain and Offset Control 20x Olympus C 1x1 3x3 Gain 123 Offset 0 20x Olympus C Create new C 2x2 4x4 Set to Factory Defaults Set to Factory Defaults Auto Exposuse Target % Satuation C ms 0 1 1 1 C 100 F 50 C 25 C 25 C 25 C 25 C 25 C 25	een band 	Ose existing:		Camera	
C Ed Qimaging 20X green band 20X Olympus C Create new Auto Exposuse Target % Saturation C 100 C 50 C 75 C 25 C 25	een band C 1x1 C 3x3 Gain 128 Offset 0 C 1x1 C 3x3 C 2x2 C 4x4 Set to Factory Defaults Auto Exposuse Exposuse Time Target % Saturation C ms s 0 1 - 1 C 100 C 50 C me 40 - 1 C 75 C 25 C us 40 - 1		-		C 1 - 1 - 1 - 1 - 1 - 1 - 1 - 1 - 1 - 1
20x Olympus Image: Constant and Constant	Image: Set in Factory Defaults Image: Set in Factory Defaults Auto Exposuse Target % Saturation Image: Set in Factory Defaults	C Ed Olmaging 20X green band			
Auto Exposuse Exposuse Time Target % Saturation C m: s 0 1 1 C 100 6 50 c ms -40 1	Target % Saturation C m s 0 1 1 C 100 6 50 G ms 40 1 C 75 C 25 C us 40 1	20x Olympus	9		
C 100 @ 50 @ me 40 +	C 100 € 50 C 75 C 25 C us40 ± C 40 ±	C Create new		-Aulo Exposure	Exposure Time
	C75 C25 C ₩ 40 ±			Target % Saturation	C m:s_0 _1 ±
C 75 C 25	cu ■ Cu _4 0 🗄				@ ms 40 +
	Set Actual: IbLActuaEx			and the second	C us 40 -
Set Actual: IbLActuaExi				Set	Actual: IbLActualEx

Figure 132. Enter the Protocol name and choose a Calibration option, which should match the microscope objective you plan to use for Scoring the Study. If no Calibrations exist then you can create one using the Calibration Wizard described in detail in the Komet User Guide. After Calibration, the Next button will show the Camera Setup panel. Adjust the settings to obtain good quality images, referring to expert advice if needed.

KGLP supports a wide range of cameras and camera setup is important for obtaining good image quality, which is central to achieving precision data from comet analysis. Image quality is also affected strongly by microscope configuration and fluorescence lamp stability and alignment. See the Komet User Guide for more details. A QC tool is available for helping to maintain consistency in these areas.

To help with Camera settings, you can place a specimen under the scope and image it by hitting the Live button. Then you can measure the effects of changing camera settings. Using factory default settings for Gain and Offset (or Brightness and Contrast) usually works well. If you have a digital camera, then using the Auto-exposure tool will help establish the correct exposure setting for good quality imaging.





Kinetic Imaging Komet 6.	0 GLP	
He Study QC View Audit Trah Setup Wizard	Wizards Plug-Ins H	Чр
Comet setup Enter Protocol name:		
SYBR Gold	_	ali
To test settings click Righ Mouse Button in Image W Then see Key Results fiel	indow.	1
<< Back Next >>		Cancel
Comet Options Processing		
Head Threshold% 10		
Smoothing Value 1	Background Height	20
Tail Break Length (um) 5	Saturation Limit	25
Tail Threshold/1000 10	Saturation Maximum	4095
Experimental		
Show Database Gallery at en	d of Slide:	
C Never	Store Full	Image
Ask each time	F Require E	quipment IDs
Scoring mode:		
C File		
IVe		
Operational		
Fluorescence	Max Head Diame	ter
Auto background	Head Symmetry	
AutoExpress on Store	Tail from Centre of	f Cell
Stop on Saturation	Interactive Tail	
	Floating Live/Op	en
🔽 Head Ellipse On	Flip on Open	

Figure 133. The Next button will take the SM to the Comet Options panel, where analysis settings are entered. To help establish the Settings, you can perform analysis on real images, by placing the Comet Region of interest (ROI) over a cell and right mouse click. KGLP will score the cell with the current settings and then you can adjust them until the indicated Head/Tail callipers correspond to the visual assessment.

Features special to KGLP options include:

- Ask Scorer if a Comet Gallery is shown at the end of each slide scored.
- Store Full Image field for the comets scored.
- Require Equipment IDs during scoring.

Show Database Gallery at end of Slide, allows either Never or Ask to be chosen. Store Full Image saves the whole image field instead of only the scored cell regions, providing review of overall specimen quality and variability as a further check on the data capture process. Require Equipment IDs allow the equipment used for scoring to be recorded for each scoring session.





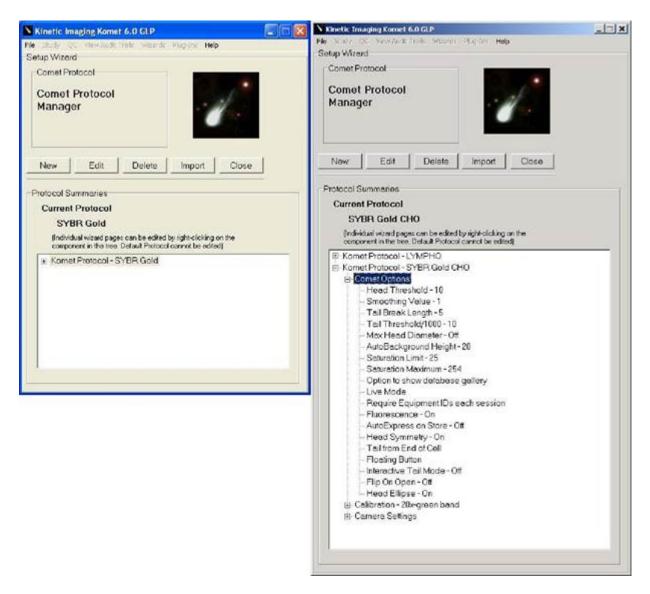


Figure 134. Once Calibration, Camera and Comet Options are setup the Comet Protocol Manager will now display the Komet Protocol Name rather than the Komet Default.

Clicking on the + symbol will open up the details of the Protocol Tree, similar to Windows Explorer function. Right click on Calibration, Camera Settings or Comet Options will provide an Edit menu option to edit the individual component of the

Protocol. The information about this Protocol will be stored in the Dataset System Audit Trail tab, once scoring the Dataset is complete. The information can be viewed and printed in the Andor Database Viewer.

.



7.8 Use Existing Study - Study Manager

Select Study	
Hello Jen Torelli. You are a Study Director and an Administrator on this system. Please select the Study for this session:	
Use Existing (please select from list):	Password Entry
Create New (please enter a name) Enter new study name here	Please enter the password for Study: Testing 123
C Review Closed Studies	
C System Setup	
Ok Cancel	OK Cancel

Figure 135. When Studies have already been setup, the SM can select one of their own studies to Review, Edit or manipulate in some other way. Access to a Study is always password protected, as shown.

Depending on how the Administrator sets up the Password Expiry, you may see this warning message, Figure 136. This allows the password to be updated and the message will not show again until a new expiration period has elapsed.

Password Expiry	
The password for this Study will expire in 1 days' time. Do you wis	h to change the password now?
	Change Password
Yes No	Change password for Study Testing 123
Change Password	Old Password:
Study password has been changed	New password:
	Confirm New Password
OK	
2	OK Cancel

Figure 136. Password Expiry warning allows password update.

Make sure to tell the Scorers the new password. Make sure that Datasets are stored in Shared or Common Access Folders.

Once you have gained access to the Study, the LMA or SM will be able to Manage Scorers, Manage Datasets and Manage Protocols.



Komet[®]

7.9 The Study Menu - Existing Studies

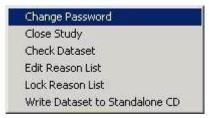


Figure 137. The Study Menu in SM mode.

SM users have full Study access capabilities allowing them to:

- Manage Study Access.
- Close a completed or compromised Study.
- Check Dataset integrity and Close it if it has been compromised.
- Manage the Reasons List (Reasons for actions which will be recorded in the audit trails).
- Write a Dataset to a formatted CD-R or CD-RW, which includes writing the Andor Database Viewer and support files for Viewing on other Windows PCs.
- Change Password

This function provides the same capability as the Password change shown in Figure 138, except the SM can instigate this at any time.

Note: All authorized Scorers should be informed of the changes.

.



7.10 Close Study

Studies can be closed for two basic reasons. Either the Study is complete and all relevant information has been recorded and communicated, or one or more of the study datasets or some other component of the Study has been compromised.

Warning	×
Are you sure that you want to 'Close' the current study? This action the application will then terminate. Restart to log onto a new study.	
Select Reason	
Please select a reason for the action:	
Close Study	
Dataset corrupted - assume opened illegally??	
Select	
Select	

Figure 138. Closing the Study prompts a warning of the actions which follow. A Reason must be specified for inclusion in the Audit Trail.

Closing the Study does not remove it from the system, but a message containing "Closed date, time, Reason and SM" is recorded in the Audit Trail and the Study no longer listed under Existing Studies. It now appears in the Closed Studies list and can only be accessed in the Review Closed Studies mode.

7.11 E-Signatures and Dataset Integrity

Datasets within a study contain the Raw Data. Each Dataset contains an Electronic Signature (esignature), which is an encrypted code word created from the entire contents of the file. If the file is tampered with in any way whatsoever, e.g. opened during scoring by another program, edited or inadvertently corrupted, then this e-signature will not match the file contents. Whenever KGLP or the Andor Database Viewer opens a Dataset the e-signature is checked and if in error you will be warned. This is an important aspect of ensuring data integrity and avoiding unauthorized changes to data – 21CFR part 11 compliance.

If Scoring is incomplete and a Dataset is found to be compromised, AND the file cannot be Repaired, KGLP will NOT allow Scoring to continue. The Dataset will have to be Recreated and Scored again.



The user will see a series of dialogs as shown in Figure 139. Successful Repair will yield a dialog indicating success.

Resume Experiment	
1	UsersiDocumentsiTesting123(Liver/Liver_RAW kdb' to resume,
Repair Create N	Resume Experiment
	Experiment Cancelled.
	ОК

Figure 139. Detection of a compromised Dataset on Resumption of Scoring leads to the options shown. Repair can be carried out if an intact Backup copy was selected for the Dataset. A New Dataset can be created and scoring started from scratch. The Scoring Session can be Cancelled or abandoned and the SM consulted. If the Scorer Cancels then the Cancel dialog is shown.

7.12 Check Dataset

Dataset integrity is checked and established whenever KGLP or Database Viewer Opens a Dataset for review or resumption of Scoring. The e-signature is checked against the file contents and a special internal encryption key in Komet GLP. The SM can initiate an integrity check in the current Study with this function.

Select the required Dataset	-
Selected Dataset: 001 Check Data Cancel	
Check Data Set Checking Dataset : 001	Dynamickomet
Please be patient this may take several minutes. Backup File Verified.	Database Verification success

Figure 140. First Select the Dataset to be checked, enter the access password, and then a Check will be activated. For large datasets with backups on network drives, this may take some time and the message box will appear, as shown. The result of the test will be delivered in a message box.

.



7.13 Scoring Datasets – Study Managers

The Study Manager is by default a Scorer in the Studies they manage. The Scoring process is identical for SDs and Scorers. To initiate scoring in an existing, open Study, the SM simply hits the Experiment button on the Komet Toolbar as shown in Figure 142 and also in Section 4 of this guide for Scorers.

III Live	Interactive	Delete Last	Express	Exgeriment	Dpen	
Background		la INA Tail DNA	Tail Length Oli	ve TM L/H	Extent TM	Experiment

Figure 141. Selecting the Experiment button on the Komet Toolbar will start preparation for scoring in the Current Study.

7.14 On Completion of a Dataset

When a Dataset has been fully scored, KGLP prompts the user to Create a Stand Alone CD. This process requires a CD-R, CD-RW or DVD-RW drive and a formatted disk installed. If the Scorer chooses to go ahead, then the Dataset is written to the preformatted CD-R, CD-RW or DVD disk along with all the support files and the Andor Database Viewer. The resultant disk will run in almost any Windows PC, and will provide a means to distribute, view, analyse and present the data on the disk. This is an ideal way to deliver the data to sponsors, archive and QA. A separate User Guide on the use of the Andor Database Viewer is provided with KGLP. The Database Viewer can be freely distributed.

Standalone CD Creation, can be activated for any Dataset, at any time, by selecting the function from the Study menu and making sure a formatted disk is in the drive.



7.15 Review Closed Studies Mode

lect Study Hello Administrator. You are a Study Administrator on this system.	/ Director and an		
Please select the Study for this sess	ion:		
C Use Existing (please select from	list):		
CO-0127-09445	select Study		
C Create New (please enter a nai	Please Select a C	losed Study	
Enter new study name here	-		
Review Closed Studies	H×04-099		
C System Setup	Study Direct M	or: IARK-LAPTOP\Administr	ator
	Number of	Password Entry	
ок с	2	Please enter the password	l for Study: HX04-001
	OK	OK	Cancel

Figure 142. Select the Review Closed Studies Mode from the Start up dialog, as shown. KGLP will display a List of Closed Studies. Select the Study you wish to review. Only Closed Studies which are managed by the current SM will be shown. As with all Study Access you will need to know the Study password for access.

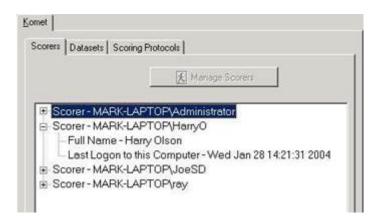


Figure 143. Once you enter KGLP, you will see the familiar information about the Study. However, while the Study is Closed you will NOT be able to access the Study Management tools.





Scorers	Datasets	Scoring Protocols
0	Komet D Data Data Back DExpe	umber of Samples - 30 umber of Slides - 2 ells per Slide - 50 ample Type - bladder
SCO	ers Datas	ets Scoring Protocols Manage Protocols
	Com Calib Calib	Protocol - LYMPHO et Options pration - 20x-green band era Settings Protocol - SYBR Gold CHO

Figure 144. You can review who scored the Study, the Dataset structures, passwords and the Folders, where the Study data was saved. You can also check Scoring Protocol settings.

7.16 View Audit Trails

File Study	View Audit Trails	Plug-Ins Help	
Current Co User: Study: Image	Study Scorers Dataset Protocol QC Setup QC Log	udy Director 4-001 4O10.TIF1	
Image		sed Study Review Mode	2

Figure 145. From the View Audit Trails dialog, the SM can access all the information relating to the Study.

Komet®



Study file	for Study: HX04-001	
Time: 8:01:03	PM	
Study Director:	WX20005DIRL\mbrowne	1.00
Password	helix3	
Paceword Channe	d: 1/31/2004 8:01:03 PM	
Number of Datase	te 0	
Study Closed:	No	
orady closed.	NO	
Amended by	WX20005DIRL\mbrowne	
Date: 1/31/2		
Time: 8:02:12		
	WX20005DIRL\mbrowne	
Password:	helx3	
	d: 1/31/2004 8:01:03 PM	
Number of Datase		
Number or Datase	Na. I	
Study Closed:	No	
study closed.	No	
and a day	WX20005DIRL\mbrowne	
Date: 1/31/2		
Time: 8:02:46		
	WX20005DIRL\mbrowne	
Password	helix3	
assword Change	d: 1/31/2004 8:01:03 PM	
Number of Datase	its: 2	
001		
002	10. · · · · · · · · · · · · · · · · · · ·	
Study Closed:	No	
current]	1. A (000000 (D) ()	
	WX20005DIRL\mbrowne	
Date: 2/16/2	JU4	
Time: 5:20:55		
	WX20005DIRL\mbrowne	
Password	helix3	
	d: 1/31/2004 8:01:03 PM	
Number of Datase	ts: 2	
001		
102	2. 2. A2	
study Closed:	Yes : Closed - Dataset corrupted - assume opened i	illegally?? : 2/16/2004 5:20:55 PM
		1
		and the second s
	Close	Print

Figure 146. The Study audit trail displays the Study audit information. It records the date, time and user who closed the Study.

7.17 Reopen the Study

The Study Menu in this mode allows you to Select another Study for Review, using the first menu item.

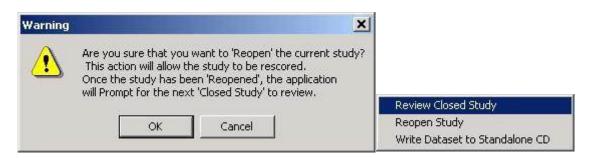


Figure 147. Choose the Reopen Study menu to Open the Study again.

You must confirm this action as the dialog shows and enter the password see Figure 148.



Password Entry	
Please enter the passw	ord for Study: HX04-001
1	
ОК	Cancel

Figure 148. Reopen requires entry of the Study Password as a security measure.

Passwor	d Expiry
?	The password for this Study will expire in 3 days' time. Do you wish to change the password now?
	<u>Y</u> es <u>N</u> o

Figure 149. If the password has expired a warning will be given. Now you can change the password for future access.

ielect Re	ason
Please	select a reason for the action:
Reop	en Study
Datase	t corrupted - assume opened illegally?? 💽
Datase	t corrupted - assume opened illegally??

Figure 150. Reopen Study will generate an entry in the Study Audit Trail and therefore requires a Reason to be entered.

Once a Study has been re-opened it is possible to Access all Study Management tools. The newly opened Study can be edited or added to exactly as it could before it was closed. It can also be used as a Template for creating a new Study.



7.18 Write Dataset to Standalone CD

Datasets can be written to a Standalone CD, using the appropriate menu, as shown in Figure 151. This allows you to redistribute the Dataset for any purpose including for example, review or training.

N	Kinetia	Imaging Komet 6.0 GLP	- 🗆 🗙
Fle	Study	View Audit Trails Plug-Ins Help	
a		iew Closed Study pen Study tor	<u> </u>
	biuuy.	e Dataset to Standalone CD	
-	mage	: KOMETDEMO10.TIF1	
		Closed Study Review Mode	9
Ir	nage	Calibration:X*0.825 : Y*0.825	
			-

Figure 151. Select the menu to Write Dataset to Standalone CD.

elect Dataset Select the require	d Dataset	
 ♥ Komet Dat ♥ Komet Dat 	aset-001	
Selected Dat	aset	
Select	Close Data	Cancel

Figure 152. Select the Dataset you wish to write.

Make sure you have a formatted CD-R or RW disk in the CD-R or RW drive. We recommend the use of UDF format for Direct CD Writing.



SECTION 8 KOMET GLP FOR SCORERS

8.1 Preparing to Score Datasets in Studies

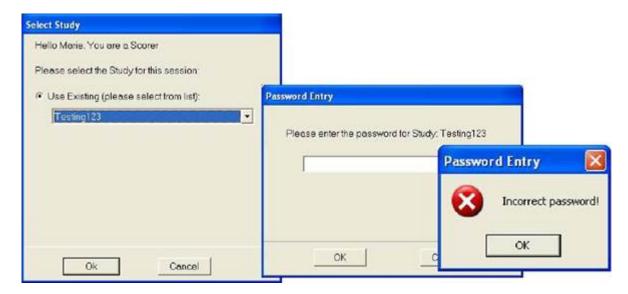


Figure 153. Start KGLP and Select a Study from the List shown. Only Studies in which you are an authorized Scorer will list. You must enter the correct password or you will receive a warning and KGLP will close.

If the Study Password has expired or has been changed you should contact your SM for details.

Once you gain access to the Study you will see the Komet Control and Image Windows. The Control Window will display a Message in Scoring Mode, in blue as follows 'Press Experiment to Start Scoring'.

III Live	Interactive	Delete Last	Express	Exgeriment	Dpen	
Background		la INA Tail DNA	Tail Length Oli	ve TM L/H	Extent TM	Experiment

Figure 154. To Start Scoring an Open Study, Select the Experiment button from the Komet Toolbar.

.





denum	
odenum	Duo denum 🔀
	The password you have entered is incorrect
	ordenum ordenum Inter password for dataset.

Figure 155. You will then be asked to Select a Dataset where your data will be stored. You must enter the correct Dataset Password or you will see a warning like the above. The password protection of Datasets allows assignment of individual Scorers to specimen

types.

If you enter a correct password KGLP will then either ask you to Select a Protocol for scoring the Dataset (new Dataset) or automatically select the one previously used.

elect the required Protocol	
Kornet Protocol - SYBR Gold Gold Syber Gold S	
Selected Protocol: SYBR Gold	

Figure 156. New Datasets require that you select a Scoring Protocol. Check with your Study Manager if you are unsure.





Comet (Options	×
1	The magnification for this Protocol is x20. Please ensure you have the correct objective selected before go	ing Live.
	OK Cancel	

Figure 157. A Magnification reminder is shown – Make sure you have the correct objective in place and make sure the microscope beam-splitter is in the correct position to send an image to the camera.

N Input Slide Information	
Please enter the ID codes for the	e next slide: —
SlideID: 001	IT Autonumber
ReplicateID: A	Autoletter
ОК	Cancel

Figure 158. The Side ID should be entered next.

Depending on the numbering scheme employed and the number of replicate slides per sample, the Autonumber or Autoletter check box, if selected, automatically updates the Replicate ID, as you move from slide to slide. You will be able to change this after scoring the slide if you make an error.

ease enter the following	identifiers for this PC setup
Microscope ID:	Oly-03-01
Camera ID:	dummy
PC Identifier:	DJXM9341

Figure 159. If the SM has selected the Require Equipment ID, you will get this dialog box.

You should at this point enter the equipment IDs which will be used for scoring the Study. Take care with these since KGLP cannot check their validity.

Each time Scoring is resumed on the Dataset you will have to enter them. After the first time these are entered you will be prompted with the previous values, which you can rectify or accept.



Autonumber will provide a postfix incrementing from your initial value e.g. 001-1, 001-2 or 001-10, 001-11; while Autoletter will postfix in the format: 003-A, 003-B etc.

8.2 Scoring Datasets

Once the Dataset has been accessed, the Scoring Protocol selected, Slide ID and the Equipment IDs entered, then you are ready to start scoring the comet slides.

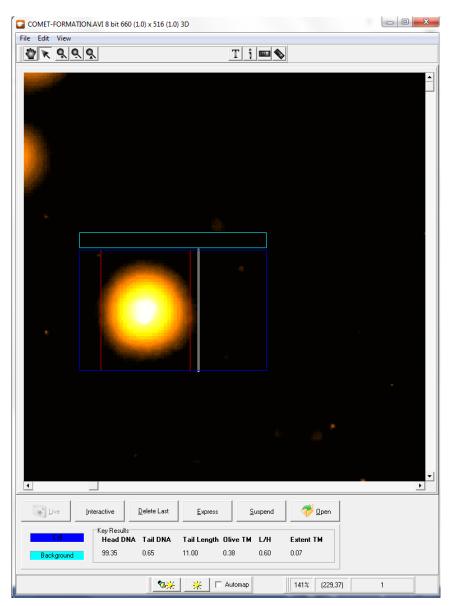


Figure 160. Now you are ready to start scoring. KGLP will automatically switch to Live (if you are scoring Live from the microscope) and initialize the camera to the correct settings specified in the Protocol. Move the ROI box over the comet and click the Right mouse button. KGLP will analyse the comet automatically and post the results to the analysis graphic in the Control panel as shown in Figure 162.

Scan the slide in a systematic raster or scrolling pattern and select fields with cells suitable for scoring. Do not choose cells by any criteria except that they are scorable, without debris in the

Komet[®]



background and avoid overlapping cells. This scanning method will remove the chance of scoring the same cell more than once, which would bias results. In a future release we will provide a motorised scanning stage option to remove this potential bias.

Important Note: If you pause in Scoring for longer than the System Timeout period (measured in minutes from 1-30), then KGLP will lock you out of Windows and show a message indicating that KGLP has locked you out. To regain access you will have to logon on to Windows again by pressing CTRL+ALT+Delete keys and perform a standard timeout logon.

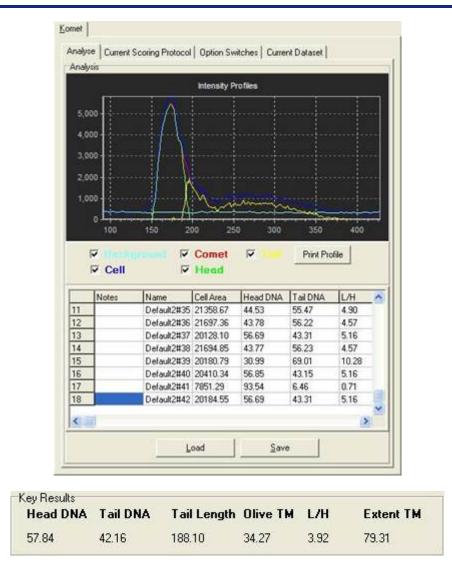


Figure 161. Results of Analysis are posted to the Control Window, including an Intensity Profile and quantification results.



8.3 Key Results and Current Context

Cell: 18	Exp. ID: 001
Data: 43	Slide ID: 001_B
Image Calibi	ation:X*1.12 : Y*1.12
Protocol:	SYBR Green (Mag: x20)

Figure 162. Key Results are posted to the Key Results area on the Komet toolbar and the progress of Scoring in the current session is shown in the Current Context area of the Control Window.

Come	Current Scoring Protocol t Protocol	option of morico [outer outer [
	YBR Gold Comet Options Calibration - 20X Green Camera Settings	Band	
Protocol C	ptons		
0			
Operatio		-	
🔽 Auto k	nal background xpress on Store IT Interac	tve Tail	
Auto k	ackground		
 Auto k AutoE Head Calibratic Apply 	ackground xpress on Store Γ Interac Ellipse On Γ Flip on	Open	Camera Settings

Figure 163. During scoring the Current Scoring Protocol can be Viewed, and, if absolutely necessary, the Camera settings can be adjusted. If the Scorer changes the camera settings, then he/she must provide a Reason for the change, which may be subject to audit.

The settings shown below the tab, Option Switches, can be changed; however, these will not affect analysis results. See the Komet[®] User Guide for details.

.



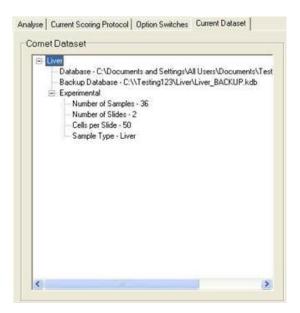


Figure 164. During scoring, the Current Dataset structure and storage location can be viewed by selecting the appropriate tab in the Control Window. These settings can only be changed by the SM.

8.4 When the Current Slide is Complete

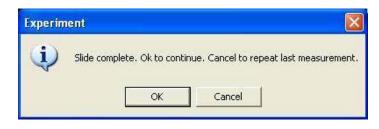


Figure 165. When you have Scored the correct number of cells for the current Slide, KGLP informs you to accept or reject the last cell.

If you accept the last cell of the slide, then KGLP proceeds with the checks needed to complete recording of all the data from that slide into the Dataset. If you reject the cell for any reason, the cell is deleted and you will be asked to provide a Reason for the deletion. You will have to re-score the cell or score another cell for inclusion into your Dataset. Rejection of a cell always requires a Reason.

If you accept the cell then the dialog in Figure 166 is shown, where you must confirm the Slide ID or correct it as shown below.

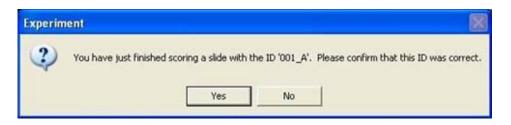


Figure 166. When you have Scored the correct number of cells for the current Slide, KGLP asks you to confirm the Slide ID. If you choose NO the dialog in Figure 167 will be shown.





N Input Slide Information	
Please enter the new ID codes	for the previous slide:
SlideID: 001	
ReplicateID: A	Autoletter
ОК	Cancel

Figure 167. To edit the Slide Information simply enter the correct values into the Slide ID and Replicate ID text boxes.

If the SM has enabled the function, you will then be asked if you wish to view the Image Gallery for this Slide, or the current Sample if more than one Slide has been scored from the Sample.

e Gallery		
Do you wish to view	the image gallery f	or this Slide?
Yes	No	
	e Gallery Do you wish to view Yes	Do you wish to view the image gallery f

Figure 168. The prompt to view the Image Gallery is shown next.

If you select Yes then an Image Gallery of comets in the current Slide or Sample is shown. See Figure 170. The Gallery can then be Viewed (and Printed if required).

N Input Slide	Information	
Please ente	r the ID codes for th	ne next slide:
SlideID:	001	-
ReplicateID	1	F Autonumber I⊽ Autoletter
	ж _	Cancel

Figure 169. The Autoletter or Autonumber will update the Replicate ID text box automatically, but the Slide ID will have to be entered manually.

Once entered click OK to start Scoring. Cancel provides an option to Suspend the current Scoring Session.



8.5 Image Gallery for Current Sample and Printing

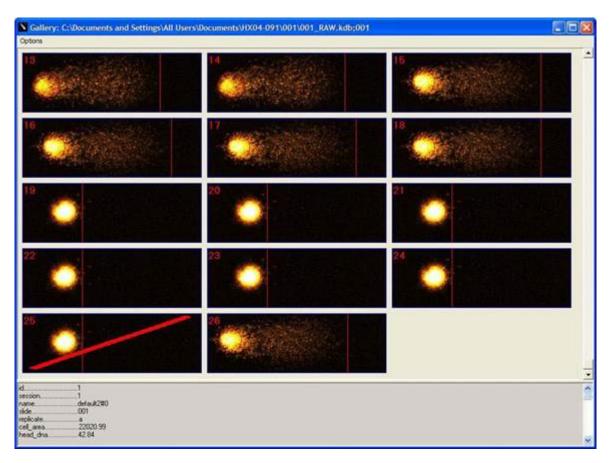
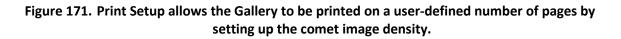


Figure 170. The Image Gallery shows an image of every cell scored in the current slide (and Replicates). Callipers (vertical red lines) show the start and end of the comet Head and Tail regions as analysed byKGLP. If 'Head Ellipse' is selected in Comet Options, then an ellipse will be shown delineating an estimate of the Head region. Deleted Cells are also recorded in the Gallery, but their omission/rejection from data analysis is indicated with a single diagonal Red Line as shown in cell 25 above.

Red numbers are shown on the individual images, indicating the data position in the current Dataset. The Options menu allows the Gallery to be printed for recording purposes. The Print Setup dialog, shown in Figure 171, allows the Scorer to select the density of images set onto a page. This allows for example, adjustment to one Sample per printed page.



Image Print Setup	
Select Layout for Images	
Total Number of Images: 100	
Number of Columns per page:	7
Number of Rows per page:	15 1
Number of Pages: 1	
ОК	ancel



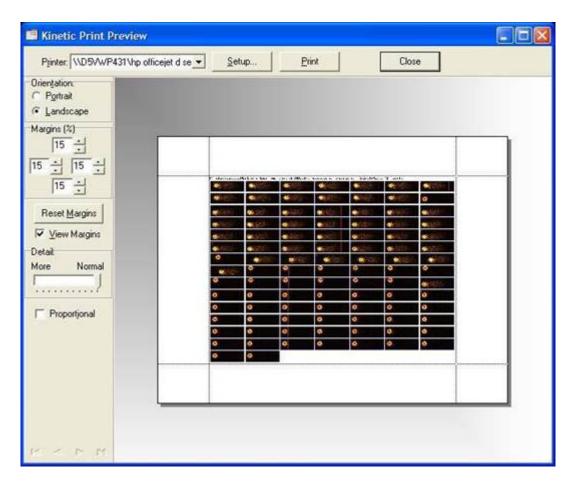


Figure 172. The Print Preview dialog allows for printer selection and setup and shows the Printer Page output prior to printing.



8.6 Deleting Last and Interactive Scoring



Figure 173. Heavily damaged cells may occasionally cause inaccurate analysis, but the Scorer can override the system in this case.

Selecting the Delete Last button on the Toolbar will mark the last cell for rejection in the Dataset and Interactive can be selected. In Interactive mode the Toolbar changes to the lower panel shown in Figure 173. The Scorer drags out a rectangular ROI holding down the left mouse button. The left edge of the ROI indicates the end of head and start of tail, while its right hand edge indicates the end of the tail. KGLP will use the left and right edges to constrain intensity analysis and re-compute the results. After Delete Last or Interactive, the user must specify a Reason for the Action from the Reason list dialog in Figure 174.

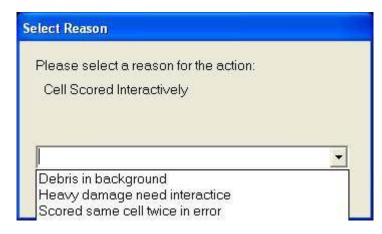


Figure 174. The Reason List provides a set of valid Reasons for specific actions, like Delete Last, Change Camera settings and Interactive. It provides an important mechanism for monitoring scoring.

Komet®

8.7 Suspending the Scoring Session

You can Suspend the Scoring Session at any time, by clicking on the Suspend button on the Komet Toolbar. Other actions which lead to Suspend are Cancelling dialogs at key points in data entry prior to the start of scoring.

Suspending a Scoring Session will make an entry in the Audit Trail as described in Figure 175. The Suspended Dataset details can be reviewed in the Dataset window the next time you score in KGLP.

Experiment		Reason
(į)	This action will Suspend your current experiment! Are you sure?	Enter reason for Suspend
	Yes No	OK Cancel

Figure 175. Scoring can be Suspended at any point. The Toolbar Suspend Button shows the Suspend Dialog. If the Scorer confirms the action then a Reason Entry box is shown, where a free form text field can be added. This information is written into the Dataset Audit Trail, along with other information about the Scoring Session, such as Start and End Time and the IDs of slides scored in the Session.

8.8 Resume Scoring of a Dataset

When you start KGLP to Resume scoring a Dataset, go through the same steps described previously to access the appropriate Study. To Resume Scoring, click the Experiment button on the Komet Toolbar and the following sequence of dialogs will be presented:

Select the required Dataset	
 S Komet Dataset - 001 Komet Dataset - 002 	
	001
	Enter password for dataset.
	OK Cancel
Selected Dataset: 001	
Select Close D	ata Cancel

Figure 176. Choose the Dataset and enter the Dataset password.





Select Protocol	
Select the required Protocol	
 Komet Protocol - EthBr Komet Protocol - SYBR Green 	
Comet Options The magnification for this Protocol	is x20. Please ensure you have the correct objective selected before going Live.
Selected Protocol: SYBR Green Select Cancel	

Figure 177. Choose the Scoring Protocol, KGLP will remind you to select the correct microscope objective to make sure the images are correctly calibrated.

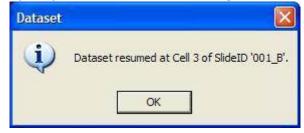


Figure 178. If scoring is Resumed mid-slide, KGLP will remind you of the Slide ID, Replicate ID and Cell number. Be sure to replace the correct slide in the correct position on the stage for scoring. You may want to record slide coordinates to ensure this.

Please enter the ID codes for the	he next slide:
SlideID: 009	-
ReplicatelD:	Autonumber

Figure 179. If Scoring Resumes after the end of a slide, you will need to enter the next Slide ID or Replicate ID when you Resume.

If the SM has set the Require Equipment IDs then make sure the correct values are entered in the Equipment ID dialog.

When KGLP Resumes Scoring the Dataset, it checks the integrity of the file(s) through the esignature. If the integrity check fails you will be given the opportunity to Repair the file(s) from the valid copy – either Raw or Backup.



KGLP also checks the Dataset access rights, by trying to write a test file to the folder(s) containing the file(s). If the folder cannot be written the following messages will be shown and you will have to gain access to the folders before scoring. This must be arranged by the SM or LMA.

Open D	atabase 🔀
į)	Cannot Open Original Database 'C:\Documents and Settings\mbrowne.WX20005DIRL\Desktop\+004-091\001\001_RAW.kdb'. This file has restriked accessed. please contact your 'Administrator'.
Komet D	atabase
8	Unable to open database: C:\Documents and Settings\nbrowne.WX20005DIRI.\Desktop\FX04-091\001\001_RAW.kdb. Experiment Cancelled OK

Figure 180. If the Scorer does not have access to the specified Dataset Folders, then KGLP will advise and will not allow scoring to proceed.

8.9 Writing Stand Alone CD - On Completion of a Dataset

When a Dataset has been fully scored, KGLP prompts the user to Create a Stand Alone CD. This process requires a CD-R, CD-RW or DVD-RW drive with a formatted disk installed. If the Scorer chooses to go ahead, then the Dataset is written to the preformatted CD-R, CD-RW or DVD disk along with all the support files and the Andor Database Viewer. The resultant disk will run in almost any

Windows PC, and will provide a means to distribute, view, analyse and present the data on the disk. This is an ideal way to deliver the data to sponsors, archive and QA. A separate User Guide on the use of the Andor Database Viewer is provided with KGLP. The Database Viewer can be freely distributed.

Kinetic Imaging Komet 6.0 GLP							
File	Study	QC V	iew Audit Trails	Wizard	ds Plug-Ins	Help	
-0	Clos Che Edit Loci	nge Pas e Study ck Datas Reason c Reason te Datas	et List	e CD	tor + Adm		

Figure 181. Standalone CD Creation, can be activated for any Dataset, any time, by selecting the function from the Study menu, making sure a formatted disk is in the CD-R, CD-RW, DVD-RW drive.



PROTOCOL FOR THE APPLICATION OF THE pH>13 ALKALINE COMET ASSAY TO THE DETECTION OF DNA DAMAGE IN MAMMALIAN CELLS

Prepared by:

Marie Vasquez, B.S. Helix3 Inc.

www.helix3inc.com

June 14, 2005 revised

The alkaline (pH>13) Comet assay, developed by N. P. Singh¹, combines the simplicity of biochemical techniques for detecting DNA single strand breaks (frank strand breaks and incomplete excision repair sites), alkali-labile sites, and crosslinking with the single cell approach typical of cytogenetic assays. The advantages of the comet technique include: (1) the collection of data at the level of the individual cell, allowing for more robust types of statistical analyses; (2) the need for small numbers of cells per sample (<10,000); (3) its sensitivity for detecting DNA damage; and (4) that virtually any eukaryote cell population is amenable to analysis. This protocol describes the methodology as used in our laboratory. Clearly, there are other methods to process samples, prepare gel slides, and analyse DNA damage; we would appreciate hearing about them.

The purpose of this protocol is to provide instruction on the application of the alkaline Comet assay to the detection of DNA damage. For recent guidelines see Tice, R.R., E. Agurell, D. Anderson, B. Burlinson, A. Harmann, H. Kobayashi, Y. Miyamae, E. Rojas, J-C Ryu, and Y.F. Sasaki (2000) The single cell/gel comet assay: Guidelines for *in vitro* and *in vivo* genetic toxicology testing. Environ. Mol. Mutagen. 35: 206-221.

¹ Singh, N.P., McCoy, M.T., Tice, R.R. and Schneider, E.L.: A simple technique for quantitation of low levels of DNA damage in individual cells. Exp. Cell Res. 175: 184191, 1988.

More information about the Comet Assay is available to all interested individuals at www.helix3inc.com or www.cometassay.com.



SECTION 9 COMET ASSAY

9.1 Preparation of Reagents

Materials:

- Dimethylsulfoxide (DMSO)
- Disodium EDTA
- Hanks Balanced Salt Solution (HBSS), (Ca $^{2+}$, Mg $^{2+}$ free)
- Phosphate Buffered Saline (PBS), $(Ca^{2+}, Mg^{2+} free)$
- Sodium Chloride (NaCl)
- Sodium Hydroxide (NaOH)
- Nucleotide stain (e.g., SYBR™ or Ethidium)
- Triton X-100
- Trizima Base

Procedure:

- A. <u>Mincing Solution: HBSS (Ca²⁺, Mg²⁺ free) with 20mM EDTA and 10% DMSO:</u> To 400 ml 1x HBSS (Ca²⁺, Mg²⁺ free), add 3.72 g EDTA and 50 mL DMSO, adjust pH to 7.0 7.5., q.s. to 500ml, store at room temperature.
- B. <u>PBS (Ca²⁺</u>, Mg²⁺ free): Dulbecco's PBS 1 L packet: add 990 mL dH₂O, adjust pH to 7.4, q.s. to

1000 ml, store at room temperature.

C. Lysing Solution: Ingredients per 1000ml: 2.5 M NaCl (146.1 g), 100 mM Na₂EDTA (37.2 g), 10 mM Trizma base (1.2 g)

Add ingredients to about 700 mL dH₂O, and begin stirring the mixture. Add ~8 g pelletized NaOH and allow the mixture to dissolve (about 20 min). Adjust pH to 10.0 using concentrated HCI or NaOH. With dH₂O q.s. to 900 mL (the Triton X-100 and DMSO will increase the volume to the correct amount), store at room temperature.



Final lysing solution:

Add fresh 1% Triton X-100 and 10% DMSO, and then refrigerate for at least 30 minutes prior to slide addition.

NOTE: The purpose of the DMSO in the lysing solution is to scavenge radicals generated by the iron released from haemoglobin when blood or tissues are used. It is not needed for other solutions or where the slides will be kept in lysing for a brief time only. Sodium lauryl sarcosinate, which we used earlier, is not included in this version.

D. Electrophoresis Buffer (300 mM NaOH / 1mM EDTA):

Prepare from stock solutions: 1. 10 N NaOH (200g/500 ml dH₂0)

2. 200 mM EDTA (14.9 g/200 ml dH₂O, pH 10)

Store both at room temperature. We prepare the NaOH and EDTA stock solutions every ~2 weeks.

For 1X Buffer (made fresh before each electrophoresis run): per litre, add 30 mL NaOH and 5.0 mL EDTA, q.s. to 1000 mL, mix well. The total volume depends on the gel box capacity. Prior to use, measure the pH of the buffer to ensure >13. Use this buffer at pH12.1 to selectively assess for strand breakage.

E. Neutralization Buffer:	0.4 M Tris - 48.5 g added to ~900 ml dH ₂ O, adjust pH to 7.5 with
	concentrated (>10M) HCI: q.s. to 1000 ml with dH_2O , store at room temperature.
F. <u>Staining Solution:</u>	Ethidium Bromide (10X Stock - 20μg/ml): add 10 mg in
	50 ml dH ₂ O, store at room temperature. For 1X stock - mix 1 ml with 9 ml dH ₂ O.
	SYBR™: Add 1µl of provided stock solution to 10 ml of
	Tris Buffer. pH 7.5 to prepare a 1:10,000 dilution. Make fresh prior to use. Staining solution is stable for several hours at room temperature and for 1-2 days when refrigerated. pH is critical for stability.

CAUTION: Handle dyes appropriately.

NOTE: SYBR[™] offers a greatly increased signal to noise ratio. However, as the stain fades rapidly upon exposure to light, the DNA must be stained just prior to analysis.



9.2 Preparation of Slides for the Comet Assay

Materials:

- Coverslips (No 1, 24 x 50 mm)
- Low Melting Point Agarose (LMP)
- Microcentrifuge Tubes
- Micropipettor and Tips
- Microscope Slides
- Microscope Slide Tray
- Normal Melting Point Agarose (NMA)
- PBS (Ca²⁺, Mg²⁺ free)
- Scintillation vials
- Frozen Ice Packs

Procedure:

- 1. Prepare 0.5% LMP (250 mg per 50 ml) and 1.0% NMA (500 mg per 50 ml) in PBS. Microwave or heat until near boiling point and the agarose dissolves. For LMP, aliquot 5ml samples into scintillation vials (or other suitable containers) and refrigerate until needed. When needed, briefly melt agarose in microwave or by another appropriate method. Place LMP vial in a 37°C water bath to cool and stabilize the temperature.
- 2. While NMA agarose is hot, dip slide up to one-half the frosted area and gently remove*. Wipe underside of slide to remove agarose and lay the slide in a tray on a flat surface to dry. The slides may be air dried or warmed for quicker drying. Store the slides at room temperature until needed; avoid high humidity conditions.
- * Procedure based on information provided in M. Klaude, S. Eriksson, J. Nygren, and G. Ahnstrom (1996) The comet assay: mechanisms and technical considerations, Mutation Res. 363, 89-86.

NOTE: Slides should be labelled prior to dipping.



The Comet Assay Laboratory Protocol

- 3. To the coated slide, add 75µl of LMP (37°C) mixed with ~10,000 cells in ~5-10µl (do not use more than 10µl). Replace coverslip and place the slide on a slide and keep cold until the agarose layer hardens (~3 to 5 minutes). Using ~10,000 cells in an area 25 by 50mm results in ~1 cell per microscope field at 250x magnification.
- 4. Gently slide off coverslip and add a third agarose layer (75μl LMP) to the slide. Replace coverslip, return to the slide tray and keep cold until the agarose layer hardens (~3 to 5 minutes).
- 5. Remove coverslip and slowly lower slide into <u>cold</u>, freshly made Lysing Solution. Protect from light and refrigerate for a minimum of 1 hour. In our experience, slides may be stored for at least 4 weeks in cold Lysing Solution without affecting the results.
- **NOTE:** The amounts indicated are based on using 24 x 50 mm coverslips. Proportional volumes can be used for coverslips differing in size. If the gels are not sticking to the slides properly, avoiding humidity and/or increasing the concentration of NMA agarose in the lower layer should eliminate the problem. Steps 3 to 5 should be performed under dim yellow lights to prevent DNA damage.



9.3 Cell Isolation

- A. <u>Whole Blood</u>: Per gel slide, mix \sim 5µl whole blood with 75µl LMPA and process accordingly.
- NOTE: A small volume of blood can be added to 1 ml of media and stored cold (refrigerated or on ice) for an extended period of time until processed. In this case, the cells must be pelleted by centrifugation and care taken to remove as much supernatant as possible before adding the 75µl LMP per 5µl of blood.
- B. <u>Isolated Lymphocytes</u>: Micro technique: Mix 20 μl whole blood with 1 ml RPMI 1640 in a microcentrifuge tube, add 100 μl Ficoll below the blood/media mixture. Spin for 3 minutes at 2000 rpm. Remove 100 μl of bottom of the media/top of Ficoll layer, add to 1 ml media and mix, spin for 3 minutes to pellet lymphocytes. Pour off supernatant, re-suspend pellet in 75μl LMPA, and process accordingly.
- C. <u>Bone Marrow</u>: Perfuse a femur (mouse) with 1 ml of cold mincing solution (HBSS with 20 mM EDTA, 10% DMSO) into a microcentrifuge tube. Remove and mix 5 μl per 75 μl LMP, and process accordingly.
- D. <u>Solid Organs</u>: Place a small piece of an organ in 1 ml of cold HBSS containing 20 mM EDTA/10% DMSO. Mince into fine pieces, remove and mix 5 10 μ l of cell suspension with 75 μ l LMP, and process accordingly.
- NOTE: For blood rich organs (e.g., liver), mince into large pieces, let settle, aspirate mincing solution, add fresh mincing solution, mince into finer pieces, remove and mix 5 μl of the cell suspension with 75 μl LMPA, and process accordingly. The purpose of the DMSO is to prevent lipid peroxidation associated with the processing of some tissue. The volume of the cell suspension to mix with 75 μl of LMP must be 10 μl or less, while the optimal cell number is ~10,000 cells per slide. Adjust the volume of mincing solution as needed.



- E. <u>Monolayer Cultures</u>: Remove the media and replace with mincing solution, scrape off cells into the mincing solution using a Teflon scraper and adjust concentration to approximately 1x10 cells/ml. Remove and mix 5 - 10 µl of the cell suspension per 75 µl LMP and process accordingly.
- F. <u>Suspension Cultures</u>: Mix ~10,000 cells in 10 μ l or less volume per 75 μ l LMP and process accordingly.
- G. <u>Viability Assessment:</u> Viability assessment is often critical for interpreting Comet data. While several methods are available, we routinely use a technique developed by G.H.S. Strauss (Non-random cell killing in cryopreservation: Implications for performance of the battery of leukocyte tests (BLT) I. Toxic and immunotoxic effects, Mutat. Res., 252, 1-15, 1991).

Materials:

- Solution A: Ethidium bromide (EB) For the stock solution, dissolve 50 mg EB in 1.0 ml 100%-EtOH, and add 100 μ l to 4.0 ml PBS. For the working solution, add 250 μ L stock solution to 9.75 ml PBS (final conc. = 0.025 μ g/ μ l), and protect from light.
- Solution B: 5-6 Carboxyfluorescein diacetate (CFDA) For the stock solution, dissolve 3 mg CFDA in 1.0 ml acetone. For the working solution, add 420 μ l stock solution to 9.58 ml PBS (final conc. = 0.125 μ g/ μ l), and protect from light.
- Solution C: Working solution: Combine A and B in ratio of 1:1. Solution can be refrigerated for up to 6 months. Protect from light.



Procedure:

- 1. Place 40 μ L of at least 10⁶ cells/mL in a microcentrifuge tube, and add 10 μ l of working solution of dual stain (5-6 carboxyfluorescein diacetate: Ethidium bromide; solution C).
- 2. Let stand for 3-5 minutes at 37°C, and then remove excess stain by washing cells with 1 ml of wash solution (tissue culture media, PBS, etc.).
- 3. Pellet cell sample, and pour off supernatant.
- 4. Repeat wash. The cells may be stored by placing pelleted cells in PBS at 4°C until ready to store.
- 5. Pellet cells, and pour off supernatant. Mix cells in remaining drop of fluid (about 10 μ l).
- 6. Place drop with cells on microscope slide, add coverslip (the smaller in size the better).
- 7. Observe cells with fluorescent microscope (FITC of Acridine orange filter combination).
- 8. Score 100 cells for the number of viable cells (green fluorescence in cytoplasm), the number of compromised cells (green fluorescence in cytoplasm, red fluorescence in nucleus), and the number of dead cells (red fluorescence in nucleus).
- NOTE: In this staining technique, a live cell will metabolize 5-6 Carboxyfluorescein diacetate to a fluorescence compound, turning the cytoplasm green. Ethidium bromide will stain the nucleus of any cell with a damaged cell membrane. Excessive exposure to UV light will kill cells, resulting in uptake of Ethidium bromide (red stain), if available. In compromised cells, the intensity of cytoplasmic stain will vary greatly depending on how metabolically competent is the cell.



9.4 Electrophoresis of Microgel Slides

Materials:

- Coverslips (No. 1, 24 x 50mm)
- Micropipettor and Tips

Procedure:

- NOTE: The procedure described is for electrophoresis under pH>13 alkaline conditions. Electrophoresis can also be conducted at pH 12.1 or under neutral conditions. Also, slides can be removed from lysis to treat with repair enzymes specific for different classes of DNA damage or with proteinase K to remove any residual protein, or can be used under non-electrophoretic conditions to detect apoptotic/necrotic cells.
- 1. After at least 1 hour at ~4°C, gently remove slides from the Lysing solution. Place slides side by side on the horizontal gel box near one end, sliding them as close together as possible.
- 2. Fill the buffer reservoirs with freshly made pH>13 Electrophoresis Buffer until the liquid level completely covers the slides (avoid bubbles over the agarose).
- 3. Let slides sit in the alkaline buffer for 20 to 60 minutes to allow for unwinding of the DNA and the expression of alkali-labile damage.
- **NOTE:** The longer the exposure to alkali, the greater the expression of alkali-labile damage.
- 4. Turn on power supply to 0.7v/cm (anode to cathode) and adjust the current to the required milliamperes (mA) by raising or lowering the buffer level.
- **NOTE:** The mA requirement is not a set value. The mA is used as a reference point that should be consistent between runs to reduce run to run variability.

Depending on the purpose of the study and on the extent of migration in control samples, electrophorese the slides for 10 to 40 minutes.

- NOTE: The goal is to obtain migration among the control cells without it being excessive. The optimal electrophoresis duration differs for different cell types. If crosslinking is one of the endpoints being assessed then having controls with about 25% migrated DNA is useful. A lower voltage and a longer electrophoresis time may allow for increased sensitivity.
- 5. Turn off the power. Gently lift the slides from the buffer and submerge in Neutralization Buffer, let sit for at least 5 minutes



- **NOTE:** Slides may be stained with 100 μl 1X Ethidium Bromide or SYBR[™] stain and scored immediately or dried as in step 6.
- Drain slides, expose them briefly to cold 100% ethanol or cold 100% methanol, and allow to dry.
 Store in a dry area.
- 7. When convenient, flood the slide with stain (\sim 300-500 µl) before placing coverslip to ensure even hydration and staining and cover with a fresh coverslip.

Before viewing slides, blot away excess liquid on the back and edges. After scoring, remove coverslip before storing.



9.5 Evaluation of DNA Damage

 For visualization of DNA damage, observations are made of fluorochrome-stained DNA using a 20-25x objective on a fluorescent microscope.

NOTE: Depending on the size of the cells being scored, other objectives can be used (i.e. 16x, 40x).

- 2. Although any image analysis system may be suitable for the quantitation of Comet data, we use a Komet analysis system developed by Andor Imaging linked to a CCD camera to quantitate DNA migration. Finally, the program calculates tail moment. Generally, 50 to 100 randomly selected cells are analysed per sample.
- 3. Statistical analysis for *in vitro* data is based on 100-150 cells per unit of exposure.
- NOTE: Although we use an image analysis system to collect the data, other approaches can be used. For example, the frequency of round cells, cells with short tails, cells with medium long tails, and cells with long tails can be counted among 100 cells and analysed. To evaluate if the assay is functioning correctly, a control and treated (ex. 50 cGy x-ray), cryopreserved population of cells (e.g., blood lymphocytes) can be used. The treated cell population can be stored as 1 mL aliquots (at a density of 100,000 cells per ml) into microcentrifuge tubes. Whenever cells are being electrophoresed, one of these "internal control" cell samples can be processed along with the study population, and analysed for DNA damage. The Comet data resulting from this population can be used to evaluate interrun variability.



9.6 Model Experiment: Human Leukocytes and Hydrogen Peroxide

This experiment consists of a comparison of basal and chemically-induced levels of DNA damage among human leukocytes. For the former, DNA migration among cells electrophoresed for either 20 or 40 minutes will be compared. For the latter, leukocytes will be exposed to H_2O_2 , an excellent inducer of DNA damage. Because of its ease of use, H_2O_2 makes for a good positive control while the Comet assay is being developed.

- 1. Transfer 20 μ l of whole blood to 1 ml of HBSS or 1 ml of HBSS containing 200 μ M H₂O₂ in a microcentrifuge tube in duplicate per treatment group.
- 2. Place samples in a refrigerator or in a 37°C incubator for 60 minutes.
- 3. Prepare 2 lysing jars (50 ml Coplin jar containing 40 ml lysing solution) and place in the refrigerator.
- 4. Label four slides for each tube, dip slides in regular agarose, and allow to dry.
- 5. At the end of the treatment period, pellet the cells, remove supernatant, re-suspend cells in HBSS, and remove 500 μ l for viability assessment.
- 6. Pellet the cells for Comet analysis.
- 7. After removing as much of the supernatant as possible, re-suspend each pellet in 300 μ l 0.5% LMP and layer 75 μ l onto each of four labelled microscope slides. Place the slides on the cold slide tray.
- 8. After allowing the agarose to solidify, add the final layer of 0.5% LMP (75 μ l). Replace the slides on the cold slide tray.
- After this layer solidifies, remove coverslips and place slides back to back vertically in a 50 ml Coplin jar containing 4 ml of cold Lysing Solution. Place the Coplin jars in the refrigerator for a minimum of 1 hour.
- 10. Transfer 2 slides from each tube within each treatment group (8 slides total) to the gel box, expose the DNA to alkali for 20 minutes, and then electrophorese for 20 minutes.
- 11. Neutralize with Tris buffer and stain with Ethidium bromide or another appropriate dye (wear disposable gloves for this step).



12. Score 50 cells per duplicate slide for SNA migration using image analysis, or score 100 cells per slide by eye, categorizing cells as undamaged (i.e., no migration), short migration, medium migration, long migration, and complete migration (i.e., no nucleus remaining).

13. Repeat steps 9 through 11 using 40 minutes of electrophoresis.

14. Compare the amount of migration per cell, the number of cells with increased migration, the extent of migration among "damaged" cells, and viability.



APPENDIX 1 INSTALLATION OF SOFTWARE

A1.1 Evaluation license

If you are evaluating Komet then you will need to obtain a license from an authorized Andor representative. In order to create the license we need to send by email the exact Computer Name (see Figures 182 & 183) on which you intend to evaluate the software. Supply this to Andor and a license can be supplied for the evaluation period. When you receive the License from Andor, place it on the Desktop of the test computer and install Komet for testing.

Without the dongle your Komet software will only run in Demo Mode. In this mode Komet will not reveal GLP functionality, but otherwise will operate normally except: you cannot acquire images, open images, apart from the Demo images supplied by Andor; and you cannot save data.



Figure 182. To obtain the Computer Name, click on the Start Menu, select Computer and click right mouse Properties



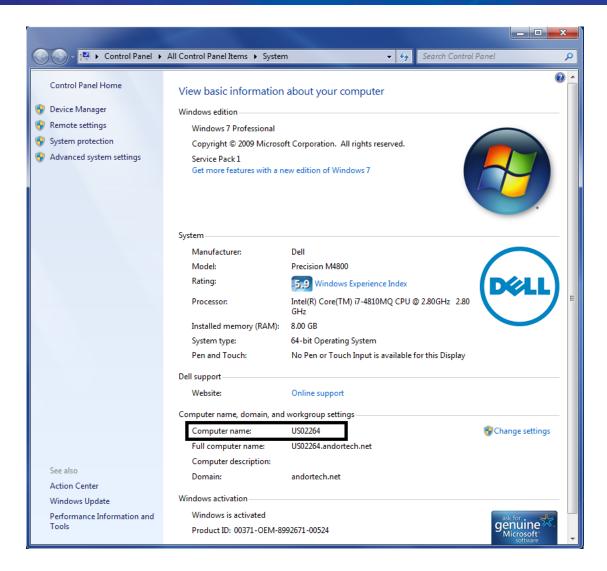


Figure 183. Obtain the Computer Name from the Computer Properties Dialogue as shown.

A1.2 Installation of Sentinel Drivers

If you have purchased Komet software or an upgrade license then you may download the installation files from the MyAndor Resource website. Before you install Komet and Database Viewer, we recommend you install the dongle or software protection key drivers. The dongle used for Komet is a Sentinel Software Protection Key and the drivers are available under **Software/Drivers and 3rd Party** as shown below. Download the file and run it to install the drivers, following the instructions provided.

.









A1.3 Initializing your Dongle with an Upgrade Code

If you have purchased Komet, the Software Protection Key you receive from Andor should have been initialized for your product. This section is aimed at users who have purchased an upgrade and already have a dongle.

There are two types of dongles. One is a parallel port dongle which is plugged into the LPT1 (printer port) and the other is a USB port dongle which can be plugged into any USB port. The LPT port dongle is now obsolete and the LPT port is not supported on modern PCs capable of running Windows 7 and later. Therefore you must obtain a new dongle as part of your upgrade. Please advise Andor if this is the case and arrange to return the old dongle.

To initialize your dongle, ensure that it is plugged into a USB port on your computer. Then follow the instructions below to install the Komet software. Once this is done Run the Configuration Manager and Open the **Software Protection** page from the menu bar. Copy and Paste into the Edit Box the 16 digit code provided and click **Initialize Your Dongle**

Your dongle will then be initialized.

A1.3 Installing your Komet Software

1) Assuming you have downloaded Komet or GLP from MyAndor: Open Windows Explorer and select the Downloads folder as shown in Figure 185. Double click on the Komet installation file (note the version number will be different from that shown). This will Run the installation.

Organize 👻 💼 Open 🛛 Share with		ew folder			=	-	
Favorites	 Name 	*	Date modified	Туре	Size		
🧮 Desktop	<mark>1</mark> 7z920-x64		9/17/2014 8:30 AM	Windows Installer	1,345 KB		
🗼 Downloads	SitplaneDe	emo80Oct14	9/25/2014 8:02 AM	LIC File	11 KB		
🖳 Recent Places	🔩 Imaris8_0_	0w64_beta0	10/6/2014 5:06 AM	Application	443,867 KB		
	1 Setup.X86.	en-us_O365ProPlusRetail_f	80c 9/30/2014 11:22 AM	Application	1,031 KB		
📜 Libraries	Setup_Kon	net_7.0.1.37	10/28/2014 12:34	Application	235,423 KB		
 Documents Music Pictures Videos 	E Support-L	ogMeInRescue	9/30/2014 11:16 AM	Application	1,493 KB		No pre availa
 Computer OS (C:) home (\\nc-dc-01) (U:) filestore (\\nc-dc-01) (Z:) 							

Figure 185. Installing from a file downloaded from MyAndor.

2) You will see this dialogue shown in Figure 186. Select Next...



Figure 186. Andor Installation Window (note that the appearance of this menu will vary with your software version).

.





To proceed with the Installation Read and indicate Acceptance of the Single User Software License Agreement shown in Figure 187.

2	Setup - Andor Komet
	License Agreement Please read the following important information before continuing.
	Please read the following License Agreement. You must accept the terms of this agreement before continuing with the installation.
	Andor Bioimaging SINGLE USER SOFTWARE LICENSE
	BY INSTALLING THE SOFTWARE YOU INDICATE YOUR ACCEPTANCE OF THESE TERMS.
	This Andor Bioimaging software (the 'Software') is copyrighted by Andor Technology. All rights are reserved. The PLIPCENASEP is grapted a LICENSE to use the coffware only.
	I accept the agreement
	< <u>B</u> ack Next > Cancel

Figure 187. Accept the EUL Agreement as shown to continue

Setup - Andor Komet	
Install Options	Kon
Select optional components for installation	
Install drivers for Andor sCMOS cameras (N	leo, Zyła)
Current selection requires at least 429.3 MB of o	lisk space.

Figure 188. Select the Andor sCMOS camera driver if you have chosen to buy a Zyla sCMOS camera with Komet.





Setup - Andor Komet	
Additional Tasks	Komet
Select the additional tasks you would like Setup to perform while installing then click Next.	Andor Komet,
Create Komet shortcuts:	
Create desktop shortcuts	
Create Start menu shortcuts	
< <u>Back</u>	Cancel

Figure 189. If you proceed with default selections then the Komet Installer will create desktop and Start menu shortcuts.

	Setup - Andor Komet
	Ready to Install Setup is now ready to begin installing Andor Komet on your computer.
	Click Install to continue with the installation, or click Back if you want to review or change any settings.
	Destination location: C:\Program Files (x86)\Andor Bioimaging
	Setup type: Full installation
	Additional tasks: Create Komet shortcuts: Create desktop shortcuts Create Start menu shortcuts
1	▼
	< <u>B</u> ack Install Cancel

Figure 190. Shows a summary before you proceed to installation.

3) The Komet Installer will ensure that all necessary steps are completed to install Andor Komet and Database Viewer (DBV) on the computer.



APPENDIX 2 HARDWARE DEVICES

A2.1 Camera Control Interfaces

Komet software supports the following cameras and acquisition devices.

Video CCD via Virtual Camera and SVid-USB2:

Many earlier Komet systems were sold with or to operate with video CCD cameras. These cameras have a video signal output delivered via BNC connector and cables. Komet formerly supported **Matrox Meteor 1 & Meteor 2 Controlled Frame Grabbers.** These frame grabbers are no longer supported. In order to provide backward compatibility for the many users wishing to upgrade to Komet 7.*n* we now support a very cost effective solution which provides an interface from the older Video source (Vid) to the modern USB 2.0 interface, known as Vid-USB interface. Andor supplies the Star Vid-USB adapter as an accessory. A photograph of the product is shown in Figure 191 below.



Figure 191. Star Video-USB interface (SVID2USB2) supported under Windows 7. A BNC adapter is supplied to install into the Yellow cable for composite video signals from the video CCD cameras.

The following video CCD cameras are supported:





- Hitachi KPM1 Monochrome PAL and NTFS Video (with integration option)
- JVC 1360BE Single Chip Colour Video
- Hamamatsu C5985
- Any other Camera providing an Analogue Video Signal.

Andor Luca S and LucaR CCD cameras



• Previous versions of Komet were sold with Luca cameras which offer excellent sensitivity and offered either 640x480 or 1kx1k pixels with 14 bit readout.

Note: Unfortunately these cameras are no longer manufactured due to obsolescence of the image sensors. However, they are still supported in Komet 7 and 7GLP.

Andor Zyla cameras



• The latest innovation in scientific CMOS cameras offers huge field of view - matching the microscope ocular. See the whole field and score more cells per image to speed through study scoring, Zyla USB plugs right into your computer's USB3 port and delivers extraordinary sensitivity and image quality.

Available in a Komet package at a highly competitive price.





QImaging Firewire Controlled Digital Camera

• QICam and similar camera such as Retiga.

The QICAM is a Firewire (IEE1394) camera that requires installation of an interface card into the computer as shown below.



Figure 192. The QImaging QICAM which provides a cost effective entry level digital CCD camera.

Any Camera running under Windows 7 can be supported.



Figure 193. Screen Capture Shortcut- Andor Virtual Capture

With Komet 7 we introduce the **Andor Virtual Camera (AVC)** which allows support of ANY camera with Komet using a screen capture technique. Any cameras capable of providing a live image window on a second monitor of the same computer can be used with Komet. The AVC is configured to acquire pixel data direct from the Live image window on the second monitor. These pixel data are used for analysis of comet images. The second monitor will cost typically around \$150 and provides a cost-effective solution for comet analysis while utilizing existing investment in hardware.

In fact, AVC also supports SVID2USB2 and other cameras that may have direct USB interfaces and are supported under the Video for Windows (VFW) or DirectShow capabilities of Windows multimedia.



A2.2 Virtual Camera and Analogue CCD Cameras

The Virtual Camera has been designed to allow users to capture images using cameras other than those currently provided by Andor, either indirectly by using screen capture (from a 3rd party application) or directly from many different types of live sources including network cameras, web cams, capture cards, TV tuners and dv cameras.

The camera provides access to capture from:

- any DirectShow or WDM compliant device
- capture cards, webcams; HD digital tuners or screen capture

	ol			1	
Video So	Integrated Webcam 💌				
Device	No Video		1		
Size	Integrated Webcam		<u></u>		
	LEAD Screen Capture (2.				
	Convert To Grey Leve	el 🔄			
v	Launch Help File on Startu	p			
		p			
	Launch Help File on Startu me Rate:	P			
		P			
		P	7		
	me Rate:		•		
			-		
	me Rate:				



When the software loads, the virtual camera will scan the system and list the available devices in the "Device" drop down control. Select the appropriate item to gain access to that device.



equisition/Auxiliary Devices	Acquisition/Auxiliary Devices
Image: State of the s	Image: Current Image: Current Image: Current Image: Current Image: Current Top
Camera	Camera
Image Control	Image Control
Video Source Device Integrated Webcam Size 640x480 Options Convert To Grey Level	Video Source Device Integrated Webcam ▼ Size 640x480 ▼ 160x120 ▼ 32x240 35x248 640x480 1230x720 1230x720 132x1080
I⊽ Launch Help File on Startup Frame Rate:	I Scuribeu I Scuribeu Fiame Rate:
Andor Virtual Camera	Andor Vitual Camera

Figure 195. Acquisition Auxiliary Menu

For devices which provide a discrete list of image sizes, then these will be listed in the "Size" drop down.

*	12			
Current	Live	Snap	Record	🗖 Stay on Top
mera				
Image Contr	ol]			
-	1			1
Video Sou	Irce LEAD Screen C			
Device	710x665		1	
Size	Options			
	Convert To (Greu Level		
	,			
_				
	Launch Help File	on Startup		
Fra	me Rate:			
T T				1
	Andor Virtua	l Camera		

Figure 196. Acquisition Auxiliary Options button

When "Screen Capture" is selected a single size will be listed, to adjust the capture window click the Options button.



Komet[®]

· · · · · · · · · · · · · · · · · · ·	indow	-Window se Finder tool	10000000	0x0070e44 CapTool (503, 397)-(542, 43;
Manual Capturing Manual Mode	apture Frames	Source out Outline Outline sha Color	source	orners
On source close:	Continue from last	position 💌	-	
On source minimize:	Continue from last	position 💌	·	
)n source area change:	Capture other win	dow 💌	-	
Capture Option:	Capture visible are	ea 🔄	-	

Figure 197. Capture Menu options

The dialog provides options for selection of the source window or specific region on the screen.

- The recommended option is to select a specific window.
- Then click and drag the "Finder tool" icon onto the window to be captured. As you move the mouse around the screen each window panel will highlight, this indicates that the window has focus.
- Release the mouse button to complete selection.

By selecting "Track window", you can automatically adjust the region to be captured. If the selected window is moved then the region will adjust appropriately. Note that size will not be adjusted.

The tracking feature is only available within a session, once iQ is closed then the window tracking will be lost. On startup, the software will remember the last position used but it cannot continue to track the window as it is moved. To re-enable tracking the window must be re-selected using the steps above.

Note: One important thing to remember is that the image view of the capture application should not be resized since this will invalidate any calibration. If a window is resized then the calibration process must be rerun. Tracking can be used so long as the window size remains constant. Also ensure that



the capture window is not occluded as this will be reflected in the screen captured image, the capture is WYSIWYG, What You See Is What You Get.

Note also that when capturing from video source, the virtual camera does not provide adequate control of the brightness and contrast. The dialog which contains these settings cannot be used while live, also settings do not get recorded for reuse in subsequent sessions. Instead it is recommended that the user adjust the settings using the Grabee software which is provided with the USB capture device. The settings will persist for the duration of the windows session, however the image should be adjusted each time Windows is restarted.

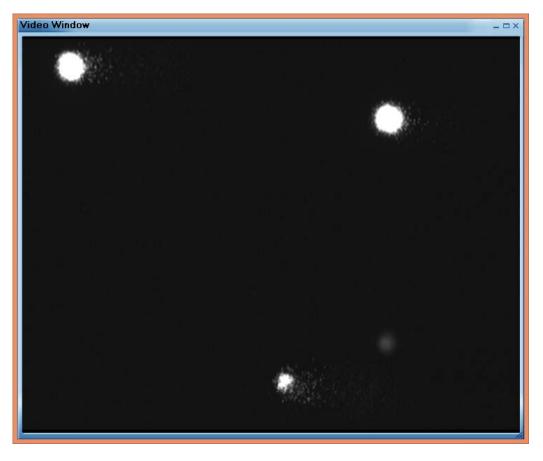


Figure 198. The Grabee Video window





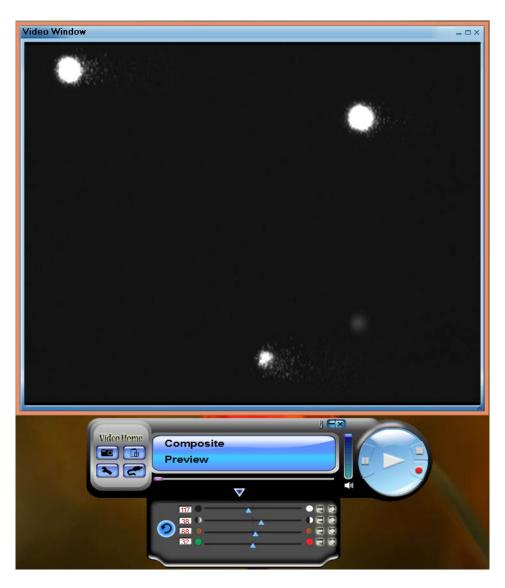


Figure 199. Grabee Video window with options



Figure 200. Grabee options for brightness and contrast adjustment

We recommend that you select a second monitor for your application space in which to show the live image window when using a third [arty camera with Virtual Camera screen capture mode.

.



A2.4 QImaging Cameras

A2.4.1 The QICam

Camera Specifications:

Air Cooled 12 bit digital CCD Camera.

Features:

Low Noise Low Dark Current Wide Dynamic Range Digital Output Super Pixel Readout (Binning) Progressive Scan Interline CCD Electronic Shutter Function

Hardware Requirements:

Firewire board with power directly connected to the Board. The Firewire board requires a free PCI slot on the computers main board and a Firewire interface cable for communication between the camera and the Firewire board in the computer.

Software Requirements:

Andor Imaging Software with the QImaging ActiveX Device Controller.






A2.4.2 The Device Tab

Selecting the Device Setup button from the Acquisition Tab opens the Acquisition /

Auxiliary Devices Window (Figure 201).

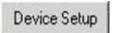

The Device Tab will automatically show the settings of the QImaging Camera connected as shown in Figure 201.

rice Camera		1
	QICamB	
12.000000000000000000000000000000000000		
	Monochrome (Cooled)	
EUTROLES .	1.68	
Firmware Version:		
Hardware Version:	÷	
	1392 × 1040	
Max Bit Depth:		
	8	
Max Binning:	4×4	
mand Status		Reset
mera is working		

Figure 201. The Device Tab.

Command Status

The Command Status section of the dialogue shows you if the camera is working (Figure 202).

Camera is	

Figure 202. Command Status.



Reset and About

The **Reset** button can be used at any time to restore the factory default settings. Selecting the Logo Icon below opens the About Dialogue Box (Figure 203). This provides information about the current software version installed on your system.



Figure 203. The Reset Button and About Dialogue Boxes.

A2.4.3 The Camera Tab

The camera setup is accessed from the **Device Setup** button on the Acquisition Tab (Figure 204).

Gain and Offset Control	
2x2 C 4x4 Gain: 1163 Offset 204	
Set to Factory Defaults	
€ 12 C 8 Exposure Time	
Target % Saturation	
C 100 C 50 C 75 C 25 Set Actual: 1ms	
mand Status	
mera is working	foca

Figure 204. The Camera Tab.





Sub Window

Sub-	Window
	Window
	656 : 494

Pressing the **Window** button with a region defined on the image in the Image Window, determines the specified X and Y dimensions of any image acquired when performing a Snap or running an Experiment (Region Defining Tools on the Image Toolbar).

Toggling the **Full Chip** button will revert the image size back to the full chip.

sub-	Window
	Full Chip
	154 : 136

Note: Full Chip image size depends upon the Window Mode and Binning values selected.

Binning

Binning	ANTING AS
C 1x1	C 4x4
	C 8x8

This is set by selecting the appropriate option button.

The default setting and the recommended setting is 2x2.

Changing the binning will alter the size of the Live image displayed in the Image

Window and also of images resulting from subsequent snaps and experiments, while still displaying the same field of view. In the above example the binning has been set to 2x2. This will result in an output image ½ the size of the original image in both the X and Y directions.

Binning is an operation to perform on-chip integration, trading spatial resolution against improved sensitivity. Charges from pixels in a rectangular neighbourhood are integrated to a single value in the output image, producing a smaller output image with proportionally improved sensitivity and reduced spatial resolution.

Smaller output images speed data transfer times, so acquisition speed also benefits.



8 Bit Acquisition

The Bit Selection should be automatically selected. Check the label on the camera to verify the bit selection.

Bit	Selecti	on
•	12	C 8

Auto Exposure

The **Set** button will calculate an exposure to give almost saturated pixels for the brightest part of the image. If a rectangular window is selected the exposure will be based on the region enclosed within the window. The camera attempts up to 5 iterations of exposure, usually achieving an acceptable exposure time. Occasionally manual intervention may be required or a second press of the **Set** button.

Target % S	aturation
€ 100 € 75	C 50 C 25
Set	

If you know your image is likely to get brighter during the course of an experiment you use the **Set** button to adjust the target saturation to 25, 50, 75 or 100 % to allow for an up to four fold increase in intensity before saturation.

Window Mode

This parameter changes the read-out on the camera chip. On performing a Snap or running an Experiment, the resultant image will be of the specified X and Y dimensions.

Wind	ow Mode	
۲	w 1392 : h 1040	
C	w 1024 : h 1024	

Warning - If the screen resolution of your Windows desktop is set to a lower value than the specified **Window Mode** parameter, you will not get an image displayed in the Image Window when in Live mode.



Gain/Offset Control

Gain:	1163	Offset: 2047
•	•	

On pressing the **Live** button on the Acquisition page, live video is displayed in the Image Window. The gain and offset of the live image can be adjusted using the **Gain** and **Offset** scroll buttons.

Note: It is recommended to keep the Gain and Offset control Set to the Factory Defaults. Use Auto **Exposure** at the desired saturation level to adjust the exposure time. If the gain and offset have to be adjusted manually, use the **Gain** control bar to increase sensitivity and the **Offset** control bar to bring up the black level.

Exposure Time

This is specified either in seconds, milliseconds or microseconds. The time each frame is exposed during the image acquisition is specified by typing a value into the text box then pressing **Enter** on your keyboard to accept the value.

0 m:s	_0 _1	÷
🕶 ms	1	- <u>-</u>
) us	0	-+

Using the Auto Exposure feature will automatically pick the best Exposure Time. The **Actual Exposure Time** displays the current exposure time in seconds.

.



APPENDIX 3 CONFIGURATION MANAGER

The **Configuration Manager** allows Komet to have Multiple User Levels. Three levels exist: Administrator, Power User and Restricted User. These levels have different status depending on whether the user is logged onto the Local Machine or a Network Domain (**Section 6**).

The Configuration Manager is only available to Administrators logged onto the Local Machine and NOT logged onto a DOMAIN.

A3.1 Run Configuration manager

To run the **Configuration Manager**, click on its icon in the Andor Folder from the Windows Start Menu. A Configuration Device Dialogue box will appear (Figure 205).

Configuration Wizard				
New	Edit	Delete	Exit	Current User- Carrie Low

Figure 205. Device Configuration Manager Dialogue Box



A3.2 Configuration Manager Menus

A3.2.1 Settings Menu

Select **Settings** from the Configuration Manager menu to set the Global Configuration Settings and to make changes to the Software Protection Key (Figure 206).



Figure 206. Settings Menu Expanded

Note: Enable Demo Version is not currently applicable to Komet.

A3.2.1.1 Settings Menu: Global Configuration Settings

When the Configuration Manager is first run there will be no current Configurations listed. You will first need to setup a configuration for your systems by selecting the **Global Configuration Settings** from the Settings menu (Figure 206). This allows you to select parameters that will affect User Access level control and Image Disk configuration.

A3.2.1.2 Multi-User Access

The **Multi-User Access** option is set by clicking in the checkbox (ON), or leaving the checkbox blank (OFF), in the Global Configurations Settings dialog (Figure 207). When **Multi-Level User Control** is **ON** the controls described will be applied to Users, according to the local machine status, decided by the Administrator. If you have **Multi-Level User Control OFF** all users have Full access.



bal Configuration Setti	inga:					
						Done
lulti-User Access						
When you select this teature, two These two levels are determine					P Multi-Level (Jser Control ON
Administrators:- FULL Access		Other Us	ers - Full Access I	to Wizerd. NO Access to Herd	ware Settings	
nage Disk Configuration						
ImageDisk Path						
Company and the second second second						
C Virtual Image Disk (Defau	ith Images ar	re NOT retained afte	er shut down (Reco	ommended for Komet Users).		
	100000		100000-20000-000	ommended for Komet Users). nded for iQ Users). An Imagel		n this drive.
Select a Drive if you want	Images to be	e retained after shut	down (Recommer		Disk folder will be created o	
Select a Drive if you want Important! An NTFS form	Images to be out hard disl	e retained after shut	down (Recommer	nded for iQ Users). An Imagel Windows Operating Syst	Disk folder will be created o	nance.
Select a Drive if you want Important! An NTFS form	Images to be nat hard disl 7 68 F	e retained after shut Ik drive which doe	down (Recomments NOT host the	nded for iQ Users). An Imagel	Disk folder will be created o	
Select a Drive if you want Important! An NTFS form Total Disk Space 9.77	Images to be out hard disl 7 G8 F	e retained after shut Ik drive which doe Free Disk Space:-	down (Recommer as NOT host the 5.72 GB	nded for iQ Users). An Imagel Windows Operating Syst	Disk folder will be created o	nance.
Select a Drive if you want Important! An NTFS form Total Disk Space- 3.77	Images to be nat hard disl 7 G8 F Ch user only h	e retained after shut Ik drive which doe Free Disk Space-	down (Recommer as NOT host the 5.72 GB wn Images.	nded for iQ Users). An Imagel Windows Operating Syst	Disk folder will be created o	nance.
Select a Drive if you want Important! An NTFS form Total Disk Space- 9.77 Private ImageDisk - Eac	Images to be nat hard disl 7 G8 F Ch user only h	e retained after shut Ik drive which doe Free Disk Space-	down (Recommer as NOT host the 5.72 GB wn Images.	nded for iQ Users). An Imagel Windows Operating Syst	Disk folder will be created o	nance.
Select a Drive if you want Important! An NTFS form Total Disk Space- 9.77 Private ImageDisk - Eac Shared ImageDisk - All	I Images to be at hard dist 7 GB F Sh user only hi users will hev	e retained after shut Ik drive which doe Free Disk Space- ias access to their ov re access to the sam	down (Recommer as NOT host the 5.72 GB wn Images.	nded for iQ Users). An Imagel Windows Operating Syst	Disk folder will be created o	nance.

Figure 207. Global Configuration Settings Dialogue Box

A3.2.1.3 Image Disk Configuration

The Image Disk Configuration (Figure 208) is accessed by selecting Global Configuration Settings item from the Settings Menu (Figure 206). As previously mentioned, all applications use a folder to store acquired image data in a RAW format. The Configuration Manager provides options for specifying the drive where this folder will be located. For best performance, we recommend a separate fast SCSI or EIDE drive, or even a RAID array for this purpose, but the Administrator can specify the drive. However, it must be accessible for full rights by ALL users when running Komet, otherwise, restricted access users will have access only to the Windows virtual memory system; the contents of which is lost whenever Komet closes down.

.



There are two options for choosing the **Image Disk Path**: Virtual Image Disk (Default setting) and Select a Drive (Figure 208).

- Virtual Image Disk A Virtual Image Disk can be selected which will force Andor to use Windows Virtual Memory System to store images. Using the Virtual Memory system will mean that ALL images on the ImageDisk are lost at power down. In contrast images stored on a "Real" Disk Folder will be retained between sessions of Komet unless deleted by the users.
- Select a Drive The path to the ImageDisk folder is configured here, when the drive is selected in the drop down box. The free disk space of the partition selected is displayed to help select an appropriate place for this folder.

Note: We strongly recommend NOT using the System Partition for ImageDisk as this can cause system operation problems when full.

In addition, the **Private Image Disk** and the **Shared Image Disk** options are applicable to the Virtual Image Disk or the Select a Drive option. However, if Virtual Image Disk is selected, regardless of whether you have chosen the Private Image Disk option or the Shared Image Disk option, all images will be lost at power down.

- Private ImageDiskThe Image Disk folder will be created on the specified drive. Each WindowsLogon User will have a folder for their own images, thereby ensuring
isolation between different users.
- Shared ImageDisk The Image Disk folder will be created on the specified drive and then All Windows Logon Users will access a shared folder, thus allowing All Komet users to share images.

Images in RAM specifies the number of frames that Komet will attempt to keep in RAM at any one time. Some image processing and visualization operations, e.g. 3D

View, ClearView & Orthogonal View in other products, may require all frames to be loaded simultaneously; therefore this setting must exceed the total number of frames being loaded. The Default Value setting is 65536.



A3.2.2 Settings Menu: Software Protection Device (Dongle)

The **Software Protection Device** enables the Komet software to run on your system. The configurations for your system have been initialized on the Dongle. There is no need to change these configurations unless new software has been purchased, at which time you will be provided with instructions, so changes can be made accordingly.

A3.2.3 File and Edit Menus

Hardware Configurations can be Created (New), Edited and Deleted. These functions are available through the Buttons on the Configuration Wizard interface or by selecting the options from the File or Edit Menus (Figures 208).

Configuration Wizard				
New	Edit	Delete		Exit
🐞 Device Configurat	ion Manager	🖀 Devi	ce Configuration M	anager
File Edit Settings Too	ls About Exit	File Edi	Settings Tools Ab	out Exit
View Configurations	ard		Edit Configuration Delete Configuration	
New Configuration Open Configuration	1		1	
	Edit		New	Edi

Figure 208. Configuration Wizard buttons, File and Edit Menus Expanded



A3.2.4 Tools Menu

The **Tools** menu is used for registering/unregistering Any ActiveX Component, registering/unregistering A Plugin Device and troubleshooting hardware devices through the File List Generator (Figure 209).

File Edit Settings	Tools	About Exit		
Configuration		jister / Unregister Any ActiveX Component jister / Unregister A Kinetic Plugin Device		
Nev	File List Generator			

Figure 209. A3.5 Tools Menu Expanded

A3.2.4.1 Tools Menu: Register/Unregister Any ActiveX Component

Select **Register/Unregister Any ActiveX Component** from the Tools menu item (Figure 209). Select the file to be registered / unregistered by defining the correct file path in the right hand panel and highlighting the appropriate file name in the left hand panel (Figure 210). The files displayed are determined by the file extension entered into the **Filter** text box. Files are registered or unregistered by clicking on the **Register** or **Unregister** buttons, respectively.

Register Done Done	MPORTANTI Eyour device driver fails to register check that any Support Files and Drivers have been installed
TANAuto on TaNaurContole on	Filter •ocx

Figure 210. Register/Unregister ActiveX Components Dialogue Box



Note: The '**regsvr32.exe'** file must reside in the same folder as the ActiveX component for the registration process to work correctly.

Select Done to exit from the Register/Unregister Any ActiveX Component menu.

Note: If the Configuration Manager finds ActiveX files with the same class name in your Andor Bioimaging folder the following dialogue in Figure 211 will be displayed.

				Se Th	ect the fix button older files will b	files found on your system. to rectify the problem . e moved to another folder lew) files will be registered.
Problem Auxiliary Devices:- Kinetic Stage Device		-			Fix1	Done
uplicate Files:-	Internal Name	Date	Time	Size		
(DYZController.ocx (DYZController2.ocx	Kinetic\KIXYZController	13/03/03 13/03/03	23:37:18 23:37:18	729.5 729.5		
ther Files Associated With						

Figure 211. Breakdown Recovery Dialogue Box

Selecting one of the problem devices that are listed will result in the duplicate files being displayed in the Duplicate Files list box. To resolve the problem press the Fix button and the older of the files will be moved to an '**OldDeviceFiles**' folder. The remaining file will then be registered.

.



A3.2.4.2 Tools Menu: Register/Unregister Plugin Components

Select **Register/Unregister Plugin Components** from the Tools menu item (Figure 212). Select the file to be registered / unregistered by defining the correct file path in the right hand panel and highlighting the appropriate file name in the left hand panel (Figure 212). The default folder for Plugin software modules is the 'Andor Bioimaging\Common File' folder. Files are registered or unregistered by clicking on the **Register** or **Unregister** buttons, respectively.

Register	Unregister	Done	IMPORITANTI If your device driv Drivers have been installed.	ver fails to register check that any Support Files and
KISOViewPlugin dl KISet2Plugin dl KICsopPlugin dl KIPperPlugin dl KIPperPlugin dl KIPperPlugin dl KISmogaPlugin dl KISsipPlugin dl KISsipVewPlugin dl KISsipVewPlugin dl				C C Phogan Files Configurator File/Active Configurator File/Active Configurator File/Active Images

Figure 212. Register/Unregister Plugin Components Dialogue Box

Select **Done** to exit from the **Register/Unregister Plugin Components** menu.



A3.2.4.3 Tools Menu: File List Generator

The **File List Generator** is a tool to help Andor Bio-imaging Division to resolve any problems you may have when updating from an older version of Komet or other

Andor product. The File List generated by this function can be sent to Andor Bio-imaging Division and lists the date and version number of every file used in running the application. To generate a suitable file list select the folder C:\ Program Files\Andor Bioimaging\Common Files and follow the instructions listed in the dialogue box (Figure 213). When requested by Andor Bio-imaging Division, send the resultant file to the technical support team along with an explanation of your problem.

e List Generator Open a Kinetic Fi Select Target Dri			_		Troublashoofing Visatio	File Version Problems.	e
C) Program Files Andor Bioimagi ConfigurationF Graphics Images	ing	STOP			To test for Kinetic Fil	from where to start your search. Ie Dates just search in the Kinetic/Common Files folder. copies of files select the Drive. (This could take some time	1)
Filename	and the second se	the second s	File Date	File Time	Error Report	Full Path	Γ

Figure 213. File List Generator Dialogue Box

.



A3.2.5 About Menu

Selecting **About** from the Configuration Manager brings up a dialogue box which contains information about Andor Bio-imaging Division (Figures 214). Press OK to close the dialogue box.

🚆 Device Configuration Manager		
<u>File E</u> dit <u>S</u> ettings <u>T</u> ools <u>A</u> bout <u>E</u> xit		
Configuration Wizard		
New Ec		Bioimaging Division
		www.andor.com
Current Configurations		www.undoi.com
vcam	Corporate Headquarters	Andor Technology USA
	7 Millenium Way	425 Sullivan Avenue
	Springvale Industrial Park Belfast BT21 7AL	Suite 3 South Windsor
apz	United Kingdom	CT 06074, USA
	Tel: +44 (0)28 9023 7126 Fax: +44 (0)28 9031 0792	Tel: +1 (860) 290-9211 Fax: +1 (860) 290-9566
	row_microscopy_support@andor.com	row_microscopy_support@andor.com
	Andor Technology Japan 4F NE Sarugakucho Building 2-7-6 Sarugaku-Cho, Chiyoda-Ku Tokyo 101-0064 Tel: +81 (3) 3518 6488 Tel: +81 (3) 3518 6489 apac_microscopy_support@andor.com	Andor Technology China Room 1213, Building B Luo Ke Time Square, No.103 Huizhongli, Beijing 100101 Tel: +86 (10) 5129 4977 Fax: +86 (10) 6445 5401 apac_microscopy_support@andor.com
	Copyright Andor Technolog	y 2006 - 2014
	ОК	

Figure 214. About Menu Expanded and the About Dialogue Box

A3.2.6 Exit Menu

Select Exit to close the Configuration Manager (Figure 215).



Figure 215. Exit Menu Expanded

A3.3 Create a New Configuration

Select **New Configuration** from the Configuration Manager File menu or select the **New** button on the Configuration Wizard interface (Figure 216).

ile Edit Settings Tools About Exit View Configurations New Configuration	Device Configuration	n Manager			
New Configuration	ile Edit Settings Tools	About Exit			
New Configuration	View Configurations		•		
		(New	Edit	Delete	Exit

Figure 216. File Menu Expanded and Wizard Buttons

The first step in setting up a New Configuration is to enter the Configuration Name in the dialogue box (Figure 217). This should reflect the devices that are to be included in the configuration set up. Press **Next** >> to continue.

Next >> Cancel Step 1 2 Enter a Name for the Configuration. It is recommended that the Name of the Configuration should give some indication of the Devices being used. It is recommended that the Name of the Configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name It is a configuration should give some indication of the Devices being used. Infiguration Name					
It is recommended that the Name of the Configuration should give some indication of the Devices being used. Integration Name Non Name 1 True	14				
onfiguation Name (Xon Jate & Time	na l	ve some indicate	ve Configuration should or		
Dicon ale 1 Texe	6 0 .	te sonne interdio	e comgaranon shouo g	o narie narie or ne	II IS I CONTINUED DE
Dicon ale 1 Texe					
Dicon ale 1 Texe					
l late & Time					onliguation Name
					iXon .
					anes.co
Current Date - 6/9/2005 8:53/23 AM					
				/9/2005 8:53:23 AM	Current Date -





The second step is to select the camera that is going to be used for your system (Figure 218). Click the required camera device and Press **Next** >> to Continue.

Select a Camera from the list.	p1234
Select the Dummy Camera Device if you do not have a Camera attached to your system.	
Ivvalable Camera Devices	
Andor Camera With fast Piezo Z and Monochromator Andor Comma	
Hamamatsu Orca I/ER 4742-95 Finewine Kinetic Dummy Camera	
Parents Contrary Contrarts	

Figure 218. Select the Camera Dialogue Box

The third step is to select the auxiliary devices that are to be included in the configuration (Figure

219).

	Second
<c back="" next="">> Cancel</c>	Step 1 2 3 4
elect the devices from the list that you wish to use in this configuration.	
sildshe Ausling Devices	
Zeine Aviocitati Device	
Kinetic Aktuator Device	
ASI Stege Device	
ASI Filter Wheel Device	
Atto KIAtto AIC Device	
Cam Hiter Wheel Device	
Conix ZMotor Device	
CRI MicroColor Device	
Leice DMR Device	
Nikon E1000 Device	
Olympus M81/BM81 Device	
Lud Stage Device	
Ludi Filter Wheel Device	
Physik Instrumente LVP2T Device	
Merzhauser Stage Device	

Figure 219. Select the Auxiliary Devices Dialogue Box



The devices are selected by enabling the appropriate check boxes. To continue Press Next >>.

The final step is to confirm the New Configuration setup. A summary of the selected camera and auxiliary devices is displayed (Figure 220). At this point the user can edit the device selection, if necessary, by returning to the previous pages in the Setup Wizard, using the **<< Back** button. Once the appropriate camera and auxiliary devices have been chosen, click the **Finish >>** button to complete the device activation procedure.

ee Back	Finish >>	Cancel	Step 1 2 3 4
	- Annonitation 		Step 12.54
Configuration Summary - Namie :-	Intern		
Name."	ixon		
Camera Devic	e:- Andor Ca	mera	
Auxiliary Devic	nes - L		
Traslicshooting			
And the second second second	d Aubiliary Device	s listed above are Registered when t	nishing the Configuration Wizard

Figure 220. Configuration Summary Dialogue Box

The **Show REGSVR32.EXE** checkbox at the bottom of the dialogue is enabled by default. When the **Finish** >> button is pressed the devices will be registered and a series of message boxes displayed indicating whether the devices have been successfully registered. These messages can be inactivated by disabling the **Show REGSVR32.EXE** checkbox. We recommend leaving the box checked to make sure all devices register correctly.

Note: To exit the Configuration Wizard at any time, ignoring changes that have been made, click the Cancel button.



A3.4 Open Configuration

Select **Open Configuration** from the File menu of the Configuration Manager (Figure 221).

Device Configuration Manager							
File	Edit	Settings	Tools	About	Exit		
Vi	iew Co	nfiguration	is e	rd			
New Configuration			1				
0	pen C	onfiguratio	n				
1							

Figure 221. Open Configuration Menu Option

The Configuration Manager automatically saves a backup version of the configuration file to an Archive folder, every time it is closed. This ensures that no configuration data is lost and allows previous configurations to be recovered, should the current configuration file become corrupted or accidentally deleted.

A catalogue of archived configuration files can be found in the **Old Configurations** list on the **Manage Configuration Archive** (Figure 222). The details of each configuration file are displayed in a Tree View structure in the right hand panel. Files are selected by clicking on the appropriate file name using the left mouse button. The Tree View is expanded by clicking on the "+" icon.

The Make a Backup Configuration File each time the Configuration Manager is Closed checkbox, shown at the top of the dialogue, remains active by default. This ensures that backup copies of configuration files are created each time the Configuration Manager is closed. The archive facility may be deactivated by disabling the checkbox.



Rewisd	Delete	Done	Current Configuration Walks a Backup Configuration File each time the Configuration Manager is Dored.
d Configuration			
te the control key to m	uli-select.	-	
kicenig 218.000 32 kicenig 218.000 33 kicenig 318.000 33 kicenig 48 3000 33 kicenig 48 3000 43 kicenig 48 3000 43 kicenig 48 3000 43 kicenig 48 3000 53 kicenig 43 3000 53 kicenig 53 3000 54 kicenig 54 3000 54 kiceni	- 450 PAN UKA - 150 AAN UKA - 141 PAN UKA - 535 PAN UKA - 535 PAN UKA - 535 PAN UKA - 535 PAN UKA - 544 AAN UKA - 544 AAN UKA - 460 PAN UKA - 460 PAN UKA - 460 PAN UKA - 150 PAN - 150 PA		

Figure 222. Manage Configuration Archive Dialogue Box

Unwanted configuration files can be removed from the **Old Configurations** listing by selecting a file and then pressing the **Delete** button.

Selecting the appropriate file and clicking the **Rewind** button activates a previously archived device configuration file.

Clicking on the **Done** button exits the **Manage Configuration Archive** dialogue.



Komet[®]

A3.5.1 Edit an Existing Configuration

Select **Edit Configuration** from the Configuration Manager Edit menu or select the **Edit** button on the Configuration Wizard interface (Figures 223).

ta d	evice Configuration Ma	anager			
File	Edit Settings Tools Abo	out Exit			
C	Edit Configuration Delete Configuration	Configuration Wizard			
T		New	Edit	Delete	Exit

Figure 223. Edit Menu Expanded and Wizard Buttons

Select the appropriate configuration file name from the **Current Configurations** list displayed in the Configuration Wizard (Figure 224). A Wizard dialogue is displayed, similar to that used during the creation of new device configuration files (**Section 2.8.8**). The edit function allows modification of the configuration name and active auxiliary and camera devices

New	Edit	Delete	Exit	Current User - Carrie Lowe	
rent Configurations					
Dummy (ionera				
X	hokromator				
FLZ					
- Won					
	rated By- Carrie Lowe				
	te Created - 2/16/2005				
	te Last Modified - 6/9/				
-	quisition Device:- And				
	Driver File:-KIDAn	dorRTAcox			

Figure 224. Current Configurations List



The available configurations are listed in a Tree View, each comprising the following information (Figure 224):

Created By	The name of the Administrator who was logged onto
	the Local Machine when the Configuration was created.
Date Created	The date the Configuration was created.
Date Last Modified	The date the configuration was last modified.
Acquisition Device	The Camera Device used in the configuration.
	The Camera Device Driver (OCX) used by the Camera
	Device can be seen by expanding this part of the tree.
Auxiliary Devices	The Auxiliary devices that are contained in the
	configuration (if any) are listed.
Auxiliary OCX Files	The Auxiliary Devices Drivers (OCXs) used by the
	Auxiliary Devices can be seen by expanding this part of
	the tree. This is useful for troubleshooting.



A3.5.2 Deleting an Existing Configuration

Select **Delete Configuration** from the Configuration Manager Edit menu or select the **Delete button** on the Configuration Wizard interface (Figures 225).

D	evice	e Configu	uration	Mana	ger			
ile	Edit	Settings	Tools	About	Exit			
0		lit Configur	ation		Configuration Wizard			
1	De	alete Confi	guration		New	Edit	Delete	Exit
				1		Lun	Delete	

Figure 225. Edit Menu Expanded and Wizard Buttons

Select the appropriate configuration file name from the **Current Configurations** list displayed in the Configuration Wizard (Figure 225). Before the selected configuration is removed from the **Current Configurations** list, the user will be required to confirm that they want to delete the selected file. The selected file name to be deleted will be displayed within the << >> brackets (Figure 226).

Press the **Yes** button to delete the configuration. Clicking the **No** button cancels out of the file deletion process.



Figure 226. Delete Configuration Dialogue Box

-



A3.6 Configuration Manager Summary

Once the Configuration Manager has been run to create a new setup, the Administrator will have to run Komet to set up all new devices used in that configuration. Hardware settings for the devices include COM ports, board selections and set-up, calibration information, filter locations and descriptions, etc. If you are upgrading from a previous version of the software, then using the upgrade utility will retain device settings and make these usable by the newer version of the software.



APPENDIX 4 KOMET MEASUREMENT DEFINITIONS

A4.1 Integrated Intensity Profiles and their Analysis

Komet analysis software is based upon the principle of integrated intensity profile analysis (Figure 227). This means that we integrate in the vertical direction (perpendicular to the direction of electrophoresis), the image intensities within the Cell and Background ROIs and work on these profiles to compute the comet parameters. The one exception is physical area, which is computed from the image data directly. We will discuss the algorithms with reference to the Density profile graphics shown below. The derived parameter definitions are listed in the following sections.

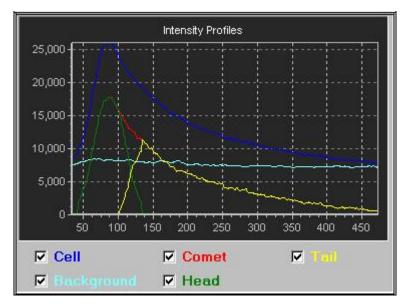


Figure 227. Intensity profile curves as presented in Komet

The profiles are computed and analysed as outlined in the following algorithmic description.

Footnote for Komet 3 Users

In Komet 4.0 & 5.0, the analysis has been integrated and both ver20 and ver22 forms of Head and Tail analysis are computed. We now refer to the ver20 Tail Moment as Olive Tail Moment, to reflect the first author of the original paper in which it was first defined.



1. Integration - Cell and Background ROIs are first integrated vertically and stored to double floating point precision, to avoid computational noise as far as possible.

2. Physical Area - based on the threshold value, T (see below) and the mean background intensity profile (MBIP), the physical area is computed by subtracting the MBIP from each line of the Cell ROI. For every pixel in the result which is greater than the threshold, T the Physical Area is incremented. The result is then scaled in X and Y by the current system scale factors, resulting from calibration.

3. Background Scaling -The background is next scaled and filtered (running average) by a factor F, where F = ((CellROIheight)/ (BackgroundROIheight)).

4. Background Offset - The background is then offset by an amount which depends on T t, the Tail Threshold setting in Comet Options, such that:

NewBack = OldBack + (T t *MaxCell/1000) where MaxCell is (max (CellProfile intensity) - max

(BackgroundProfile)).

5. Filtering - Cell and Background profiles are then filtered or smoothed by a running average filter with a range, R = (1+2*S), where S is the smoothing factor setting in Comet Options - a larger value means more smoothing.

6. Profile Subtraction - The Background profile is then subtracted from the cell profile, to create the difference profile, which is used for the comet head and tail computations.

7. Difference Profile Analysis - The difference profile is analysed for i) comet integrated intensity, ii) centre of gravity (mean), iii) maximum (mode), iv) standard deviation, v) skewness, vi) distributed moment and vii) total length. These are defined in the next section. Breaks (zero values) in the profile do not terminate analysis, unless there is a contiguous run of breaks which exceed the TailBreakLength setting in Comet Options.

8. Fit Head Profile to Difference Profile - Komet then attempts to fit a head profile to the difference profile data. This is achieved, by finding the peak intensity or mode, moving from the left-most extent of the ROI limits. The start of the head is identified as the first profile point which is above the Head Threshold setting in Comet Options. The mode is used as a fulcrum point to identify the head extent and to create the right hand side of the Head profile.

Komet[®]



If the **Head Symmetry** checkbox in Comet Options is checked then the Head profile is made symmetrical (if possible) and the data to the right of the Head Centre (Comet Mode) will be set to the minimum of either the left hand data value or the actual data. Selecting the minimum means that the profile will be symmetrical where physically possible, but will not over-estimate the head intensity.

If **Head Symmetry** is not checked then the data to the right of the Head centre will be set to the actual values over the range.

This means that if there is an apparent skew of DNA (intensity) towards the tail which you wish to detect as tail then you should select Head Symmetry.

On the other hand if you wish to discount the skewed intensity distribution then do not select Head Symmetry. Skewed Head intensity will not be included in the tail and this strategy may be useful in minimising false positive scores.

Once identified, the new head profile is used to compute i) head integrated intensity, ii) centre of gravity (mean), iii) maximum (mode), iv) standard deviation, v) skewness, vi) distributed moment and vii) head length.

9. Produce Tail Profile - The head profile is subtracted from the difference profile to produce the tail profile.

10. Tail Distribution Analysis - The tail distribution is then analysed for i) tail integrated intensity, ii) centre of gravity (mean), iii) maximum (mode), iv) standard deviation, v) skewness, vi) distributed moment and vii) tail length. Breaks (zero values) in the profile do not terminate analysis, unless there is a contiguous run of breaks which exceed the TailBreakLength setting in Comet Options.

11. Remaining Parameters Computed - The other derived parameters are then computed as outlined below.

.



A4.2 Komet Analysis - A Mathematical Description

Another way of viewing the analysis performed by Komet is to describe the process mathematically, rather than by algorithm. This approach may lead to a clearer understanding of the underlying process. We have restricted the description to the salient points without trying to include all of the algorithmic detail in mathematical notation.

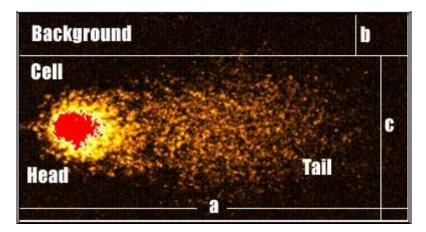


Figure 228. A typical damaged comet cell showing analysis Cell and Background ROIs

The image in Figure 228 shows a cell region of interest (ROI) and the comet cell clearly defined within it as well as a background ROI created directly above the cell. The extent of the cell ROI is a pixel in the x direction and c pixels in y and the extent of the background ROI is a pixels in x and b pixels in the y direction, as shown.

We define the Cell image I c as a function of i j, where i and j are pixel indices in x and y respectively -

 I_{C} (i, j) and the Background image as I_{b} (i, j). From this everything else follows the description below, which can be compared to the algorithmic description for those interested.

1) First we compute the intensity profiles, P $_{\rm C}$ for cell and P $_{\rm b}$ for background, by integrating in the y direction over the ROIs as follows:

$$P_{c}(i) = \sum_{j} I_{c}(i, j) i = 1 \dots a; j = 1 \dots c and P_{b}(i)$$

= $\sum_{j} I_{b}(i, j) i = 1 \dots a; j = 1 \dots b$



2) We then compute the background corrected profile P $_{bc}$ (i)

$$P_{bc}(i) = P_{c}(i) - (P_{b}(i) . c/b)$$
 $i = 1 a$

3) The Comet profile is then computed from this by applying the Tail threshold, T_t, which is specified in parts per thousand:

C (i) = P $_{bc}$ (i) - (max (P $_{bc}$ (i)).T $_{t}$ /1000) i = 1 a

A4.2.1 Head and Tail Discrimination

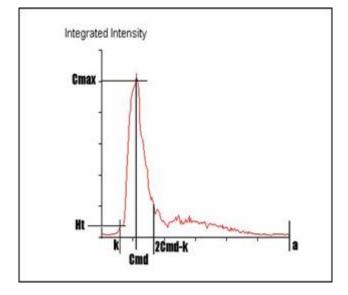


Figure 229. A background corrected comet profile with key features annotated

Discrimination of the comet Head and Tail proceed after the corrected comet profile C(i) has been computed (Figure 229). However, it must be stressed that this discrimination can be reliably undertaken only if the comet cell shows a certain "normal" morphology. This is true whether the analysis proceeds by conventional 2D image analysis or by intensity profile analysis. In Komet we require that the damage level is low enough such that the head region exhibits the brightest part of the comet intensity profile. In severely damaged cells, or necrotic or apoptotic cells this may not be the case and the user will generally over-ride the automatic analysis and define the Tail region with the mouse, as documented elsewhere in this guide.



However, assuming that analysis can proceed normally, we have included the option of "Head Symmetry" to achieve maximum flexibility in analysis. To aid in visualisation of this process refer to the intensity profile in the figure above.

The system creates the head intensity profile, as follows:

1) Detect start of head profile pixel, k

where first is a function which detects the first occurrence of the condition, H $_{\rm t}$ is Head threshold value and a is the cell ROI extent in x as defined previously.

2) If the user has selected *Head Symmetry* in Comet Options then:

where the min function takes the minimum of the two

parameters.

Note that when the Comet profile to the right of the Comet mode shows lower values than the equivalent values to the left of the mode, the min function ensures that Head intensity profile is not over-estimated.

3) If the user has **not** selected *Head Symmetry* in Comet Options then:

H(i) = C(i); $i = k \dots 2Cmd-k$, where Cmd is comet mode



Once the Head profile H(i) has been computed, then computing the Tail profile, T(i) is simply a matter of subtracting the Head profile from the Comet profile as defined here:

T(i) = C (i) - H(i); i = k a , where a is ROI limit T(i) = 0; i = 1 ... k-1.

A4.2.2 Ensuring Intensity Profile Equivalence

To ensure consistency of intensity analysis between the three resultant profiles we finally reassign the comet profile to be equal to the sum of the Head and Tail profiles. This rejects the intensity profile at the front of the comet, which is considered below the Head threshold and ensures that the detected Head and Tail intensity profiles combine to be exactly equivalent to the discriminated comet intensity profile. Thus we have:

Comet, Head and Tail parameters are then computed from the intensity profile as described in Section A4.4.



A4.3 Komet Measurements and Parameter Definition

It is important to differentiate between derived Global Comet parameters, in which we generally use the definitions first outlined by Olive and co-workers, and the individual component measurements on head, tail and comet. The first are global descriptions of the data, while the latter are specific measurements of the raw data. Hence we have adopted the notation which indicates these differences i.e. Tail Length and Tail Moment are now used to describe the Comet global parameters, while Tail Extent and Tail Distributed Moment are used for the component intensity profile parameters.

We have also implemented a number of new measures which have proven to be used in the literature. In the current version of Komet we have harmonised the calculations and added calculation of the inertia as defined by Hellman (1995) and the Comet Moment has been recently redefined by Kent *et al.* (1995).

.



A4.4 Comet Parameters

These are computed from the data in C (i), the Comet intensity profile, using parameters defined in this appendix.

Preliminary definitions - Optical Intensity - or population $P = \sum C(i)$.

Comet.Mode = Mode*IMAGE.Xres. Mode is the position of the peak intensity in the distribution, Cmd as defined previously in this Appendix.

Comet.Mean, Cm =

Profile Centre of gravity = $(\sum (C(i)*i)/P)$ *IMAGE.Xres.

Comet.StdDev = $((\sum (C(i)*i^2)/(P-1)) - ((P*Cm^2)/(P-1))) * (IMAGE.Xres).$

Comet.Extent = (RightEdge- LeftEdge)*IMAGE.Xres,

where RightEdge and LeftEdge are found by intensity searching,								
controlled	by	Tail	and	Head	Thresholds.	In	Figure	229,
a=RightEdg	a=RightEdge and k=LeftEdge.							

Comet.Distributed Moment	= (∑(C(i)*(i-k))/∑C(i)) * IMAGE.Xres	i = k a
	where k is defined in Figure A3.3	
Comet.Inertia	= (Σ(C(i)*((i-k) * IMAGE.Xres) ²)/ΣC(i))	i = k a
Comet.Skew	= ∑(((i-k)-Cm)³*C(i))	i = k a
Comet.CoefVar	=(Comet.StdDev/Cm)*100.0. CoefVar is the profile Coefficient of Variation	n.
Comet.Optical Intensity	= ∑(C(i)).	



A4.5 Head Parameters

These are computed from H(i), the Head intensity profile.

Preliminary definitions - Optical Intensity = population $P = \sum H(i)$.

Head.Mode = Mode *IMAGE.Xres.

Mode is the position of the peak intensity in the distribution, which is equal to Cmd in a "normal" comet.

Head.Mean, Hm =

Profile Centre of gravity = $(\Sigma(H(i)*i)/P)$ *IMAGE.Xres.

Head.StdDev = $((\sum (H(i)*i^2)/(P-1)) - ((P*Hm^2)/(P-1))) * (IMAGE.Xres).$

Head.Extent = (RightEdge-LeftEdge)* IMAGE.Xres

where RightEdge and LeftEdge are found by intensity searching and controlled by Head Threshold. In Figure 229, RightEdge=(2Cmd-k) and LeftEdge=k

Head.Distributed Moment	= (Σ(H(i)*(i-k))/ΣH(i)) * IMAGE.Xres	i = k a
	where k is defined in Figure A3.3.	
Head.Inertia	= (Σ(H(i)*((i-k) * IMAGE.Xres) ²)/ΣH(i))	i = k a
Head.Skew	= ∑(((i-k)-Hm)³*H(i))	i = k a
Head.CoefVar	= (Head.StdDev/Hm)*100.0.	
С	coefVar is the profile Coefficient of Variation.	
Head.Optical Intensity	= ∑(H(i)).	



A4.6 Tail Parameters

These are computed from T(i), the Tail intensity profile.

```
Preliminary definitions - Integrated Intensity = population P = \sum T(i).
```

Tail.Mode = Mode * IMAGE.Xres.

Mode is the position of the peak intensity in the tail distribution.

Tail.Mean, Tm =

Profile Centre of gravity = $(\Sigma(T(i)*i)/P) * IMAGE.Xres.$

Tail.StdDev = $((\sum (T(i)*i^2)/(P-1)) - ((P*Tm^2)/(P-1))) * (IMAGE.Xres).$

Tail.Extent = (RightEdge-LeftEdge)* IMAGE.Xres

where RightEdge and LeftEdge are found by intensity searching and controlled by Tail Threshold. In Figure 229, RightEdge=a and LeftEdge=(2Cmd-k)+1.

```
Tail. Distributed Moment = (\sum (T(i)^*(i-k)) / \sum T(i))^* IMAGE.Xres i = k \dots a
```

where k is defined in Figure A3.3.

Tail.Inertia	= (∑(T(i)*((i-k) * IMAGE.Xres) ²)/∑T(i))	i = k a
--------------	--	---------

Tail.Skew = $\sum(((i-k)-Tm)^{3*}T(i))$ i = k a

Tail.CoefVar = (Tail.StdDev/Tm)*100.0.

CoefVar is the profile Coefficient of Variation.

Tail.Optical Intensity $= \sum (T(i)).$

.



A4.7 Global Comet Parameters

Results will vary depending on the selection of Comet Options, and particularly with selection/deselection of the Head Symmetry option.

Head%DNA = (Head.OptInten /(Head.OptInten + Tail.OptInten))*100. Tail%DNA = 100 - Head%DNA.

Comet.OptInten = Head.OptInten + Tail.OptInten - where OptInten is the Optical Intensity.

Tail Length = Tail Extent.

(TailfromCenter) Tail Extent + Head Extent/2

Tail/Head (LoH) = Tail.Extent / Head.Extent.

Olive Tail Moment=(Tail.mean - Head.mean)*Tail%DNA/100. Extent Tail Moment=Tail Length * Tail%DNA/100.



A4.8 Conclusions

The data produced by Komet are extensive and fall into two main categories. On the one hand we have the derived parameters for Global Comet description, which are most commonly used for data analysis and comparison with other systems. On the other hand we have the pure geometric/densitometric measurements on all of the component profiles, which provide a detailed quantitative description of the profiles and from which the interested user can derive many powerful methods of Comet evaluation. The distributed moment is a case in point, which supplies a profile description which weights intensities by distance from the leading edge of the profile. The final result is normalised by the integrated intensity to reduce the effects of staining density. While we will continue to improve the analysis and handling of data, the raw Komet results produced are unlikely to change for the foreseeable future.



APPENDIX 5 KOMET REPEATABILITY MEASUREMENTS

A5.1 Measurements Made

Two different sources of image were used in the repeatability tests:

A previously stored Comet data file image was used as a standard image to check the reproducibility of making the same measurements and to investigate the effect of varying the size of the interactive box containing the Comet.

Live grabbing of a Comet image. This was used to test the reproducibility of results when the same image was grabbed between each measurement and to test the impact of choosing to average over three images when grabbing (the Integrations option in the Camera Settings dialogue box).



A5.2 Summary of Results

Case 1: Standard Comet data file image, interactive Comet box fixed.

This case reproduced the same conditions for every measurement -the same image with the same interactive definitions of the Comet and background boxes. Not surprisingly the same results were obtained each time, so they are not reproduced here.

Case 2: Standard Comet data file, interactive Comet box redefined.

In this case the box containing the Comet was redefined between each measurement. Sometimes the box size was changed; sometimes its position was moved. The Comet box always fully contained the Comet; only the amount of background enclosed with the Comet was varied.

Name	Physical Area	Head DNA	Tail DNA	L/H
Min	33244	55.7661	42.4881	1.74032
Max	37260	57.5119	44.2339	1.96171
Mean	34838.32	56.8621	43.1379	1.834889
SD	1030.57	0.421149	0.421149	0.052862

The table below shows the parameter values for 50 sets of measurements:

Tail Moment	Tail Length	Comet Mean	Comet Mode	Comet Std. Dev.
37.6898	214.413	133	168.59	53.9968
42.1458	227.559	133	173.138	60.1916
39.46107	219.9273	133	170.4005	56.59249
1.051832	3.2798	0	1.074515	1.609348

Comet Radius	Comet Skew	Comet Kurtosis	Comet Variance	Comet Opt. Dens.
92.5997	2.84297	0.804464	32.0016	8004.58
98.1782	4.11351	1.10086	34.7652	8253.08
95.36046	3.44513	0.953734	33.20724	8097.371
1.335263	0.41391	0.095184	0.747578	60.66702



Head Mean	Head Mode	Head Std. Dev.	Head Radius	Head Skew
130	130	19.1569	58	0
130	130	19.249	62	0
130	130	19.21126	59.96	0
0	0	0.026954	1.261033	0

Case 2: Standard Comet data file, interactive Comet box redefined.

(contd.)

Head Kurtosis	Head Variance	Head Opt. Dens.	Tail Mean	Tail Mode
0	14.7361	4730.28	176	218.707
0	14.8069	4746.43	176	225.279
0	14.77789	4739.701	176	221.4636
0	0.020737	4.111566	0	1.63982

Tail Std. Dev.	Tail Radius	Tail Skew	Tail Kurtosis	Tail Variance
41.4826	125.707	2.71143	0.546978	18.9471
50.5136	132.279	4.72481	1.12124	22.4227
45.51865	128.4636	3.640203	0.839433	20.54679
2.50157	1.63982	0.666072	0.188034	0.983595

Tail Opt. Dens.
3502.71
3754.9
3596.16
60.96388



Case 3: Live grabbing of an image, interactive Comet box size fixed.

The Camera Settings dialogue parameters were used to obtain the best image of a Comet. Once this was established images were repeatedly grabbed under the same conditions using the Live/Store button in the Comet toolbar. Each time, the same fixed Comet box was used on the raw grabbed image for measurement.

The table below shows the parameter values for 75 sets of measurements. Some parameters are missing -a reminder to check what parameters are being recorded before you start!

Name	Physical Area	Head DNA	Tail DNA	L/H
Min	29805	85.8919	11.1881	1.2495
Max	29897	88.8119	14.1081	1.30886
Mean	29862.84	86.96873	13.03127	1.274574
Std. Dev.	21.37085	0.846886	0.846886	0.01527

Tail Length	Comet Std. Dev.	Comet Radius	Comet Skew	Comet Kurtosis
239.904	39.7602	101.114	3.07068	0.409898
253.919	42.086	101.986	3.30262	0.53217
245.5229	41.1532	101.5844	3.156709	0.45787
4.04759	0.710035	0.206929	0.071249	0.037422

Comet Opt. Dens.	Head Radius	Head Opt. Dens.	Tail Skew	Tail Kurtosis
7271.46	96	6387.7	5.541	-1.88718
8236.31	97	7428.73	9.45833	-1.11116
7892.204	96.31081	6971.181	6.942444	-1.38014
309.4808	0.465985	281.7465	1.347708	0.237614

Tail Opt. Dens.
881.977
1147.6
1044.368
76.58095



Case 4: Live grabbing of an image with averaging, interactive Comet box size fixed.

As Case 3 except that in this test each image measured was the average of three grabbed images. This is set up using the Integrations box in Camera Settings dialogue box. As before, once this was established images were repeatedly grabbed under the same conditions, using the Live/Store button in the Comet toolbar. Each time the same fixed Comet box was used on the averaged image for measurement.

The table below shows the parameter values for 75 sets of measurements:

Name	Physical Area	Head DNA	Tail DNA	L/H
Min	28969	90.3627	1.6752	1.23217
Max	29282	98.3248	9.63728	1.62508
Mean	29067.4	92.16932	7.830681	1.588514
Std. Dev.	42.03891	1.004191	1.004191	0.052815

Tail Moment	Tail Length	Comet Mean	Comet Mode	Comet Std. Dev.
0.968657	162.647	278	281.088	24.1959
7.36831	218.716	286	288.909	34.1881
6.12743	210.2369	283.52	287.5473	33.06493
0.801866	7.363791	2.896223	0.997541	1.180842

Comet Radius	Comet Skew	Comet Kurtosis	Comet Variance	Comet Opt. Dens.
60.0877	3.68791	0.530478	8.60797	527731
67.9093	4.14549	0.874687	11.8335	1500000
66.54732	3.833171	0.848555	11.49773	1373436
0.997599	0.073387	0.039085	0.379911	119205.2

Head Mean	Head Mode	Head Std. Dev.	Head Radius	Head Skew
281	281	23.2855	65	0
284	284	26.22	68	0
282.1733	282.1733	25.29016	66.17333	0
0.665224	0.665224	0.398288	0.665224	0



Case 4: Live grabbing of an image with averaging, interactive Comet box size fixed (contd.)

Head Kurtosis	Head Variance	Head Opt. Dens.	Tail Mean	Tail Mode
0	8.25726	562351	345	339.824
0	9.23241	1450000	349	364.358
0	8.962424	1330965	347.0267	360.2318
0	0.125288	107691.7	0.90005	2.859324

Tail Std. Dev.	Tail Radius	Tail Skew	Tail Kurtosis	Tail Variance
24.4721	104.824	2.74936	-1.95847	6.74929
43.572	138.358	10.3227	-0.98245	12.822
25.76361	132.1785	8.086101	-1.51472	7.156872
2.158124	5.246722	1.02159	0.19062	0.686396

Tail Opt. Dens.
9580.98
143437
114200.2
18224.61



APPENDIX 6 TROUBLESHOOTING

A6.1 GENERAL PROBLEMS

A6.1.1 No Image appears when I click on Live

Things to try....

Check the **beamsplitter** is in the camera position on the microscope. Check that the camera is getting **power**.

Check that a cell is in **focus** at the **centre** of the microscope **field of view**. Click on **Store** to Stop Grabbing.

Click on the Live button.

Depending on what type of camera you have, try some of the following. Check the **Contrast** and **Brightness** settings.

Check the **Offset** and **Gain** settings. Check **COM port** settings.

Check **Binning** for digital cameras. Check the **Exposure** Time.

Press the **AutoMap** button below the Image Window.

Help!

If you have installed the imaging card yourself then refer to the **Imaging Card Installation guide** in Appendix A. If this is not helpful or if no image appears, contact your supplier for further advice.

A6.1.2 No Image appears when I select On-CCD

If you are using a Matrox Meteor 2 imaging card, when you select On-CCD in the Camera Settings dialogue you may initially get a blank image in the Image Window. This problem can be easily resolved by moving the Exposure Time scroll bar. Once you have done this, the image will be redisplayed.

Note: If you set a high exposure time it may slow your system down significantly. It is recommended that you set the exposure time to as low a value as possible before you exit from Komet.



A6.1.3 Driver Problems

It is recommended that you install the latest **Windows Service Pack**, otherwise you may experience driver problems. This is included on the Komet Installation CD-ROM.

In the Windows Explorer double click on the file **Update.exe**, and follow the on-screen instructions. Reboot your computer following the installation.

A6.1.4 User Name System Set-up Lost

When you start your computer, on opening Windows you are asked to enter a

Windows User Name.

When you start Komet you are asked to enter a Komet UserID.

On exiting from Komet the current system set-up is saved to the Registry. This system set-up is specific to **both** the **Windows User Name** and the **Komet UserID**.

If you wish to start Komet with this system set-up at a future date you must ensure that you enter the same **Windows User Name** on starting Windows and the same **Komet UserID** on starting Komet. E.g. If you log on to Windows with the User Name "Administrator" and you then start Komet with the UserID "Genghis Khan", to reload the system set-up for a future session of Komet you must again log on to Windows with the User Name "Administrator" and then start Komet with the UserID "Genghis Khan". If either the Windows User Name or the Komet UserID is not identical to their previous values the system set-up will not be reloaded and Komet will start using default values.

A6.1.5 Saturation Detection with the Meteor or Meteor 2

The Matrox Meteor full range is limited to 254, not 255. To detect the camera saturation, you should therefore set the Saturation Maximum parameter in the Comet Options dialogue to 254.

A6.1.6 Regions Resizing and Disappearing

If running Komet with the Hamamatsu Orca 1 camera, when you resize or move regions on the image you may see the region disappear, or grow extremely large. If this happens, open the Comet Options page and edit the value for the Saturation Maximum to 4095.

This may be caused by a non 12-bit image being opened in the Image Window.